


# Operating Manual

## Cryo-Compact Circulators

### The *HighTech* Series

CF31

CF41

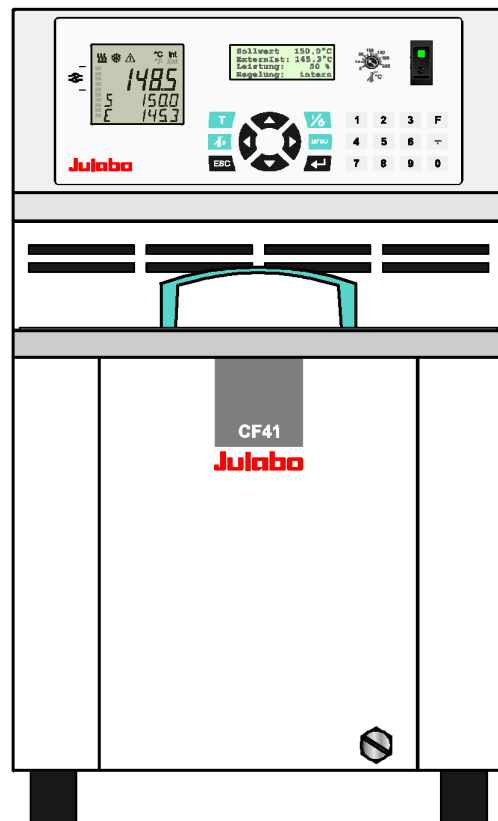

Distributed by:



ADVANCED APPLIED TECHNOLOGIES

Contact Us:

Irl Ph: 01 4523432

UK Ph: 08452 30 40 30

Web: [www.carlstuart.com](http://www.carlstuart.com)

Email: [info@carlstuart.com](mailto:info@carlstuart.com)

Printed in Germany

Changes without prior notification reserved

1.951.4871 BE0 01/07

**Julabo**  
Innovative Temperature Technology

JULABO Labortechnik GmbH  
77960 Seelbach / Germany

☎ +49 (0) 7823 / 51-0

📠 +49 (0) 7823 / 24 91

✉ [info@julabo.de](mailto:info@julabo.de)

🌐 [www.julabo.de](http://www.julabo.de)

---

## Congratulations!

You have made an excellent choice.

JULABO thanks you for the trust you have placed in us.

This operating manual has been designed to help you gain an understanding of the principles of operating and possibilities of our Cryo-Compact Circulators. For optimum utilization of all functions, we recommend that you thoroughly study this manual prior to beginning operation.

## Quality Management System



### The JULABO Quality Management System:

Development, production and distribution of temperature application instruments for research and industries conform to the requirements according to DIN EN ISO 9001:2000.

Certificate Registration No. 01 100044846

## EC Conformity



The products described in the operating instructions conform to the requirements of the following European guidelines:




Low voltage regulations with respect to legal harmonization of the member countries concerning electric devices for use within certain voltage limits.

EMC guideline with respect to legal harmonization of the member countries concerning electromagnetic compatibility.

**Julabo**

JULABO Labortechnik GmbH  
Eisenbahnstr. 45  
77960 Seelbach / Germany

## TABLE OF CONTENTS

EC Conformity .....	2
1. Use according to intended purpose.....	5
2. Operator responsibility – Safety recommendations .....	6
2.1. Handling .....	6
2.2. Use.....	6
2.3. Disposal .....	7
3. Technical specifications .....	8
3.1. Warning functions and safety installations .....	10
4. Safety notes for the user.....	11
4.1. Safety recommendations.....	11
5. Unpacking and checking .....	12
6. Installation .....	12
6.1. Temperature application to external systems .....	13
6.2. Tubing .....	15
7. Operating controls and functional elements .....	16
8. Operating procedures .....	20
8.1. Bath fluids .....	20
8.2. Power connection.....	21
8.3. Filling.....	22
8.4. Switching on / Selecting the language .....	23
9. Manual operation .....	23
9.1. Start - Stop .....	23
9.2.  Setting the temperatures .....	24
9.3.  Safety installations, warning functions .....	25
9.3.1. Excess temperature protection.....	25
9.3.2. Early warning system, low level protection .....	28
10.  Menu functions.....	29
10.1. Configuration.....	30
10.2. Control .....	34
10.3. Start of a profile .....	38
10.3.1. Edit after start.....	40

---


10.3.2. Interrupting a profile.....	42
10.3.3. Interruption after a power failure .....	42
10.3.4. Termination of a profile .....	43
10.4. Integrated programmer .....	44
10.5. Analog inputs/outputs .....	47
10.5.1. REG+E-PROG – temperature recorder outputs .....	49
10.5.2. REG+E-PROG - programmer input.....	50
10.5.3. STAND-BY input.....	52
10.5.4. ALARM output.....	53
10.6. Limits.....	54
10.7. Serial interface.....	56
10.8. ATC Absolute Temperature Calibration, 3-point calibration.....	57
10.9. Setting the pump pressure.....	60
11. Troubleshooting guide / Error messages.....	61
12. Electrical connections.....	64
13. Remote control.....	66
13.1. Setup for remote control .....	66
13.2. Communication with a PC or a superordinated data system .....	66
13.3. List of commands.....	67
13.4. Status messages .....	70
13.5. Error messages .....	71
14. JULABO Service – Online remote diagnosis .....	72
15. Installation of electronic module .....	73
16. Maintaining the cooling performance.....	73
16.1. Draining .....	74
17. Cleaning / repairing the unit.....	75
18. Warranty conditions .....	76

## 1. Use according to intended purpose

JULABO Cryo-Compact Circulators have been designed for temperature application to specific fluids in a bath tank. The units provide pump nozzles for temperature application to an external system (loop circuit).



Setpoint : 150.00°C  
 InternRef : 140.03°C  
 Power : 60%  
 Control : Intern

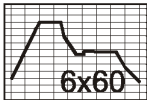



**ICC**

**TCF**

**ATC<sup>3</sup>**

**RS232**  
**RS485**

**Pt100**

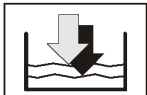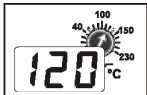

**SMART PUMP**

- ☑ The Cryo-Compact circulators are operated via the splash-proof keypad. The implemented microprocessor technology allows to set and to store different values that can be indicated on the VFD COMFORT-DISPLAY and LCD DIALOG-DISPLAY. Three menu keys facilitate adjusting setpoints, warning and safety functions and menu functions.
- ☑ The integrated programmer allows storing and running temperature and time-dependent processes.
- ☑ The control electronics including “ICC - Intelligent Cascade Control” automatically adapts the heat supplied to the thermal requirements of the bath.
- ☑ The TCF - Temperature Control Features allow the user to have access to all important temperature control parameters. This means: Full control on the control mode and the chance to manually adjust or adapt control to the specific application.
- ☑ Absolute Temperature Calibration (ATC3). With the 3-point calibration an offset is adjusted at three temperatures to ensure an accurate temperature pattern at the selected spot in the bath over the full temperature range.
- ☑ Electrical connections:  
 The serial interface, switchable from RS232 to RS485, allows modern process technology without additional interface.  
 Connection for Pt100 external sensor for external temperature measurement and control.  
 Alarm output for external alarm message or control of JULABO refrigerating baths or solenoid valve (cooling water).  
 The electronic module (option) provides 3 further analog connections (alarm input, standby input, recorder output, programmer input).
- ☑ The excess temperature protection conforming to IEC 61010-2-010 is a safety installation independent from the control circuit. This protection can be indicated and set on the LCD DIALOG-DISPLAY.
- ☑ The early warning system for low level signals that bath fluid needs to be refilled before the low level protection conforming to IEC 61010-2-010 causes a complete shut-down of the main functional elements.
- ☑ The pump capacity (electronically adjustable via the motor speed) enables to adapt to varying conditions for internal and external temperature applications..



JULABO Cryo-Compact circulators are not conceived for direct temperature application to food and luxury articles or pharmaceutical and medico-technical products. Direct temperature application means: Unprotected contact of the object with the bath medium (bath fluid).

## 2. Operator responsibility – Safety recommendations

The products of JULABO Labortechnik GmbH warrant a safe operation if installation, operation and maintenance is carried out according to common safety regulations. This section informs you about potential dangers that may arise from operating the Cryo-Compact Circulator and also mentions the most important safety precautions.

The operator is responsible for the qualification of the personnel operating the units.

The operator should be constantly informed about the dangers involved with their job activities as well as preventive actions.

Make sure all persons expected to carry out operation, installation and maintenance of the unit read and understand the safety information and operating instructions.

When using hazardous materials, the Cryo-Compact Circulator may only be operated by persons that are absolutely familiar with these materials and the Cryo-Compact Circulator. These persons must be fully aware of possible risks.

If you have any questions concerning the operation of your unit or the information in this manual, please contact us!

<b>Contact</b>	JULABO Labortechnik GmbH Eisenbahnstraße 45 77960 Seelbach / Germany	☎ +49 7823 51-0 ☎ +49 7823 2491	✉ <a href="mailto:info@julabo.de">info@julabo.de</a> 🌐 <a href="http://www.julabo.de">www.julabo.de</a>
----------------	--	------------------------------------	--

### 2.1. Handling

You received a product conceived for industrial use. Nevertheless, avoid strikes to the housing, vibrations, damages to the keypad foil (keys, display) or contamination.

Make sure the product is regularly checked for proper condition. Regularly check (at least every 2 years) the proper condition of the mandatory, warning, prohibition and safety labels.

Take care that the mains supply features a low impedance to avoid any negative affects on the instrument being operated in the same mains.

This unit is designed for operation in a controlled electromagnetic environment. This means that transmitting devices (e.g. cellular phones) should not be used in the immediate vicinity.

Magnetic radiation may influence other units with components susceptible to magnetic fields (e.g. a monitor). We recommend to keep a minimum distance of 1 m.

Permissible ambient temperature: max. 40 °C, min. 5 °C.

Permissible relative air humidity: 50 % (40 °C).

Do not store in an aggressive atmosphere. Protect from contaminations. Do not expose to sunlight.

#### Appropriate Operation

Only qualified personnel is authorized to perform configuration, installation, maintenance and repairs of the recirculating cooler. Untrained personnel should be instructed by trained personnel.

### 2.2. Use




Only use recommended materials (bath fluids). Only use non-acid and non corroding bath fluids.

The bath can be filled with flammable materials. Fire hazard!



There might be chemical dangers depending on the bath medium used.

Observe all warnings for the used materials (bath fluids) and the respective instructions (safety data sheets). Insufficient ventilation may result in the formation of explosive mixtures. Only use the unit in well ventilated areas (see page 12).

When using hazardous materials, **the user must** attach the enclosed safety labels to the front of the unit so they are well visible:

Warning label W09: Colors: yellow, black		Danger area. Attention! Observe instructions. (operating manual, safety data sheet)
Mandatory label M018 Colors: blue, white		Carefully read the user information prior to beginning operation Scope: EU
Semi S1-0701 Table A1-2 #9		Carefully read the user information prior to beginning operation Scope: NAFTA

Particular care and attention is necessary because of the wide operating range. There are thermal dangers: Burn, scald, hot steam, hot parts or cold parts and surfaces that can be touched.

Warning label W26: Colors: yellow, black		Hot surface warning. (The label is put on by JULABO)
Warning label W017: Colors: yellow, black		Low temperature warning. (The label is put on by JULABO)

Observe the instructions in the manuals for instruments of a different make that you connect to the Cryo-Compact Circulator, particularly the respective safety recommendations. Also observe the pin assignment of plugs and technical specifications of the products.


### 2.3. Disposal

The Cryo-Compact Circulator contains a so-called back-up battery that supplies voltage to memory chips when the unit is switched off. Do not dispose of the battery in domestic waste!

Depending on battery regulations in your country, you might be obliged to give back used or defect batteries to gathering places.

The product may be used with oil as bath fluid. These oils fully or partially consist of mineral oil or synthetic oil. For disposal, observe the instructions in the safety data sheets.

This unit contains the refrigerants R404A or R134a – at this time considered not to have any negative effects on the ozone layer. However, during the long operating period of the unit, disposal prescriptions may change. So only qualified personnel should take care of disposal.

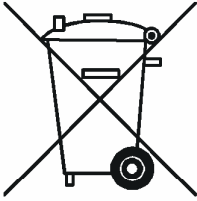	<p><u>Valid in EU countries</u>            Directive 2002/96/EC of the European Parliament and of the Council of 27 January 2003 on waste electrical and electronic equipment (WEEE).            This directive requires electrical and electronic equipment marked with a crossed-out trash can to be disposed of separately in an environmentally friendly manner. Contact an authorized waste management company in your country.            Disposal with household waste (unsorted waste) or similar collections of municipal waste is not permitted!</p>
---	--

### 3. Technical specifications

		<b>CF31</b>	
Working temperature range	°C	-30 ... +200	
Temperature stability	°C	±0.02	
Temperature selection		digital	
via keypad		indication on LCD DIALOG-DISPLAY (°C/°F)	
remote control via personal computer		indication on monitor	
Temperature indication:		VFD COMFORT-DISPLAY	
Resolution	°C	0.01	
Pump pressure stage		indication on VFD COMFORT-DISPLAY	
Temperature control		ICC - Intelligent Cascade Control	
Working temperature sensor		Pt 100	
Safety temperature sensor		Pt 100	
Absolute Temperature Calibration			
at internal control	°C	±3	
at external control	°C	±10	
Heater wattage (at 230 V)	kW	2,0	
Heater wattage (at 115 V)	kW	1,0	
Cooling capacity	°C	<u>20    0    -20</u>	
Medium ethanol	kW	0.32   0.25   0.15	
Cooling compressor		1-stage	
Refrigerant		R134a	
Cooling machine		Air-cooled	
Electrical connections: see page 10			
Electronically adj. pump capacity		1 ... 4 stages	
Flow rate min ... max. at 0 bar	l/min	22 ... 26	
Pressure max. at 0 liter	bar	0,7	
Suction max at 0 liter	bar	0,4	
Bath opening (WxL)	cm	16x3	
Bath depth	cm	14	
Filling volume	liters	2.0 ... 3.0	
Overall dimensions (WxDxH)	cm	24x46x40	
Weight	kg	33	
Ambient temperature	°C	5 ... 40	
Protection class according to IEC 60 529		IP 21	
Mains power connection 230 V/50 Hz	V/ Hz	207-253 / 50	
Current input (at 230 V)	A	11	
Mains power connection 230 V/60 Hz	V/ Hz	197-242/60	
Current input (at 230 V)	A	12	
Mains power connection 115 V/60 Hz	V/ Hz	103-127 / 60	
Current input (at 115 V)	A	14	

All measurements have been carried out at: rated voltage and frequency ambient temperature: 20 °C  
 Technical changes without prior notification reserved.



		<b>CF41</b>	
Working temperature range	°C	-40 ... +200	
Temperature stability	°C	±0.02	
Temperature selection via keypad		digital indication on LCD DIALOG-DISPLAY (°C/°F)	
remote control via personal computer		indication on monitor	
Temperature indication:		VFD COMFORT-DISPLAY	
Resolution	°C	0.01	
Pump pressure stage		indication on VFD COMFORT-DISPLAY	
Temperature control		ICC - Intelligent Cascade Control	
Working temperature sensor		Pt 100	
Safety temperature sensor		Pt 100	
Absolute Temperature Calibration			
at internal control	°C	±3	
at external control	°C	±10	
Heater wattage (at 230 V)	kW	2,0	
Heater wattage (at 115 V)	kW	1,0	
Cooling capacity	°C	20 0 -20 -30	
Medium ethanol	kW	0.47 0.4 0.28 0.12	
Cooling compressor		1-stage	
Refrigerant		R404A	
Cooling machine		Air-cooled	
Electrical connections: see page 10			
Electronically adj. pump capacity		1 ... 4	
Flow rate min ... max. at 0 bar	l/min	22 ... 26	
Pressure max. at 0 liter	bar	0,7	
Suction max at 0 liter	bar	0,4	
Bath opening (WxL)	cm	19x3	
Bath depth	cm	19	
Filling volume	Liter	4.0 ... 5.5	
Overall dimensions (WxDxH)	cm	28x46x46	
Weight	kg	42	
Ambient temperature	°C	5 ... 40	
Protection class according to IEC 60 529		IP 21	
Mains power connection 230 V/50 Hz	V/ Hz	207-253 / 50	
Current input (at 230 V)	A	13	
Mains power connection 230 V/60 Hz	V/ Hz	197-242/60	
Current input (at 230 V)	A	13	
Mains power connection 115 V/60 Hz	V/ Hz	103-127 / 60	
Current input (at 115 V)	A	18	

Alle Daten beziehen sich auf Nennspannung und Nennfrequenz Umgebungstemperatur 20 °C  
Technische Änderungen vorbehalten.

## Technical specifications

---

### Electrical connections:

Computer interface	RS232 / RS485
External sensor	Pt100

<b>Optional:</b> (Order No. 8900100 Electronic module with analog connections)	
Programmer input -100 °C to 400 °C = 0 - 10 V or 0 - 20 mA or 4 - 20 mA	
Input for the signal of a flow meter or external manipulated variable	
Temperature recorder outputs	0 - 10 V (0 V = -100 °C, 10 V = 400 °C)
	0 - 20 mA (0 mA = -100 °C, 20 mA = 400 °C)
	4 - 20 mA (4 mA = -100 °C, 20 mA = 400 °C)
Standby input	for external emergency switch-off
Alarm output	for external alarm signal

### 3.1. Warning functions and safety installations

Excess temperature protection	adjustable from 0 °C ... 220 °C
Low liquid level protection	float switch
Early warning system for low level	float switch
Classification according to DIN 12876-1	class III
Alarm message	optical + audible (permanent)
Warning message	optical + audible (in intervals)
Overload protection	for compressor and pump motor
High temperature warning function	optical + audible (in intervals)
Low temperature warning function	optical + audible (in intervals)
Supervision of working sensor	plausibility control
Reciprocal sensor monitoring between working and safety sensors	difference >25 °C

#### Environmental conditions according to IEC 61 010-1:

Use only indoor.

Altitude up to 2000 m - normal zero.

Ambient temperature: +5 ... +40 °C (for storage and transportation)

Air humidity:

Max. rel. humidity 80 % for temperatures up to +31 °C,

linear decrease down to 50 % relative humidity at a temperature of +40 °C

Stromversorgung: entspricht der Schutzklasse I, VDE 0106 T1





Over voltage category II

Pollution degree 2




**Not for use in explosive atmosphere**

## 4. Safety notes for the user

	<p>In addition to the safety warnings listed above, warnings are posted throughout the manual. These warnings are designated by an exclamation mark inside an equilateral triangle. „Warning of a dangerous situation (Attention ! Please follow the documentation).“ The danger is described according to an alarm keyword. Read and follow these important instructions.</p>
	<p><b>Warning:</b> Describes a possibly highly dangerous situation. If this is not avoided, serious injury and danger to life could result.</p>
	<p><b>Caution:</b> Describes a possibly dangerous situation. If this is not avoided, slight or minor injuries could result. A warning of possible damage can also be contained in the text.</p>
	<p><b>Notice:</b> Describes a possibly harmful situation. If this is not avoided, the product or anything in its surroundings can be damaged.</p>

### 4.1. Safety recommendations

Follow the safety recommendations to prevent damage to persons or property. Further, the valid safety instructions for working places must be followed.

	<ul style="list-style-type: none"> <li>• Only connect the unit to a power socket with earthing contact (PE – protective earth)!</li> <li>• Place the instrument on an even surface on a pad made of <b>non-inflammable</b> material.</li> <li>• Do not stay in the area below the unit.</li> <li>• Make sure you read and understand all instructions and safety precautions listed in this manual before installing or operating your unit.</li> <li>• Observe the fire point of the bath medium used. The excess temperature protection should be set at least 25 °C below the fire point.</li> <li>• Never operate the unit without bath fluid in the bath.</li> <li>• Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the fluid.</li> <li>• Prevent water from penetrating into the hot bath oil.</li> <li>• Exercise caution when emptying hot bath fluids! Check the temperature of the bath fluid prior to draining (by switching the unit on for a short moment for example).</li> <li>• Employ suitable connecting tubing. Make sure that the tubes are securely attached.</li> <li>• Never operate damaged or leaking equipment.</li> <li>• Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit.</li> <li>• Always empty the bath before moving the unit.</li> <li>• Never operate equipment with damaged mains power cables.</li> </ul>
---	---



- Some parts of the bath cover and the pump connections may become extremely warm during continuous operation. Therefore, exercise particular caution when touching these parts.




- Some parts of the bath cover and the pump connections may become extremely cold during continuous operation. Therefore, exercise particular caution when touching these parts.

## 5. Unpacking and checking

Unpack the unit and accessories and check for damages incurred during transit. These should be reported to the responsible carrier, railway, or postal authority, and a request for a damage report should be made. These instructions must be followed fully for us to guarantee our full support of your claim for protecting against loss from concealed damage. The form required for filing such a claim will be provided by the carrier.

## 6. Installation

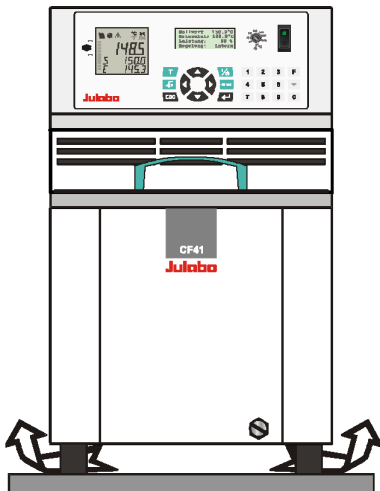

- Place the unit on an even surface on a pad made of **non-flammable** material.
- Cooling machine, pump motor and electronics produce intrinsic heat that is dissipated via the venting openings.! Never cover these openings!
- Be sure that the flow of ventilation can exit under the instrument.
- Keep at least 20 cm of open space on the front and rear venting grids
- The place of installation should be large enough and provide sufficient air ventilation to ensure the room does not warm up excessively because of the heat the instrument radiates to the environment. (Max. permissible ambient temperature: 40 °C). With regard to a disturbance in the cooling loop (leakage), the guideline EN 378 prescribes a certain room space to be available for each kg of refrigerant.
  - > For 0.48 kg of refrigerant R404A, a room space of 1 m<sup>3</sup> is required.
  - > For 0.25 kg of refrigerant R134a, a room space of 1 m<sup>3</sup> is required.Model CF41 with  
0.16 kg filling quantity of refrigerant R404A = 0,33 m<sup>3</sup> volume  
Model CF31 with  
0.15 kg filling quantity of refrigerant R134a = 0.6 m<sup>3</sup> volume

- Do not set up the unit in the immediate vicinity of heat sources and do not expose to sun light.
- Before operating the unit after transport, wait about one hour after setting it up. This will allow any oil that has accumulated laterally during transport to flow back down thus ensuring maximum cooling performance of the compressor.

The following questions shall help to recognize possible dangers and to reduce the risks to a minimum.

- Are all tubes and electrical cables connected and installed?  
Note: sharp edges, hot surfaces in operation, moving machine parts, etc.
- Do dangerous steams or gases arise when heating?  
Is an exhaust needed when working?
- What to do when a dangerous substance was spilled on or in the unit?  
Before starting to work, obtain information concerning the substance and determine the method of decontamination.

## 6.1. Temperature application to external systems

The Cryo-Compact Circulator is used for temperature application to external systems (loop circuit)



**Caution:** Securely attach all tubing to prevent slipping.



**Notice:** Flood hazard!


If the liquid levels in the Cryo-Compact Circulator bath and the external system are at different heights, overflowing must be prevented after the power has been turned off.

### Return flow safety device

For this reason, shut-off valves can be integrated in the loop circuit.

Order No.	Description
8 970 456	Shut-off valve (suitable up to +90 °C)
8 970 457	Shut-off valve (suitable up to +200 °C)


**Temperature application to external, closed systems**

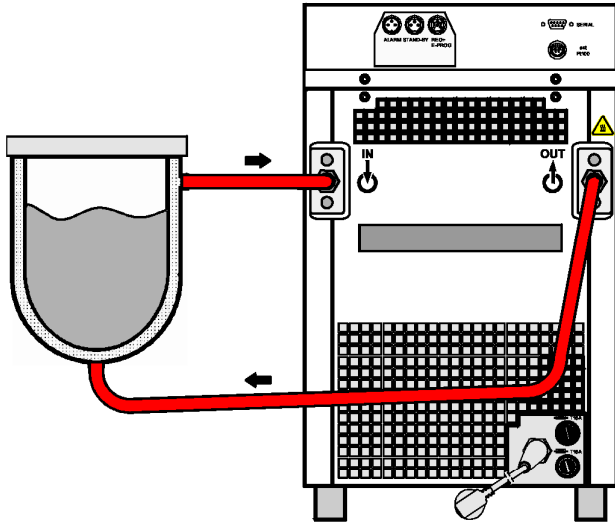

- Unscrew the M16x1 collar nuts on the pump connectors with a 19 mm (3/4") wrench and remove the sealing disks. Using the collar nuts, screw on the tubing connection fittings (for tubing 8 mm or 12 mm in diameter) delivered with the unit and tighten firmly. (Pressure pump: 14, suction pump: 15)
- Push on the tubings, and secure with tube clamps.
- Attach the tubing to the connectors of the external closed system, e.g., an instrument with a pressure-resistant temperature jacket or a temperature coil, and fasten with tube clamps to prevent slipping.

Tubing see page 15

**Temperature application to external, open systems**

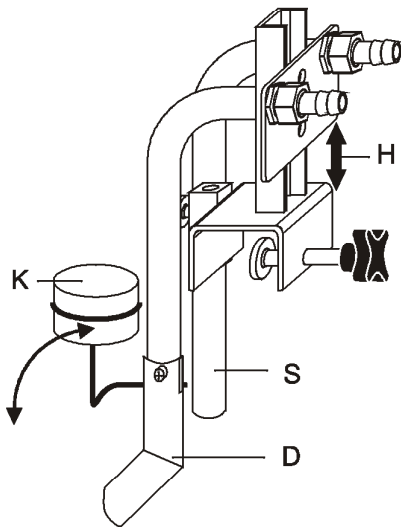

S = Suction pump connection  
 D = Pressure pump connection  
 K = Float  
 H = Height adjustment

The Cryo-Compact Circulator is equipped with both a pressure and suction pump for external temperature application in open systems. Differing flow rates of the pressure and suction pumps should be compensated. To maintain a constant liquid level, the JULABO „D+S“ Level Adapter is recommended for the external bath tank. The flow rate of the pressure pump will be then regulated by a built-in float device. The liquid level may be changed by a height adjustment on the „D+S“ Level Adapter.

Accessory: „D+S“ Level Adapter Order No. 8 970 410

**Important:**

- ⓘ The liquid level should be equal in the internal and external baths (absolute height).
- ⓘ If you take out samples (for example Erlenmeyer flasks) from the external bath, turn the Cryo-Compact Circulator off with the Start/Stop key.

## 6.2. Tubing



### **Warning:** Tubing:

At high working temperatures the tubing used for temperature application and cooling water supply represents a danger source.

A damaged tubing line may cause hot bath fluid to be pumped out within a short time.

This may result in:

- Burning of skin
- Difficulties in breathing due to hot atmosphere



### **Caution:**

Safety recommendations

- Employ suitable connecting tubing.
- Make sure that the tubing is securely attached.
- Avoid sharp bends in the tubing, and maintain a sufficient distance from surrounding walls.
- Regularly check the tubing for material defects (e.g. for cracks).
- Preventive maintenance: Replace the tubing from time to time.

### Recommended tubing:

Order No.		Suitable for
8930008	1 m CR <sup>®</sup> -tubing 8 mm inner dia. (-20 ... +120°C)	CF31, CF41
8930012	1 m CR <sup>®</sup> -tubing 12 mm inner dia. (-20 ... +120°C)	CF31, CF41
8930108	1 m Viton <sup>®</sup> tubing 8 mm inner dia (-50 °C bis 200 °C)	CF31, CF41
8930112	1 m Viton <sup>®</sup> tubing 12 mm inner dia (-50 °C bis 200 °C)	CF31, CF41

### Tubing insulation

8930410	1 m Insulation, 14 mm inner dia..	CR <sup>®</sup> -tubing 8 mm inner dia
8930412	1 m Insulation, 18 mm inner dia.	Viton <sup>®</sup> tubing 12 mm inner dia.

### Tube clamps

8970480	2 Tube clamps, size 1	CR <sup>®</sup> -tubing 8 mm inner dia
8970481	2 Tube clamps, size 2	Viton <sup>®</sup> tubing 12 mm inner dia.

### Metal tubing, flexible, triple insulated


8 930 209	0.5 m	2 fittings M16x1 female	-100 °C ... +350 °C
8 930 210	1.0 m		
8 930 211	1.5 m		
8 930 214	3.0 m		

### Metal tubing, flexible, insulated




8 930 220	0.5 m	2 fittings M16x1 female	-50 °C to +200 °C
8 930 221	1.0 m		
8 930 222	1.5 m		
8 930 223	3.0 m		

## 7. Operating controls and functional elements

Front view CF41 / CF31

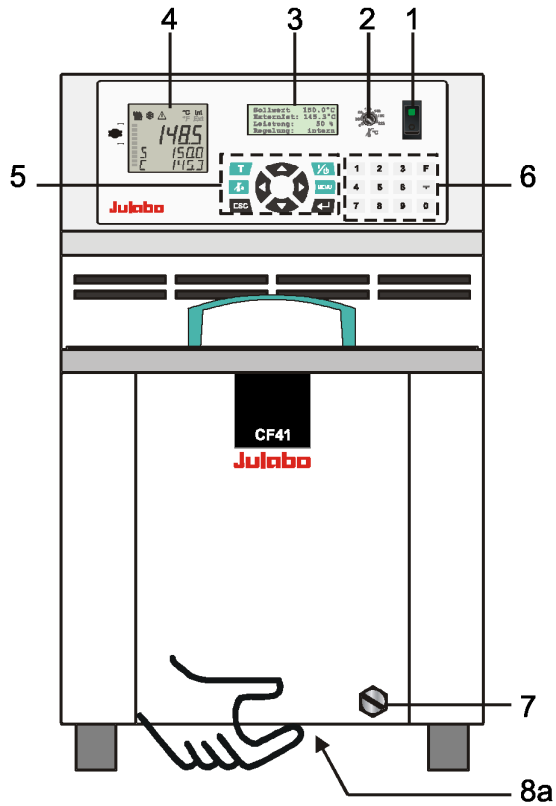

Rear view CF41/ CF31

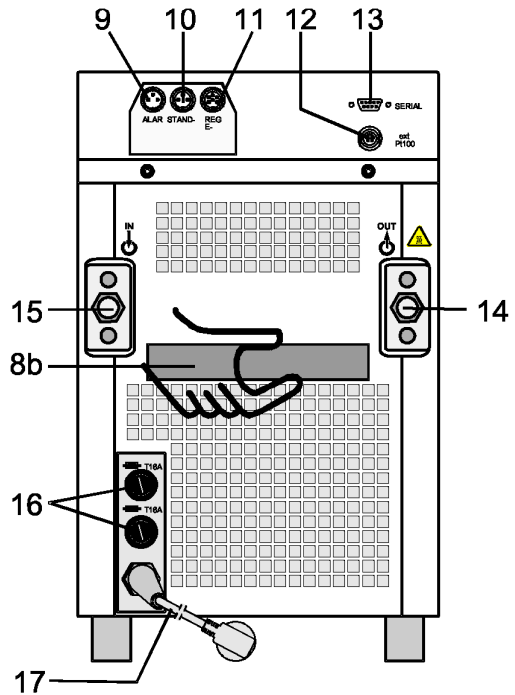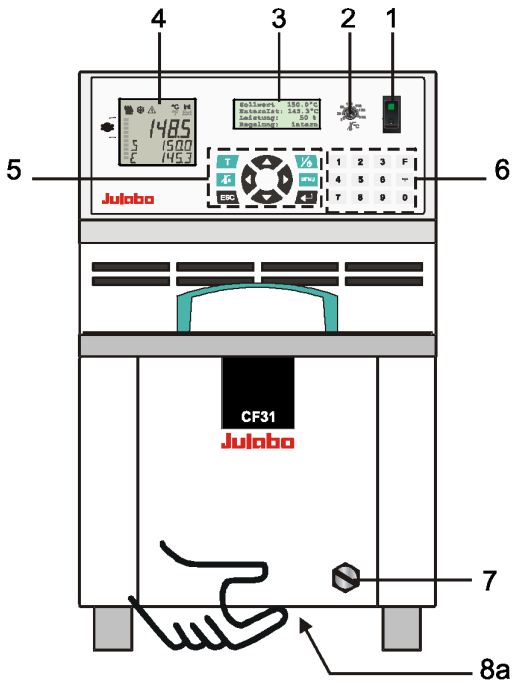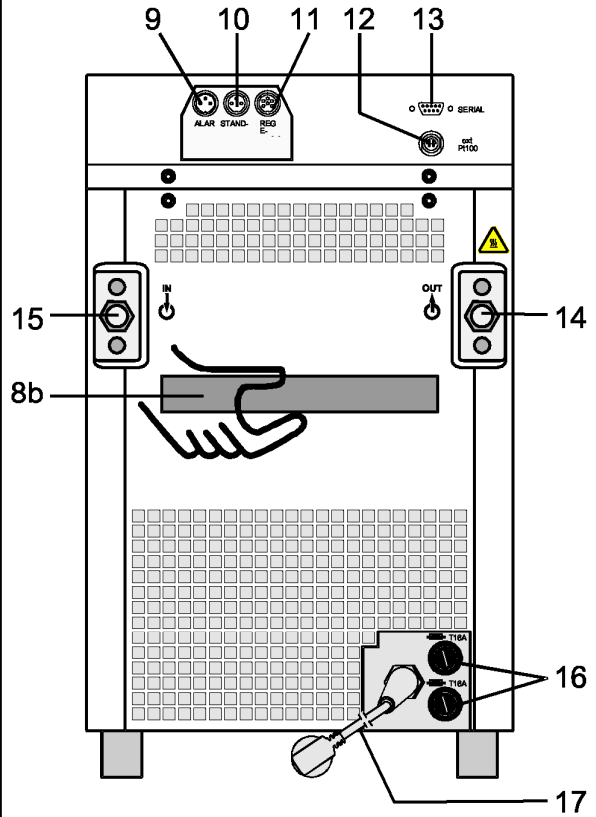





1

Mains power switch, illuminated

2



Adjustable excess temperature protection according to IEC 61010-2-010

3

Setpoint1:	150.00°C
IntAct :	148.53°C
Power :	80 %
Control :	intern

**LCD DIALOG-DISPLAY**

Line 1: Setpoint and origin of setpoint programming  
(Key / RS232 or RS485 / ext. Pt100 / EProg)


Line 2: Actual value - internal or external,  
identical to line 1 of the VFD-COMFORT-DISPLAY

Line 3: Heating capacity in %.

Line 4: Control type: internal / external

RS232 :	150.00°C
ExtAct :	145.30°C
Power :	80 %
Control :	extern

4

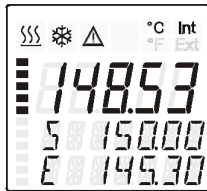

**VFD-COMFORT-DISPLAY**

Header: Control indicators see sections 4.1 and 4.2

Line 1: Actual value internal or external  
The display is depending on the selected control mode in the menu > Control < (internal or external).

Line 2: Working temp. setpoint, constantly S xxx.xx

Line 3: Actual value (E = external or I = internal)  
Alternating with the display in line 1

Use the keys ▼ ▲ to indicate further values in line 3. However, the functions of these keys are different with the programmer started.

- PI Capacity in % - with manipulated variable set to >control<\*
- PS Capacity in % - with manipulated variable set to >serial<\* or >eprog<\*

- H Heater capacity in Watts
- U Mains voltage Volts
- F Flow rate in liters/minute  
(providing EPROG input set to >Flowrate<)

\*see >Configuration> page 30

4.1



Control indicators in the header:

Heating / Cooling / Alarm /

**R**emote control

4.2




Control indicators in the header:

Temperature indication **I**nternal or **E**xternal actual value

Temperature indication in °C or °F (see page 33)

4.3

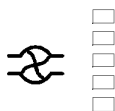

Display for the adjusted pump pressure stage in the -OFF- mode.

Display for the effective pump pressure stage (rotation speed) after start.

Four stages, adjustable via the  button, in the menu >PUMP<.

Operating controls and functional elements

5 Keypad 1



Start / stop key



Key for selecting the working temperature - Setpoint 1, 2, 3



Key for selecting the warning and safety values



MENU key - for selecting the menu functions



Cursor keys (left or right)



Edit keys (increase or decrease)



Enter key  
1) Store value / parameter  
2) Next lower menu level



Escape key  
1. Cancel entries  
2. Return to a higher menu level

6 Keypad 2



Numeral keypad: numerals 0 to 9



minus / decimal point



F key: Shows the help text in the LCD DIALOG-DISPLAY when alarm messages appear.




7 Drain port


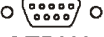




8a Handle: front



8b Handle: rear

Rear view

Option: Electronic module	Order No. 8 900 100
The Cryo-Compact Circulator automatically recognizes the connected electronic module.	
9	 Alarm output (for external alarm signal)
10	 Standby input (for external emergency switch-off)
11	 Programmer input and temperature recorder output
	REG+E-PROG

- 12  Socket for external measurement and control sensor  
or external setpoint programming  
ext Pt100
- 13  Interface RS232 / RS485: remote control via personal computer  
SERIAL
- 14 Pump connection  OUT - pressure pump, M16x1
- 15 Pump connection  IN – suction pump, M16x1 (return)
- 16  Mains fuses, T16A
- 17  Mains power cable with plug

## 8. Operating procedures

### 8.1. Bath fluids



**Caution:**

Carefully read the safety data sheet of the bath fluid used, particularly with regard to the fire point!

If a bath fluid with a fire point of  $\leq 65$  °C is used, only supervised operation is possible.

**Water:**

The quality of water depends on local conditions.

Ferrous water can cause corrosion - even on stainless steel.

Chloric water can cause pitting corrosion.

**Mixture water -glycol:**

Strictly observe the safety data and handling instructions from the manufacturer.

The proportion of water might evaporate by and by. Check the mixing ratio regularly and refill water if necessary.

**Recommended bath fluids:**

Bath fluids	Temperature range	Flash point	fire point
water	5 °C ... 80 °C		
mixture water/glycol	-30 °C ... 50 °C		
JULABO Thermal G	-30 °C ... 80 °C		
JULABO Thermal M	+40 °C ... +170 °C	>280 °C	>305 °C
JULABO Thermal H	+20 °C ... +250 °C	>270 °C	>360 °C
JULABO Thermal H5S	-50 °C ... +105 °C	>110 °C	>130 °C
JULABO Thermal H10S	-20 °C ... +180 °C	>175 °C	>210 °C
JULABO Thermal H20S	+0 °C ... +220 °C	>230 °C	>270 °C
JULABO Thermal H250	+60 °C ... +200 °C	>225 °C	>250 °C

Order No. 10 liters	Bath fluids	Order No. 5 liters
8 940 124	JULABO Thermal G	8 940 125
8 940 100	JULABO Thermal M	8 940 101
8 940 102	JULABO Thermal H	8 940 103
8 940 106	JULABO Thermal H5S	8 940 107
8 940 114	JULABO Thermal H10S	8 940 115
8 940 108	JULABO Thermal H20S	8 940 109
8 940 116	JULABO Thermal H250	8 940 117

**Notice:**

Please contact JULABO before using other than recommended bath fluids. JULABO takes no responsibility for damages caused by the selection of an unsuitable bath fluid.

Unsuitable bath fluids are liquids which e.g.

- are very highly viscous (much higher than 70 mm<sup>2</sup> /s at the respective working temperature)
- have corrosive characteristics or tend to cracking.

ATTENTION: The maximum permissible viscosity is 70 mm<sup>2</sup> /s.

**V Caution:**

The temperature controlling i.e. of fluids in a reactor constitutes normal Cryo-Compact Circulator practice.

We do not know which substances are contained within these vessels.

Many substances are:

- inflammable, easily ignited or explosive
- hazardous to health
- environmentally unsafe

i.e.: **dangerous**

**The user alone is responsible for the handling of these substances!**

## 8.2. Power connection

**Caution:**

Only connect the unit to a power socket with earthing contact (PE – protective earth)! We disclaim all liability for damage caused by incorrect line voltages!

Check to make sure that the line voltage matches the supply voltage specified on the identification plate.

### 8.3. Filling




**Notice:**

Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the liquid.

**Guideline:**

A volume change of 12 % per 100 °C temperature variation is to be considered.

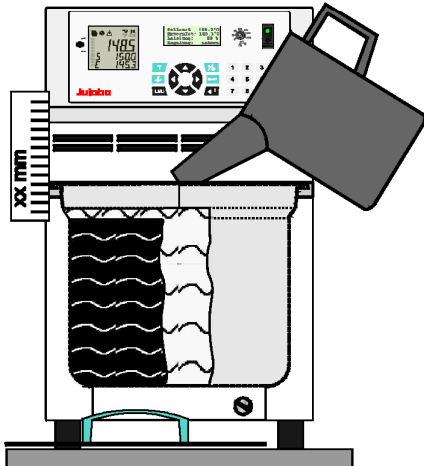

Take care that no liquid enters the interior of the Cryo-Compact Circulator .

- ① Connect the tubing from the external system to the pump connectors and check for leaks




Respect instructions from page 12 to page 15!

- ① Check to make sure that the drain tap (7) is closed.

**Recommendation:**


For filling, use for example an measuring jug with nuzzle.

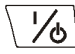- ① Recommended maximum filling level with water as bath fluid: 30 mm below the tank rim
- ① Recommended maximum filling level with bath oils: 40 mm below the tank rim
- Turn the mains switch (1) on (Switching on - see page 23)
- Switch on unit by pressing the start/stop key .
- Tempering fluid is pumped into the externally connected system. Refill fluid up to recommended filling level.

**Important:**

- ① When using a bath fluid, the change in volume in case of change in temperature has to be respected. Fill in a little amount of bath fluid only so that the low level alarm is not triggered.
- ① Low level alarm is triggered at the following liquid level:
  - CF31 75 to 80 mm below the tank rim
  - CF41 80 to 85 mm below the tank rim
- ① When reaching the working temperature, check the liquid level. If the cooling coil is not completely covered with bath fluid, refill it.

## 8.4. Switching on / Selecting the language

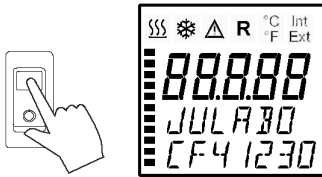

```
**** JULABO ****
*** HighTech ***
** Compact **
* Version 1.02-41 *
```

```
-OFF-
S 150.00
I 24.60
```

### Switching on:

- The Cryo-Compact Circulator is turned on and off with the mains switch. The integrated control light will illuminate to indicate that power has been applied.

**i** During the self-test all segments of the VFD-Info-Display and the LCD DIALOG-DISPLAY light up.

Then the type of unit and the software version is indicated.

(example: **JULABO CF41 230** Volt / **Version 1.02-41.**)

The display „OFF“ or „R OFF“ indicates the unit is ready to operate.

**i** The Cryo-Compact Circulator enters the operating mode activated before switching the Cryo-Compact Circulator off:

**keypad control mode** (manual operation) or

**remote control mode** (operation via personal computer).

### Selecting the language:

There are two options for the language of the LCD DIALOG-DISPLAY:

German or English. Select the desired language in the menu

>Configuration< under the submenu >Language/Sprache<.

```
Configuration
Off-Mode : pump off
>Language : english
Reset : no
```

Press the respective keys in the following order:



[▼ up to submenu >Language/Sprache <]



2x

## 9. Manual operation

### 9.1. Start - Stop

```
24.60
S 150.00
E -----
```

**Start:** Press the start/stop key .

The actual bath temperature is displayed on the VFD COMFORT-DISPLAY.

The LCD DIALOG-DISPLAY informs about adjustments and conditions on the Cryo-Compact Circulator (see example on the left).

**Stop:** Press the start/stop key .

The VFD COMFORT-DISPLAY indicates the message „OFF“.

**Autostart:** see chapter 10.1. Configuration

The Autostart function enables the start of the Cryo-Compact Circulator directly by pressing the mains switch or using a timer.

## 9.2. Setting the temperatures



Factory setting:

```
Setpoints
>Setpoint1: 20.00°C
Setpoint2: 37.00°C
Setpoint3: 70.00°C
```




Press the  key to call up the menu for temperature selection.

3 different working temperatures are adjustable. Their values are freely selectable within the operating temperature range.


 This setting may be carried out with the circulator being in the Start or Stop condition!



### Example: Setting working temperature „Setpoint 3“


```
Setpoints
Setpoint1: 20.00°C
Setpoint2: 37.00°C
>Setpoint3: 70.00°C
```

1. Press the  key.  
The LCD DIALOG-DISPLAY indicates the valid settings.
2. Use the keys   to select Setpoint 3.  
(One segment of the line blinks)

**Change the value to 85 °C.**

```
8 5 
Setpoints
Setpoint1: 20.00°C
Setpoint2: 37.00°C
>Setpoint3: 85.00°C
```

3. Use the numeral keypad to enter 8 and 5.  
Then press enter  to store the value.  
The value remains visible on the LCD DIALOG-DISPLAY for about 10 seconds, or press  to update the display immediately.

 In the >Start< condition this value is immediately used for controlling the working temperature.  
The indication on the VFD COMFORT-DISPLAY is updated.  
The heater control indicator blinks.



```
25.83
S 85.00
E -----
```




**Notice:** See **SetMax:** and **SetMin:**  
in chapter 10.6. Limits on page 54

### Example: Selecting the working temperature




- Press the  key. Use the keys   to select setpoint 1, 2 or 3 and press enter .

 The Cryo-Compact Circulator uses the new working temperature value for temperature control.




### 9.3. Safety installations, warning functions

#### 9.3.1. Excess temperature protection

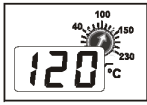



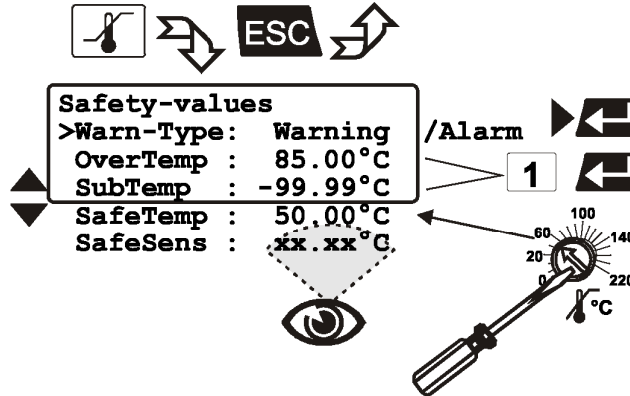

Settings for the excess temperature protection according to IEC 61010-2-010 and for the high and low temperature warning functions are made




in a menu that is called up with the key .





Legend:


<b>1</b>		Setting via keypad
		Select parameter
		Select menu item
		Change menu level scroll down / scroll up



1. Press the key .  
The LCD DIALOG-DISPLAY shows the valid settings (see example above).
2. Use the keys   to select the menu item.


**Settings:**

**Warn-Type:** Select the parameter with the key  and press enter  (warning or alarm).

**OverTemp and SubTemp:** Use the numeral keypad to enter the value and press enter .

**SafeTemp:** Set the new cut-out value using a screwdriver via the LCD DIALOG-DISPLAY. (Example: 50 °C)

Example:

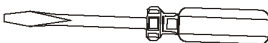

The value remains visible on the LCD DIALOG-DISPLAY for about 10 seconds, or press **ESC** to update the display immediately.

**Recommendation:**

Set the excess temperature protection >SafeTemp< at 5 to 10 °C above the working temperature setpoint.



**Warning:**



The excess temperature protection >SafeTemp< should be set at least 25 °C below the fire point of the bath fluid used.

In the event of wrong setting there is a fire hazard!

We disclaim all liability for damage caused by wrong settings!

**Warn-Type:** >Warning< or >Alarm<

For the two menu items >OverTemp< and >SubTemp< choose between a warning message being signaled or a complete shutdown of the main functional elements such as heater and circulating pump being effected.

**OverTemperature: SubTemperature:**

If for a sensitive temperature application task adherence to a working temperature value >Setpoint< is to be supervised, then set high and low temperature warning values.

In the example below, the >Setpoint< of 85 °C is surrounded by the values >OverTemp< 87 °C and >SubTemp< 83 °C. The electronics immediately registers when the actual temperature attains a temperature out of the limits and it follows a reaction according to what is set in the menu item >Warn-Type<.

**Warn-Type: >Warning<**

An audible warning (interval tone) sounds and a message appears on the VFD COMFORT-DISPLAY.

**OverTemp**

XXXXXX  
WARNING  
CODE 03

**SubTemp**

XXXXXX  
WARNING  
CODE 04

Press **F** for help on the LCD.

Actual temperature  
above high tempera-  
ture value.

Actual temperature  
below low tempera-  
ture value.

**Warn-Type: >Alarm<**

A complete shutdown of heater and circulating pump is effected. An audible alarm (continuous tone) sounds and a message appears on the VFD COMFORT-DISPLAY.

**OverTemp**

XXXXXX  
ALARM  
CODE 03


**SubTemp**

XXXXXX  
ALARM  
CODE 04

Press **F** for help on the LCD.

Actual temperature  
above high tempera-  
ture value.

Actual temperature  
below low tempera-  
ture value.

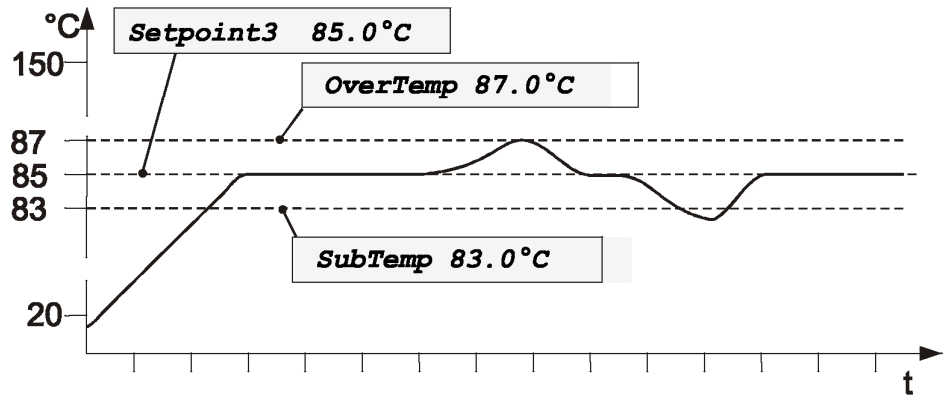

ⓘ The warning functions are only triggered when the actual bath temperature, after start from the „OFF“ or „rOFF“ mode, lies within the set limits for 3 seconds.

**SafeTemperature:** Indicated is the cut-out value set with a screwdriver on the excess temperature protection device.

Setting range:  
20 °C ... 220 °C

This safety installation is independent of the control circuit. When the temperature of the bath fluid has reached the safety temperature, a complete shutdown of the heater and pump is effected.

The alarm is indicated by optical and audible signals (continuous tone).

The following error message appears on the VFD COMFORT-DISPLAY:

```
XXXXXX
ALARM
CODE 14
```

Press **F** for help on the LCD

```
Working temperature
exceeds the adjusted
safety temperature.
```



**Notice:**


Check the safety installation at least twice a year!

- Excess temperature protection according to IEC 61010-2-010  
With a screwdriver turn back the adjustable excess temperature protection until the shut-down point (actual temperature).



**SafeSens:** Indicated is the temperature value of the safety sensor.

9.3.2. Early warning system, low level protection

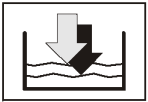

(patented)


This low level protection is independent of the control circuit and is divided in two sections.

1. Switch in stage 1 recognizes a critical fluid level 😊.

An audible warning (interval tone) sounds and a message appears on the VFD COMFORT-DISPLAY

<b>XXXXXX</b> <b>WARNING</b> <b>CODE 40</b>	Press <b>F</b> →	Niveau level warning Critical fluid level Replenish the bath tank with bath fluid
---	------------------	--

Refill bath fluid!

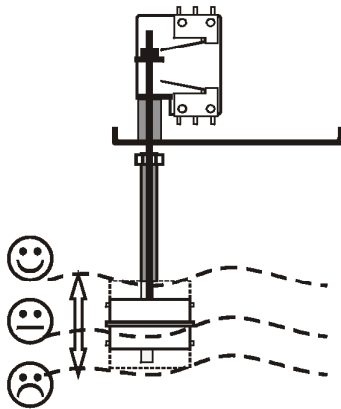

2. Switch in stage 2 recognizes a low fluid level 😞.

If stage 2 of the low level protection device (according to IEC 61010-2-010) is triggered, a complete shutdown of the heater and circulating pump is effected.

A continuous alarm tone sounds and a message >ALARM< >CODE 01< appears on the VFD COMFORT-DISPLAY.

<b>XXXXXX</b> <b>ALARM</b> <b>CODE 01</b>	Press <b>F</b> →	Insufficient liquid level or float is defective.
---	------------------	--

Turn off the unit with the mains switch, refill bath fluid and turn the unit on again!

**Important:** Check the safety installations from time to time.


	<p><b>Warning:</b></p> <p>For refill always use the same bath fluid type that is already in the bath.                  Bath oils must not contain any water contaminants!                  Explosion hazard at higher temperatures!</p>
--	---

	<p><b>Notice:</b></p> <p>Check the safety installation at least twice a year!                  Low level protection according to IEC 61010-2-010                  To check the function of the float, it can be manually lowered with a screwdriver for example.</p>
--	--

## 10. Menu functions


The term „menu functions“ refers to adjustments such as

- Configuration of the Cryo-Compact Circulator
- Internal and external control with adjustable control parameters
- Start menu for the integrated programmer
- Integrated programmer; 6 profiles with 60 sections each
- Configurable inputs and outputs; analog recorder outputs, input for external programmer or manipulated variable or flow rate
- Determination of limits for all important setting ranges and capacity variables
- Online communication, with adjustable interface parameters
- Absolute temperature calibration; sensor calibration
- Electronically adjustable pump capacity

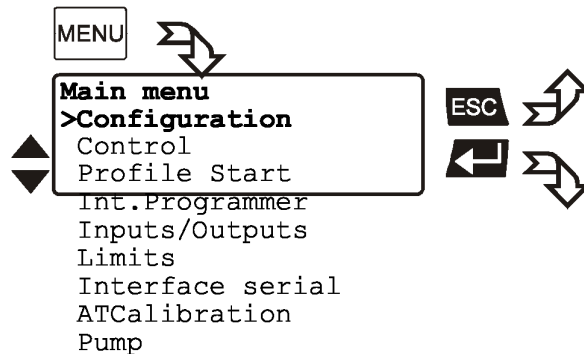


- i** If the electronic module is not fitted, the item >inputs/outputs< will be missing in the menu.

Legend

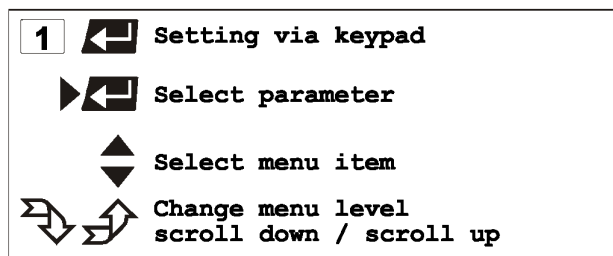

1. If you press  , use the keys   to move in menu level 1.
2. If the desired menu item is highlighted on the LCD DIALOG-DISPLAY (>), press enter  to change to menu level 2.

## 10.1. Configuration

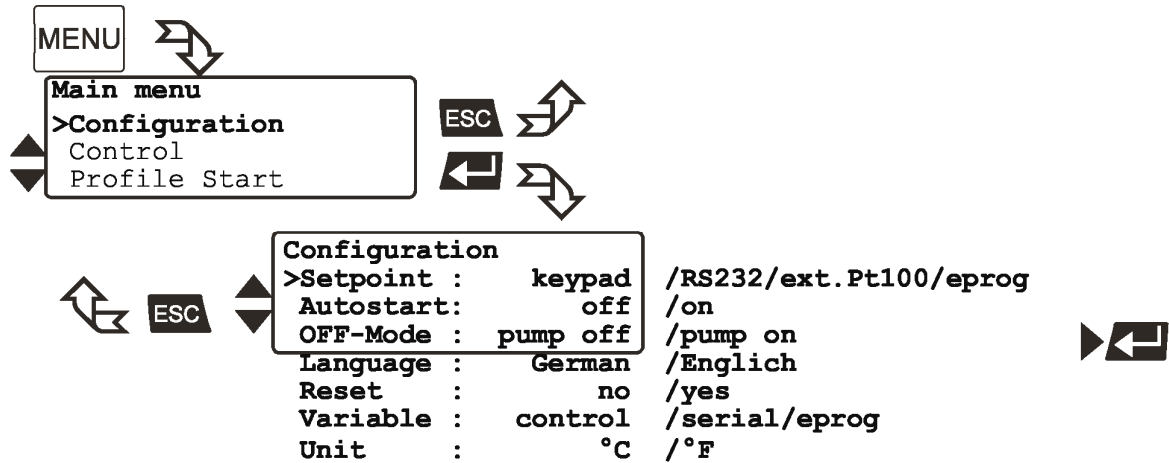

1. Use the up/down keys to select the desired menu item.
2. Use the left/right keys to select the parameter (parameter blinks).
3. Press enter to store the new parameter.

### Setpoint

The Cryo-Compact Circulator provides four possibilities for setpoint setting. The selected mode is indicated in line 1 on the LCD DIALOG-DISPLAY.

#### Possible parameters:

```
Setpoint1: 150.00°C
IntAct   : 148.53°C
Power    : 80 %
Control  : intern
```

**keypad** – Setpoint setting via keypad or via the integrated programmer.  
(factory setting)

```
RS232    : 150.00°C
ExtAct   : 145.30°C
Power    : 80 %
Control  : extern
```

**RS232 or RS485** - Setpoint setting via the serial RS232/RS485 interface through a PC or superordinated data system.

In the header of the VFD COMFORT-DISPLAY an „R“ illuminates. It indicates that remote control mode is set.



**i** RS232 or RS485: see menu > Serial interface < page 56.

```
-OFF-
S  xxxx
I  xxxx
```

#### **Important:**

Connect the Cryo-Compact Circulator to a PC using an interface cable. Check the interface parameters of both interfaces (on Cryo-Compact Circulator and PC) and make sure they match.  
(see 13.1. on page 66 )

```
Pt100    : 150.00°C
ExtAct   : 145.30°C
Power    : 80 %
Control  : extern
```

**ext.Pt100** – Setpoint setting via the analog socket „ext. Pt100“ using an external temperature sensor or an appropriate voltage/current source.

<b>eprog</b>	:	50.0°C
<b>IntAct</b>	:	24.7°C
<b>Power</b>	:	80 %
<b>Control</b>	:	intern

**eprog** – Can only be adjusted when an electronic module with analog connections is used (option).  
Setpoint setting via the analog interface REG+E-PROG through an external programmer.

**Important:**

Connect the external programmer to the Cryo-Compact Circulator via the socket REG+E-PROG.

In the menu >Inputs/Outputs< set the parameter >EPROG< and the input variables (see page 50).

The E-Prog input can only be used either under menu item >**Setpoint**< or under menu item >**Variable**< (page 32).

### Autostart

Possible parameters:

**on** - AUTOSTART on

**off** - AUTOSTART off (Factory setting)

**Note:**

The temperature system has been configured and supplied by JULABO according to N.A.M.U.R. recommendations. This means for the start mode, that the unit must enter a safe operating state after a power failure (non-automatic start mode). This safe operating state is indicated by „ **OFF** „, resp. **R OFF** on the VFD-COMFORT-DISPLAY. A complete shutdown of the main functional elements such as heater and circulating pump is effected simultaneously.

The values set on the Cryo-Compact Circulator remain stored, and the unit is returned to operation by pressing the start/stop key (in manual control mode). In remote control mode, the values need to be resent by the PC via the interface.

Should such a safety standard not be required, the AUTOSTART function (automatic start mode) may be activated, thus allowing the start of the instrument directly by pressing the mains power switch or using a timer.

ⓘ The AUTOSTART function can only be used, if setpoint setting is carried out via >keyboard< or >eprog< or >ext. Pt100<.



**Warning:**

For supervised or unsupervised operation with the AUTOSTART function, avoid any hazardous situation to persons or property.

The instrument does no longer conform to N.A.M.U.R. recommendations.

The safety and warning functions of the instrument should always be used to their fullest capacity.

### **OFF-Mode**

Normally the circulating pump is switched via the start/stop signal. However, if circulation should be maintained also for the -OFF- condition, the parameter **>pump on<** needs to be set.

Possible parameters:

**pump off** (factory setting)

**pump on**

### **Language**

There are two options for the language of the LCD DIALOG-DISPLAY: German or English.

Possible parameters:

**deutsch / englisch German / English**

### **Reset**

Use this to reset all values to factory setting (except date and time).

Possible parameters:

**yes**

**no** (factory setting)

### **Variable** - actuating variable

The variable corresponds to the extent to which the heater or cooling machine of the Cryo-Compact Circulator is controlled. Heat or cold is applied to the bath according to this variable. If this happens with the control electronics of the Cryo-Compact Circulator, called **>control<** in this particular case, the bath temperature is exactly heated and maintained constant at the adjusted setpoint.

① Programming of variables for the parameters **> serial <** or **> eprog <** is only accepted, if the unit is in Start mode.

Possible parameters:

**control** – The internal control electronics of the Cryo-Compact Circulator controls the heater and cooling machine. Self-tuning is possible.

**serial** – The heater or cooling machine receives the control signal via the serial interface. Self-tuning is not possible.

**eprog** - The heater or cooling machine receives the control signal via the E-Prog input. Self-tuning is not possible. (option).

#### **Important:**

Set the parameter **>EPROG<** and the input variables also in the menu **>Inputs/Outputs<** (see page 50).

#### **Note:**

The E-Prog input can only be used either under menu item **>Setpoint<** (page 30) or under menu item **>Variable<** (see above).



**Warning:**

The working temperature range of the Cryo-Compact Circulator is determined during configuration. If set to >Control<, this range cannot be exceeded.

If set to > **serial** < and > **eprog** <, heat or cold is applied to the bath without control.

The permissible maximum temperature can be exceeded. The user has to take adequate precautions for temperature control.

Materials, such as gaskets or insulations for example, may be damaged or destroyed, if the permissible maximum temperature is exceeded.

The safety and warning functions of the instrument should always be used to their fullest capacity. (See page 25)

**Unit**

Temperature values can be displayed in the unit °C or °F.

With this change all adjusted temperature values are converted and displayed into the new unit.

Example:

```
Setpoints
>Setpoint1: 20.00°C
Setpoint2: 37.00°C
Sezpoint3: 70.00°C
```

°C → °F ↓

```
Setpoints
>Setpoint1: 68.00°F
Setpoint2: 98.60°F
Setpoint3: 158.00°F
```



°F

```
-OFF-
S  XXXX
I  XXXX
```


Possible parameters:

°C (factory setting)

°F

## 10.2. Control

The Cryo-Compact Circulator is conceived for internal and external temperature control. Switching is carried out in this submenu. Depending on what is set, only the respective set of parameters is indicated.

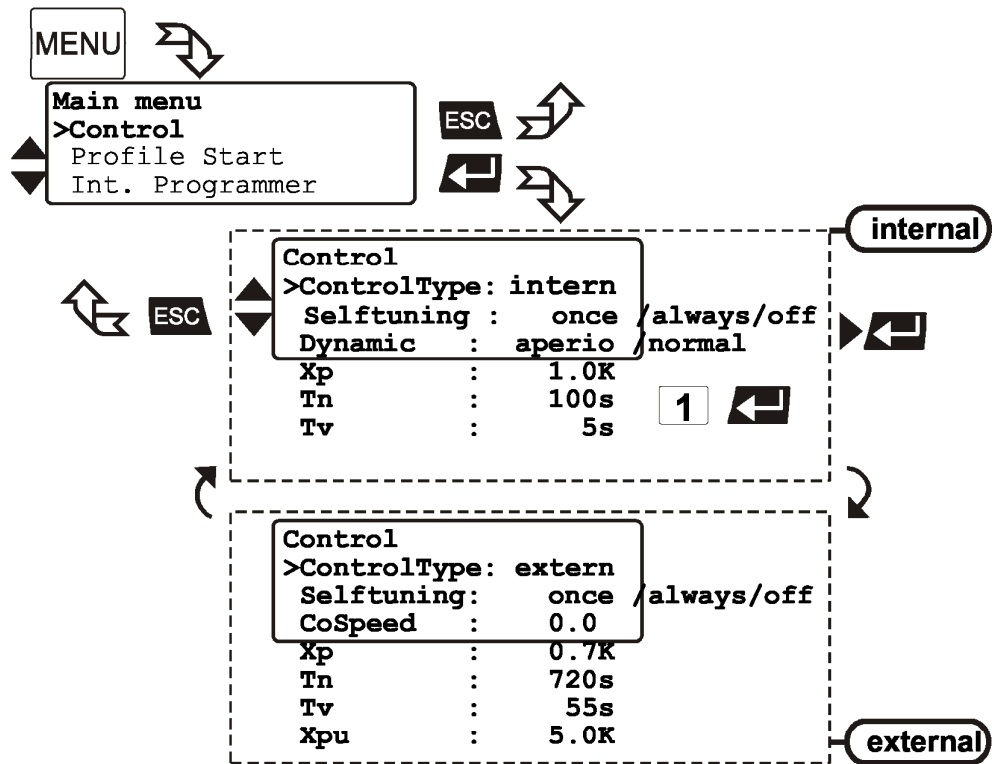

1. Use the keys to select the menu item.
2. With the keys select the parameter (parameter blinks)  
or  
use the numeral keypad to enter the numbers (number segment blinks).
3. Press enter to store the new parameter or value.

Example: 0.8




```

-OFF-
S xx.xx
E xx.xx
    
```

**ControlType:** Can only be adjusted in the -OFF- condition

Possible parameters:

**intern** internal temperature control  
**extern** external temperature control with external Pt100 sensor

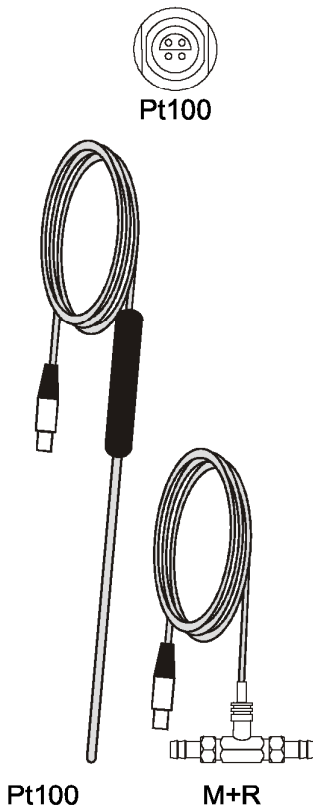**IMPORTANT:**

Additional measures for external temperature control.

- ❗ For external control and temperature measurement an external Pt100 sensor must be connected to the socket (7) on the rear of the Cryo-Compact Circulator .

Order No.	Description	Material	cable
8981003	200 x 6 mm dia.,	stainless steel	1.5 m
8981005	200 x 6 mm dia.,	glass	1.5 m
8981006	20 x 2 mm dia.,	stainless steel	1.5 m
8981010	300 x 6 mm dia.,	stainless steel	1.5 m
8981013	600 x 6 mm dia.,	stainless steel/Teflon	3 m
8981014	1200 x 6 mm dia.,	stainless steel/Teflon	3 m
8981103	Extension cable for Pt100 sensor		3.5 m
8981020	M+R in-line Pt100 sensor		

- ❗ The M+R in-line Pt100 sensor is a flow sensor and can be installed in the loop circuit.
- ❗ Sensor calibration of the external Pt100 sensor is carried out in the menu >ATCalibration<, with the > Status< set to >no< (see page 57).
- ❗ Suggested adjustments for external temperature control: AreaUp/ AreaDown and IntMax / IntMin see chapter >Limits< on page 54.
- ❗ External temperature control does not allow for setpoint setting via the socket „ext. Pt100“ (see page 30).

**Notice:**

Place the external sensor into the bath medium and securely fix the sensor.


**Selftuning:**

When performing an selftuning for the controlled system (temperature application system), the control parameters Xp, Tn and Tv are automatically determined and stored.

**Possible parameters:****off - no selftuning.**

The control parameters ascertained during the last identification are used for control purposes.

**once – single selftuning (factory setting)**

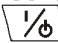The instrument performs a single selftuning of the controlled system after each start with the start/stop key  or after receiving a start command via the interface.


**always – continual selftuning**

The instrument performs a selftuning of the controlled system whenever a new setpoint is to be reached.

Use this setting only when the temperature application system changes permanently.

**Dynamic:**


This parameter affects the temperature pattern only in case of **internal** control.

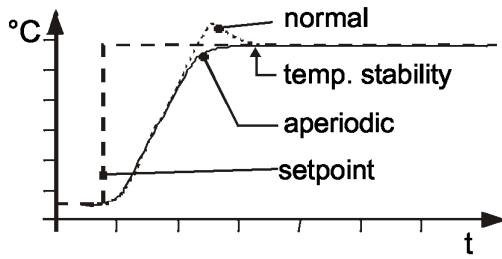

**Possible parameters:**

**normal** –Allows for reaching the setpoint faster, but overshooting of up to 5 % is possible.

**aperiodic** – (factory setting)  
Target temperature is attained without overshooting.

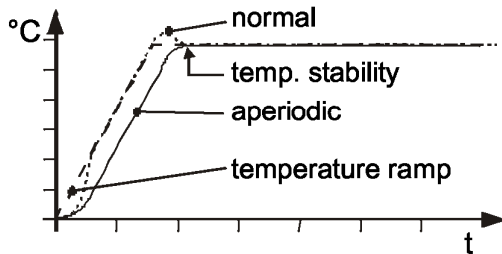


① With both adjustments an adequate temperature stability is reached after approximately the same time.

**CoSpeed:**

This parameter affects the temperature pattern only in case of **external** control.

**Possible parameters:** 0 ... 5

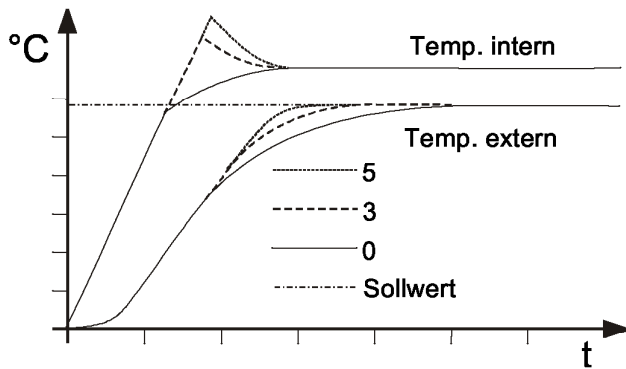

During selftuning, the control parameters  $X_p$ ,  $T_n$  and  $T_v$  of a controlled system are automatically determined and stored. Depending on the controlled system, time for tuning can be unequally longer. This controller layout allows protection of sensitive objects requiring temperature application.

As soon as a co-speed factor is set, it is considered for calculating the control parameters. As shown in the diagram, tuning times become shorter the higher the co-speed factor is, but overshooting can happen in the internal system.

**Control parameters – intern / extern**

The control parameters preset in factory are in most cases adequate for achieving an optimum temperature pattern for the samples requiring temperature application.

Each parameter may be manually set via the keypad if necessary, to allow optimum control performance.

Setting range:  
internal / external  
0.1 ... 99.9 K

**Proportional range >Xp<**

The proportional range is the range below the selected temperature value in which the control circuit reduces the heating power from 100 % to 0 %.

Setting range:  
internal / external  
3 ... 9999 s

**Resetting time >Tn< (Integral component)**

Compensation of the remaining control deviation due to proportional regulation. An insufficient resetting time may cause instabilities to occur. Excessive resetting time will unnecessarily prolong compensation of the control difference.

Setting range:  
internal / external  
0 ... 999 s

**Lead time >Tv< (Differential component)**


The differential component reduces the control settling time. An insufficient lead time will prolong the time required to compensate for disturbance effects and cause high overshooting during run-up. An excessive lead time could cause instabilities (oscillations) to occur.

Setting range:  
0.1 ... 99.9 K

**Proportional range >Xpu<**





The proportional range Xpu of the cascaded controller is only needed for external control.

**Optimization instructions for the PID control parameters:**

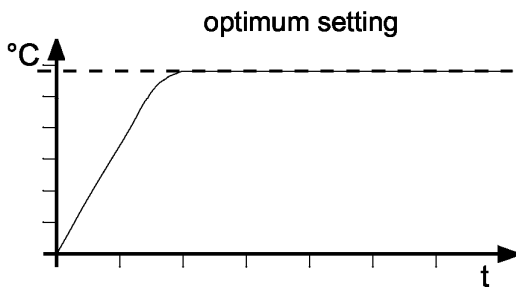


The heat-up curve reveals inappropriate control settings.

**Inappropriate settings may produce the following heat-up curves:**

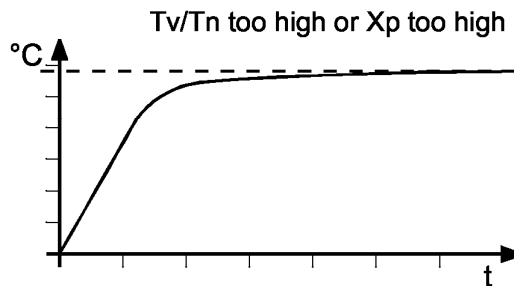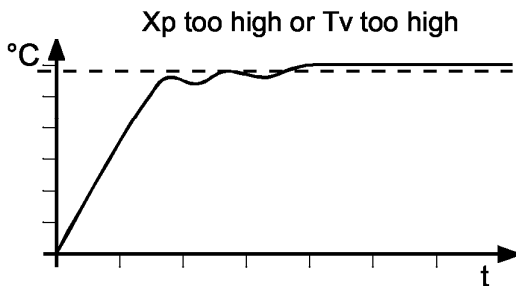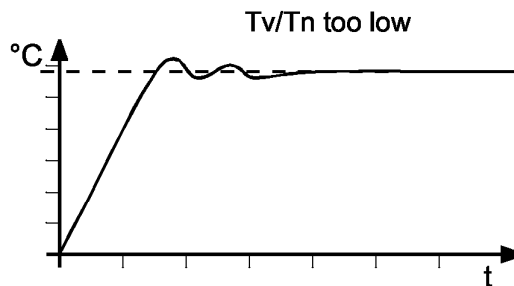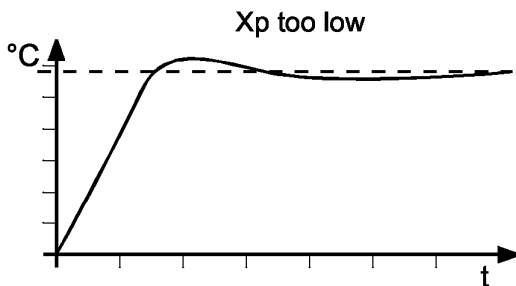

### 10.3. Start of a profile

The start menu of the integrated programmer allows calling up and defined starting of one of six previously stored temperature profiles. The profiles are started manually or via the integrated timer.

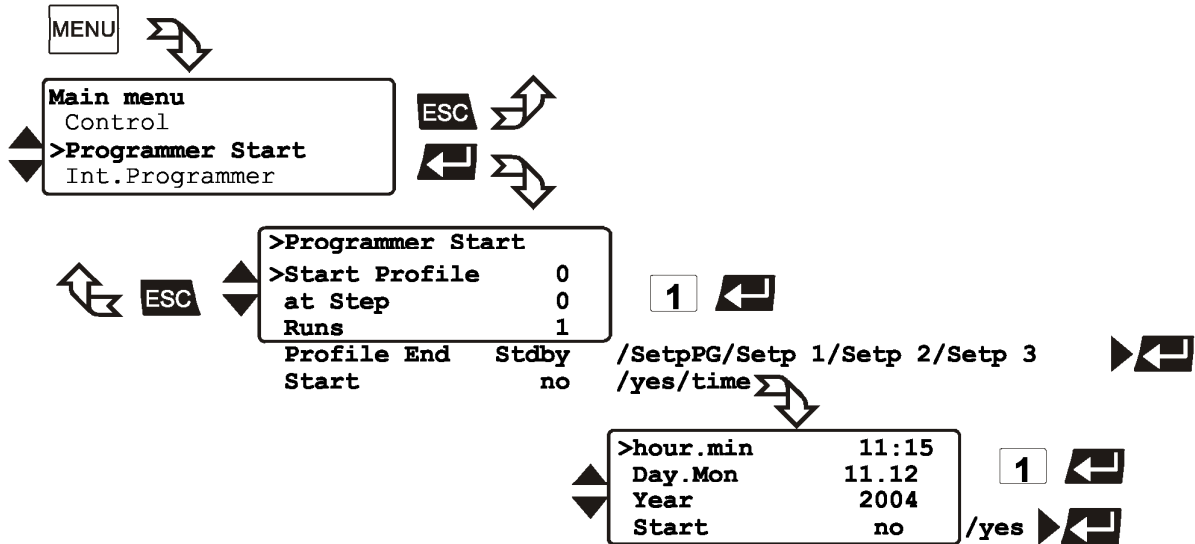

Example: 11:15



1. Use the keys to select the menu item.
  2. With the keys select the parameter (parameter blinks) or use the numeral keypad to enter the numbers (number segment blinks).
  3. Press enter to store the new parameter or value.
- ① If the parameter **time** is selected, a further submenu opens for setting the start time.

Possible parameters:

**Start Profile 1 to 5**

Select the profile that should be started.

**at Step 1 to 60.**

Select the section with which the profile should be started.

**Runs 1 to 99**

Select the number of profile repetitions.

**Start no / yes** ⇒ (manual start)

or

**time** ⇒ (via integrated timer)

**Profile End Stdby / SetpPG / Setp 1 / Setp 2 / Setp 3**

This parameter is used to determine if the unit switches to the -OFF- condition at the end of a profile, or if temperature control is continued and the working temperature setpoint that is be used in such case.

Examples:

```

-OFF-
S 25.00
E -----
    
```



```

Setpoint3: 80.00°C
IntAct    : 80.00°C
Power     : 2%
Control   : intern
    
```

```


Setpoint
Setpoint1: 25.00°C
>Setpoint2: 58.00°C
Setpoint3: 85.00°C
    
```

```


Setpoint2: 58.00°C
IntAct    : 60.00°C
Power     : 0%
Control   : intern
    
```

```

60.00
S 58.00
E -----
    
```




**Stdby** – If the end of the profile is reached, the Cryo-Compact Circulator switches to the -OFF- condition. The VFD display indicates the working temperature setpoint, that was displayed in line 2 already before the program was started (example: S 25.00 °C ).

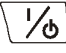**SetpPG** – (Setpoint of the ProGrammer) At the end of the profile, the setpoint of the last program section is written to the position of setpoint 3 (example: 80.00 °C) and indicated in line 1 of the LCD display. The Cryo-Compact Circulator operates and controls this temperature until the Start/Stop key  is pressed or a new profile is started.


**Setp 1, Setp 2, Setp 3** – („Setp“ short for **Setpoint**)

Before starting the programmer, a temperature value is set for one of

the three setpoints in the  menu.

(example: Setpoint 2 58.00 °C)

The Cryo-Compact Circulator operates and controls this

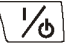temperature at the end of the profile until the Start/Stop key  is pressed or a new profile is started.

**Start no / yes** ⇒ (manual start)  
 or  
**time** ⇒ (via integrated timer)

When selecting the parameter **time**, a new submenu is called up for entry of the start time.



A flashing segment indicates that a start time needs to be entered.

**hour.min** Start time  
**Day.Mon** day and month  
**Year** year

Set each entry with enter .

**Start no / yes**

A flashing line indicates that the parameter „yes“ needs to be entered.

Press the key  to select the parameter and press enter .

```

Starttime: 06:00
Date      : 15:12
*** wait ***
Acttime   02:34:45
    
```

The temperature system switches to waiting mode and a flashing line „wait“ appears on the LCD DIALOG-DISPLAY. The start time and actual time are permanently indicated on the display.

**Indication after successful start:**

The started programmer displays the actually calculated setpoint in line 1. The value changes until the final temperature of the section is reached (temperature ramp).

Examples:

```
Setp. :    125.8°C
IntAct:    124.7°C
Prof. :    4 Step: 12
Remain:    00:03:45
```

```
Setp. :    125.8°C
IntAct:    124.7°C
Prof. :    4 Step: 12
All :    00:03:45
```

```
Setpoint :    125.8°C
IntAct :    124.7°C
Prof :    4 Step:12
S12:150:00 h:m 01:10
```

```
Edit
>Edit Profile:    4
Step:            12
Setp:            150.00°C
Time [hh:mm]    01:10
```

1<sup>st</sup> line: Setpoint of the programmer

2<sup>nd</sup> line: Actual temperature value  
 at **internal control** = IntAct: xxx.xx  
 at **external control** = ExtAct: xxx.xx

3<sup>rd</sup> line: Selected profile and the actual section (step).

4<sup>th</sup> line: With the keys ▼ ▲ the following information can be displayed in line 4.

- Remaining time of the actual section (standard display)
- Remaining total time: Profile x number of repetitions  
**(All: Days/hours:minutes:seconds)**

**Particularitys:**

1. If a total time of more than 999 days is calculated, only the hours are displayed on the unit.  
**(All: hours:minutes:seconds)**
2. If the time period of a section is edited after start, then the remaining total time can no longer be displayed.

- The data of a section as stored under „Edit Profile“.

Use the keys ◀ ▶ to successively indicate the sections of a profile. Unused sections or incorrect data are skipped.

**10.3.1. Edit after start**

A JULABO programmer of generation II allows editing of a running profile. A programmer of the previous generation requested the profile to be interrupted, edited and then started again.

**Start and quit the edit mode:**






```
Edit
>Edit Profile:    1
Step:            6
Setp:            100.00°C
Time [hh:mm]    01:05
```

- After start, press the MENU key to enter the submenu „Edit Profile“.

Press **ESC** to quit the edit mode.



**Deleting a complete section:**

- Use the keys   to move to line 2.
- Use the keys   to choose the desired section.
- Press the  key to delete the complete section.

```

Edit
>Edit Profile:      1
**Step is deleted **
  Setp:             °C
  Time [hh:mm]
    
```

If the desired section is already chosen, the section is deleted even if the cursor > is in line 1, 3 or 4.

Deletion is indicated in the respective line of the LCD display:  
 \*\* Step is deleted \*\*




```

Edit
>Edit Profile:      1
** Step is active **
  Setp:             100.00°C
  Time [hh:mm]    01:05
    
```

The section that is currently processed cannot be deleted. If you still try it, the message **\*\* Step is active \*\*** appears on the LCD display.

**Changing a section:**










① The values in lines 3 and 4 can be changed, even on the section that is currently processed.

- In the input window, select the menu item with the keys  .
- Use the numeral keypad to enter the numbers and press enter .

```

Edit
  Step:              6
>Setp:              83.00°C
  time [hh:mm]     01:25
    
```

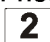Examples:

**Setpoint:** target temperature 83 °C     
**Time** time period 1 hour:25 minutes  
     


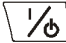
**Particularity for changes in a section that is currently in process.**


- ① If the target temperature is changed, the slope of the ramp is automatically calculated on basis of the remaining time of the section.
- ① If the time period is changed, the slope of the ramp is automatically calculated on basis of the new total time of the section.


### 10.3.2. Interrupting a profile

```

Setp. :    125.80°C
IntAct:    124.70°C
Prof. :    4 Step:12
*** pause ***
    
```

-  Press the start/stop key  to interrupt or continue a profile. The setpoint and time period set for the corresponding section are thus stopped at the values presently achieved. The instrument is put on hold and the message „pause“ flashes on the LCD DIALOG DISPLAY.

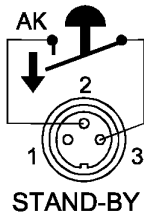


-  A profile can be interrupted or restarted by an external emergency shut-off.
 

**Important:**  
For this, in the menu >Inputs/Outputs< the menu item >Ext. StBy < must be set to >activ< (see page 52) and in the menu >Configuration< >Autostart< needs to be set (see page 31).

```

Setp. :    xxx.xx°C
IntAct:    xxx.xx°C
Prof. :    4 Step:12
*** STAND-BY ***
    
```

-  Setpoint control and timer are interrupted by breaking the contact „AK“. The instrument is put on hold and the message „Stand-By“ flashes on the LCD DIALOG-DISPLAY.

 **Notice: This is not an actual emergency switch-off.**




**Warning:**

Following a power interruption, it would be possible in this condition for the instrument to restart automatically. The safety and warning functions of the instrument should always be used to their fullest capacity. See **Warning** page 31.

### 10.3.3. Interruption after a power failure


There is every possibility to operate the programmer safe to power failures.

**Important:**  
For this in the menu >Configuration< >Autostart< needs to be set (see page 31).

-  If the AUTOSTART function is activated, the programmer starts again at a point approx. 20 seconds before the interruption took place. However, an uncontrolled change of the bath temperature happened.

```

Starttime: 06:00
Date :    15:12
*** wait ***
Acttime 02:34:45
    
```

-  Power failure when the instrument is on hold if start is effected via the built-in timer:  
If the starting time is not yet reached, the programmer is put on hold again.  
If the starting time is exceeded, the programmer starts immediately.

### 10.3.4. Termination of a profile

```

Setp : 125.80°C
IntAct: 124.70°C
Prof. : 4 Step:12
Remain: 00:03:45
    
```



```

Programmer Start
>Start Profile      x
at Step            x
Runs               1
Start              no
    
```

① A profile can be terminated by pressing **ESC**. The programmer switches back to the Start menu.

- Press **ESC** again to quit the menu or use the keys **▼▲** to remain in the Start menu.

The start of another temperature profile can now be prepared if necessary.

Example:

```

Setpoint1: 150.00°C
IntAct : 83.24°C
Power : 0 %
Control : intern
    
```

```

-OFF-
S 150.00
E -----
    
```

#### Interruption after a power failure


In case of a power failure, the reaction of the Cryo-Compact Circulator is the same as when switched off and on again with the mains switch.

The LCD DIALOG-DISPLAY gives information on the adjustments and conditions of the Cryo-Compact Circulator .

The VFD COMFORT-DISPLAY indicates "OFF".

## 10.4. Integrated programmer

The integrated programmer allows any desired temperature program sequences to be realized. Such a temperature sequence is called profile. A profile consists of individual sections defined by duration (t:) and target temperature. Target temperature is the setpoint (T:), that is achieved at the end of a section. The programmer uses time and temperature difference values within a section to calculate a temperature ramp.

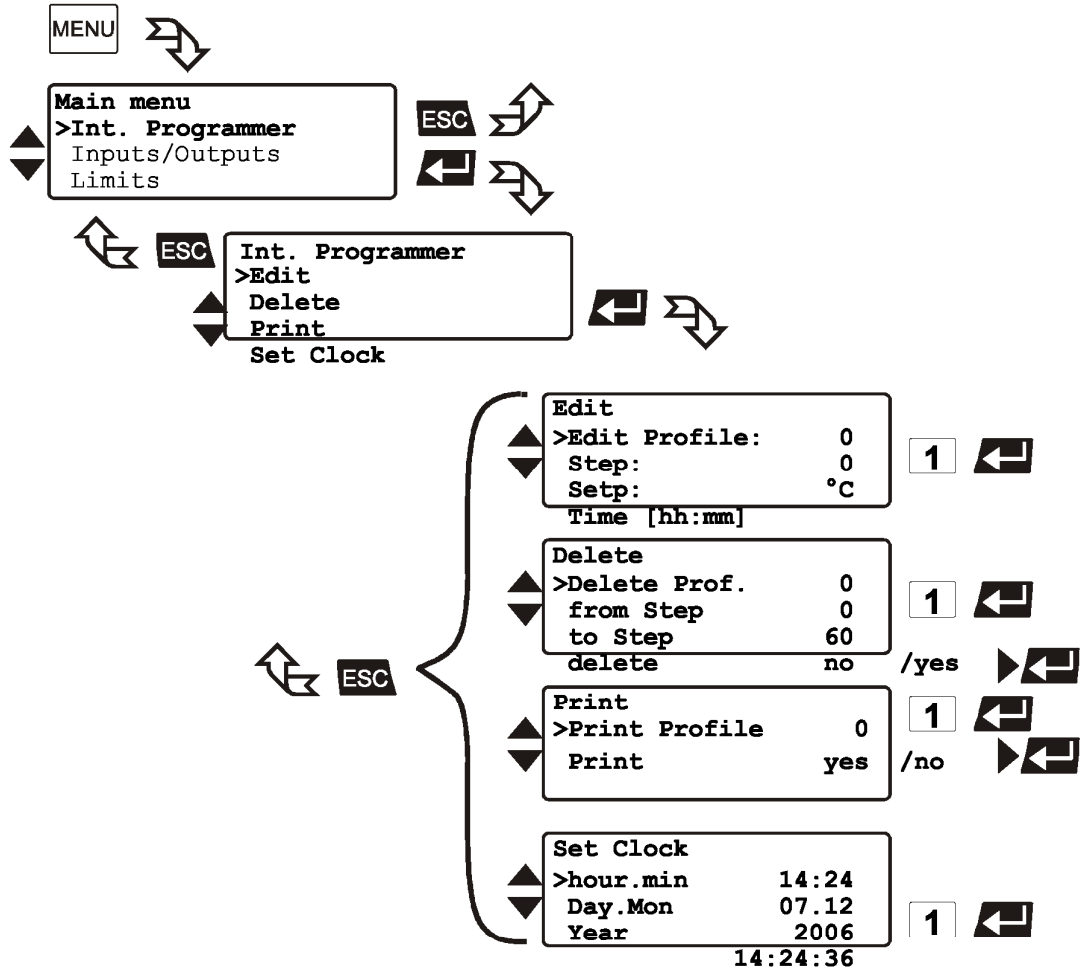

1. Use the keys ▼ ▲ to select the desired menu item in menu level 2 on the LCD DIALOG-DISPLAY (➤) and press enter to change to menu level 3. An input window opens for each menu item.
2. In the input window, select the menu item with the keys ▼ ▲.
3. Use the keys ◀ ▶ to select the parameter (parameter blinks) or use the numeral keypad to enter the numbers (number segment blinks).
4. Press enter to store the new parameter or value.

```

Int. Programmer
>Edit
Delete
Print
Set Clock
    
```

- Edit**            Compile profiles  
                      Display sections
- Important:**    Also a running profile can be changed.  
                      See „Edit after start“ page 40
- Delete**           Delete sections
- Print**             Print a programmed profile
- Set clock**        Set the real time on the instrument

**Compile profiles:**

A flashing segment indicates that a number needs to be entered. Under submenu „Edit Profile“ enter a profile number. Six profiles may be stored (nos. 0 to 5).

Example:

```

Edit
>Edit Profile:    1
Step:             6
Setp:            100.00°C
Time [hh:mm]    01:05
    
```

**Edit Profile:** Profile No.1            **1**

Then programme the desired values for each section. Use the keypad to set section number, target temperature and time period. Set each entry with enter .

**Step:**            Section No. 6                            **6**

**Setpoint:**      target temperature - 100 °C        **1 0 0**

**Time**             time period - 1 h:05 minutes        **0 1 - 0 5**


When the program is running, only sections having complete information for target temperature and time period are considered. It makes sense, to leave out section numbers in the profile, in order to use them later for corrections in the profile

**Important:**

If a time of 00:00 is set for a profile, the profile is continued with the next section only after the setpoint temperature ( $\pm 0.2$  °C) is reached.

Example:

Step	(No.)	1	3	5	6	9	11	14
Setpoint	(°C)	50	50	100	100	80	80	20
Time	(h.m)	00:08	00:02	00:10	01:05	00:10	00:05	00:30

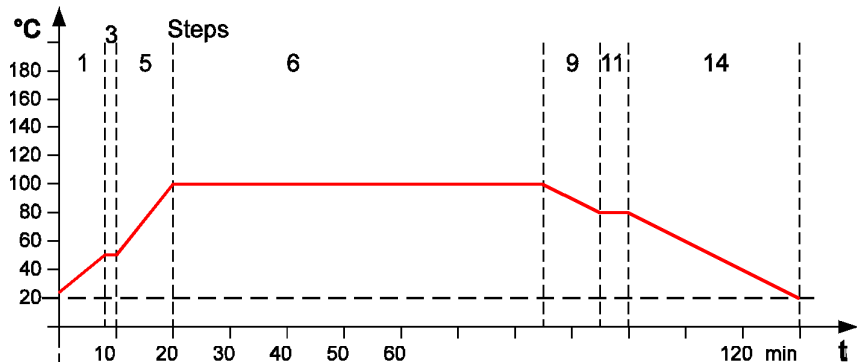

**Display sections (Step):**

```

Edit
>Edit Profile:    1
Step:            3
Setp:           50.00°C
Time [hh:mm]    00:02
    
```

1 Use the keys to select the submenu „Step“, enter the desired number **3** and press enter

or

2 If the submenu „Step“ is selected, section by section can be shown in ascending order by the key .

**i** Each previously set value is displayed.

Example:

```

Delete
>Delete Prof.    4
from Step       8
to Step        34
delete
                yes
    
```



```

Delete
>Delete Prof.    4
from Step       8
***** delete *****
    
```

**Delete:**

A flashing segment indicates that the respective profile number needs to be entered in which one or more consecutive sections are to be deleted.

Example:

Delete step 8 to step 34 in profile 4.

```

4  8  3 4   yes 
    
```

Line 4 indicates the deletion.

Example:

```

Print
>Print Profile  2
Print          yes
    
```



```

Print
>Print Profile  2
***** print *****
    
```

**Print:**

Each profile may be printed via the serial interface for control or documentation.

A flashing segment indicates that the number of the profile to be printed needs to be entered.

Example: Print profile 2

```

2  yes 
    
```

Printing is indicated in line 2.

**Setting clock**

The integrated clock allows starting a profile at any date and time.

The clock is preset in the factory.

**i** Lines 1 to 3:


Check for correctness of the preset date and time and correct if necessary.

The time is displayed permanently in line 4.

```

Set Clock
>hour.min      14:24
Day.Mon        07.12
Year           2006
                14:24:36
    
```

## 10.5. Analog inputs/outputs

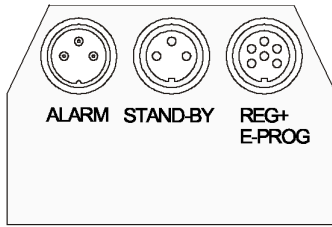


**i** In order to use the analog inputs and outputs, the Cryo-Compact Circulator must be equipped with the electronic module available as option.

Order No. 8 900 100 Electronic module

This submenu enables setting of the input and output values for the programmer input and the temperature recorder outputs of socket REG+E-PROG.





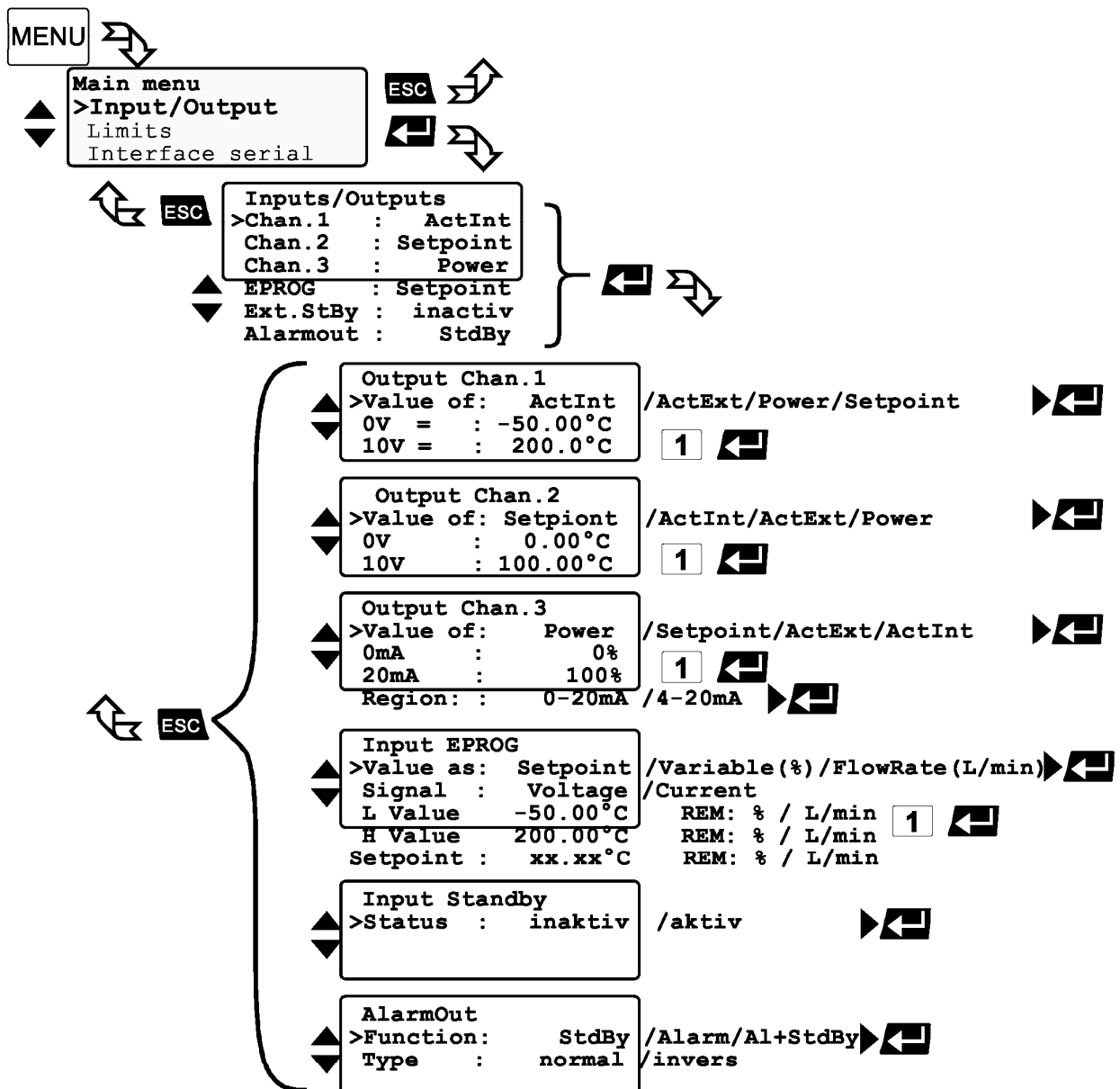





The >STAND-BY< input and the >ALARM< output are configurable.

**i** If the electronic module is not fitted, the item >inputs/outputs< will be missing in the menu.



Menu level 2

Inputs/Outputs	
>Chan.1	: ActInt
Chan.2	: ActInt
Chan.3	: Setpoint
EPROG	: Setpoint
Ext.StBy	: inactiv /activ
AlarmOut	: StdBy

1. Use the keys   to select the desired (➤) menu item in menu level 2 on the LCD DIALOG-DISPLAY.
2. Use the keys  and  to store the desired parameter or press enter  to enter menu level 3. An input window opens for each menu item.
3. In the input window in menu level 3 select the menu item with the keys  .
4. Use the keys   to select the parameter (parameter blinks) or use the numeral keypad to enter the numbers (number segment blinks).
5. Press enter  to store the new parameter or value.


Menu level 2

<b>Channel 1</b>	voltage output for recorder (V)
<b>Channel 2</b>	voltage output for recorder (V)
<b>Channel 3</b>	current output for recorder (mA)
<b>EPROG</b>	external programmer input
<b>External standby</b>	Standby input (for external emergency switch-off)
<b>Alarm output</b>	Output for external alarm signal



10.5.1. REG+E-PROG – temperature recorder outputs

REG+E-PROG (11): Outputs of the connector

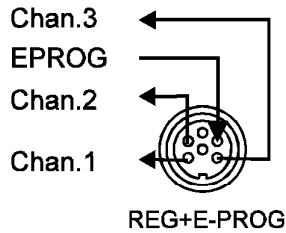

1. First define the desired output value for channels 1 to 3:

- ActInt** internal actual temperature value (bath temperature)
- ActExt** ernal actual temperature value (external sensor)
- Power** periodic or intermittent heating or cooling
- Setpoint** active setpoint temperature  
(Setpoint1, 2, 3,/ integ.r. programmer /external programmer)

2. Then select the display size for channels 1 to 3:

**Voltage outputs channels 1 and 2**


Assign the voltage values of 0 V to the lowest and 10 V to the highest necessary temperature or power rating required as an output value (°C / %).

**Current output channel 3**

Assign the current values 0 mA or 4 mA to the lowest and 20 mA to the highest temperature or power rating required as an output value (°C / %).

3. The current output (chanell 3) offers 2 ranges for selection:  
0 to 20 mA or 4 to 20 mA

**Examples:**

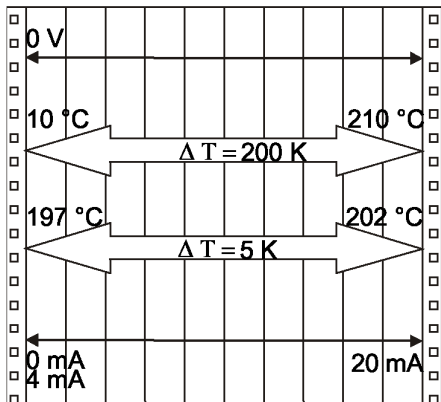

Chan. 1	
>Value of :	ActInt
0V	= 10°C
10V	= 210°C

lowest temperature value: 10 °C  
 highest temperature value 210 °C  
 Fig. shows 200 °C scaled to paper width  
 rise: 50 mV/°C

lowest temperature value: 197 °C  
 highest temperature value: 202 °C  
 Fig. shows 5 °C scaled to paper width  
 rise: 2000 mV/°C

### 10.5.2. REG+E-PROG - programmer input

#### Menu >Configuration<

```

Configuration
>Setpoint :      eprog      Keypad/RS232/extPt100
Autostart:      off
Standby  :      pump off
Language  :      english
Reset    :      no
Variable  :      Control /serial/eprog
    
```

#### Menu >Inputs/Outputs<

```

Input EPROG
>Value as: Flowrate      Setpoint/Variable
Signal  : Voltage
L Value : 0.0L           (Liters/Minute)
H Value : 50.0L
Flowrate: xx.xL
    
```

#### E-PROG - Input

Setting needs to be carried out, if

1. setpoint programming is to be made via an external programmer.  
For this, in the menu >Configuration< first set the menu item >Setpoint< to >eprog<.
2. the heater variable should be controlled via an external control pulse.  
For this, in the menu >Configuration< set the menu item > Variable.< to >eprog<.
3. the signal of an external flow meter should be registered.

- ⓘ The E-Prog input can only be used either under menu item >Setpoint< or under menu item > Variable <. If the input is neither occupied by >Setpoint< or > Variable <, the signal of a flow meter can be connected.

- ⓘ First set menu >Configuration<, then menu >Inputs/Outputs<.

- Connect the external programmer to socket (11) REG+E-PROG of the Cryo-Compact Circulator .


#### Selecting the signal:

The programmer (E-PROG) input of the Cryo-Compact Circulator can be matched to the output signal of the external programmer.

**Voltage**            voltage input  
**Current**            current input


#### „L Value“ - Setting the LOW value: (See below ↻)

Adjust and set the lowest desired working temperature on the programmer (e.g. 0 °C).

Enter the same value on the Cryo-Compact Circulator by pressing the appropriate buttons on the keypad and press enter  to set.

#### „H Value“ - Setting the HIGH value: (See below ↻)

Adjust and set the highest desired working temperature on the programmer (e.g. 300 °C).

Enter the same value on the Cryo-Compact Circulator by pressing the appropriate buttons on the keypad and press enter  to set.

#### Example:

```

Input EPROG
>Value as: Setpoint
Signal  : Current
L Value : 0.0°C
H Value : 300.0°C
Setpoint: 50.0°C
    
```


#### Example:

- Set a temperature of 50 °C on the external programmer!  
The value adjusted and set on the external programmer is displayed in line 4 of the LCD DIALOG-DISPLAY for control purposes.

```
Eprog : 50.0°C
IntAct : 24.7°C
Power : 80 %
Control : intern
```

❗ After returning the LCD display to standard display by pressing **ESC** this value is displayed in line 1 (Example: Eprog : 50.00 °C).

This EPROG input enables the use of different voltage and current values as program parameters.

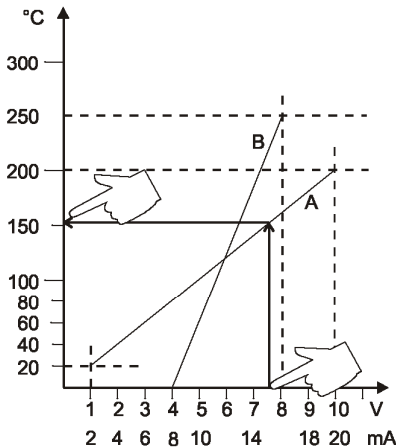

- **L Value“ - Setting the LOW value:** (See below ↻)
  - 1) Adjust and set the lowest desired value on the voltage or current source resp. (Example A: 1 V).
  - 2) Assign a lower temperature threshold value to this adjusted voltage/current value by pressing the appropriate buttons on the keypad of the instrument (Example A: 20 °C ) and set by pressing enter ↵.
- **„H Value“ - Setting the HIGH value:** (See below ↻)
  - 1) Adjust and set the highest desired value on the voltage or current source resp. (Example A: 10 V).
  - 2) Assign an upper temperature threshold value to this adjusted voltage/current value by pressing the appropriate buttons on the keypad of the instrument (Example A: 200 °C) and set by pressing enter ↵.
- Return to the standard display by pressing **ESC**.

❗ Example B in the diagram serves to illustrate that the end point values are freely selectable.

```
Input EPROG
>Value as: Setpoint
Signal : Voltage
L Value : 20.0°C
H Value : 200.0°C
Setpoint: 152.0°C
```


**Example out of diagram A:**


- Adjusting the voltage source for an output of 7.6 V!
- Line 5 of the LCD DIALOG-DISPLAY shows the externally set setpoint value. The instrument calculates this value from the rise angle of the two pre-decided end points (in example A: 7.6 V correspond to an external setpoint temperature of 152.0 °C ).



```
Eprog : 152.00°C
IntAct : 50.76°C
Power : 100 %
Control : intern
```

After returning the LCD display to standard display by pressing **ESC**, this value is displayed in line 1 (Example: EPROG 152.00 °C).

 **Notice:**  
If this adjustment is not correctly performed at two different points, the setpoint setting will be incorrect.

 **Important:**  
The usable temperature range between > L Value < and > H Value < is limited to the configured working temperature range of the Cryo-Compact Circulator resp. of the unit combination (working temperature range see technical specifications on page 8)


### 10.5.3. STAND-BY input

#### Ext. StBy (10)

(for external emergency switch-off)  
(Connector see page 65)

#### Possible parameters:

- inactiv** - standby input is ignored
- activ** - standby input is active

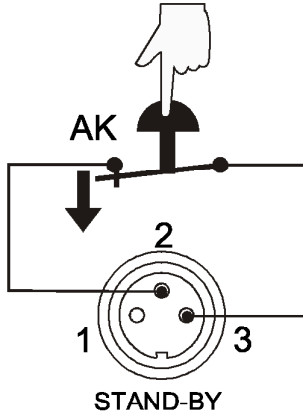


#### Activate the standby input:

1. Under menu item **>Ext. StBy<**, set the parameter to **>activ<**.
2. Connect an external contact 'AK' (e.g. for emergency switch-off) or an alarm contact of the superordinated system.

In case the connection between Pin 2 and Pin 3 is interrupted by opening the contact 'AK', a complete shutdown of the circulating pump and heater is effected, and the unit enters the condition „E OFF“.

As long as the contact remains open, line 4 of the LCD DIALOG-DISPLAY flashes and displays the message „STAND-BY“.

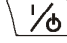
If the contact is reclosed, the instrument returns to the standby state and „E OFF „ is displayed. Press  to start.


```
Setpoint1: 150.00°C
IntAct   : xxx.xx°C
Power    :   xx %
STAND-BY
```

```
E OFF
S xx.xx
E -----
```

#### **ⓘ Additional tips for using the STANDBY input:**

The standby function can be used in conjunction with the AUTOSTART feature (see page 31).


1. If the Autostart function is NOT turned ON, the standby input is used as described above.
  2. If the Autostart function is enabled, the instrument will revert back to the original method of entering the setpoint (i.e. keypad, RS232, Analog input, etc.).
- Entering the setpoint with the keypad. e.g.   
As described above, a bipolar shut-down is accompanied by the „STAND-BY“ display and the OFF state. The programmable controller starts again when the contact is reclosed. The temperature of the bath fluid changed during the STAND-BY state.
  - Entering the setpoint with the programmer (see pages 42). The display „STAND-BY“ appears. The setpoint value and the time are both held at the current value. The temperature of the bath fluid will be held constant at this temperature. The programmer continues once the contact is reclosed.

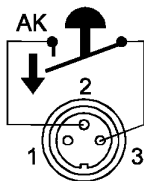

```
Setpoint : xxx.xx°C
IntAct   : xxx.xx°C
Profil   : 4 Step:12
*** STAND-BY ***
```

 **Notice: this is not an actual shutoff feature.**

10.5.4. ALARM output

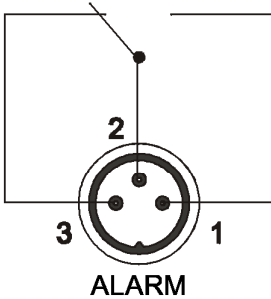

**Alarm output (9)**

(for external alarm signal)

This socket is a potential-free change-over contact. With the adjustments in the menu >Inputs/Outputs< all operating conditions can be signaled without having to change the pin assignments.

Menu >Inputs/Outputs<

```

AlarmOut
>Function:      StdBy /Alarm/Al+StBy
Type           :      normal /invers
    
```

Signification of the terms under menu item >Function< :

The Cryo-Compact Circulator is in condition

>StandBy< or >Alarm<

```

- OFF -
S xx.xx
E -----
    
```

```

XXXXXX
ALARM
CODE xx
    
```


For >Type: normal< pins 2 and 3 are connected in any case according to the selected >Function<.

For >Type: invers< pins 2 and 1 are connected in any case according to the selected >Function<.

- Switching capacity    max. 30 W / 40 VA
- Switching voltage    max. 125 V~/–
- Switching current    max. 1 A

## 10.6. Limits

The menu >Limits< allows for setting the minimum and maximum values for all important setting ranges and capacity variables.

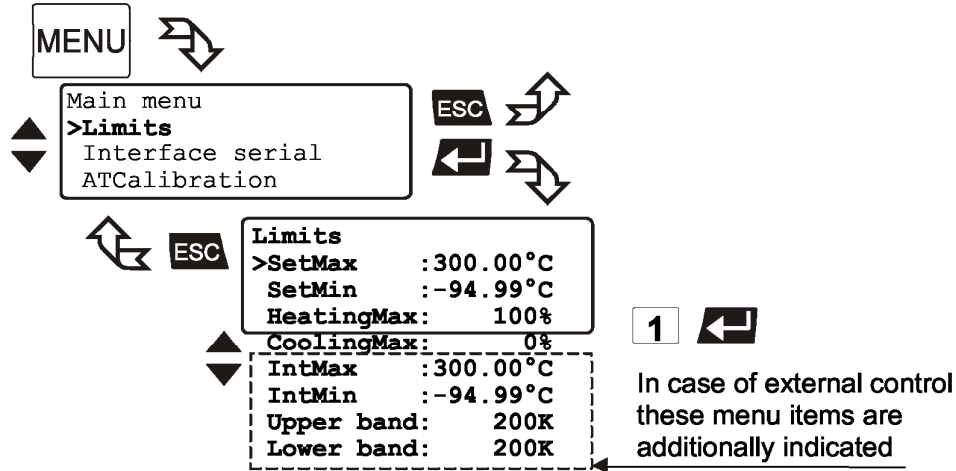

**SetMax:** and **SetMin:** Maximum and minimum setpoint

Restriction for the adjustable temperature range.

Setting range: see working temperature range

The limitation of the operating temperature range effects the temperature

setting under the menu called up with the **T** key - Setpoint 1, 2, 3.

It is possible to adjust only working temperatures that lie within the limit range set here.

Existing settings for Setpoint 1, 2, 3 and also for >OverTemp< and >SubTemp< (see page 25) are automatically defered within the limit range.

**HeatingMax:** and **CoolingMax:**

Adjusted maximum heating / cooling.

Setting range: 0 to 100 % in steps of 1 %

Heating and cooling powers of the Cryo-Compact Circulator are adjustable. 100 % corresponds to the values in the technical specifications of the equipment.

Factory setting: **CoolingMax**

CF31 100 % (fixed value)

CF41 0 to 100 % in steps of 1 %

**IntMax:** and **IntMin:** Limit setting  
 Maximaler und minimaler Sollwert im internen Bad.  
Setting range: see working temperature range 8


The limits IntMax and IntMin are only valid for external control. IntMax and IntMin are used to limit the expected internal bath temperatures to any upper and lower values. The temperature controller cannot exceed these limits even if it would be required for reaching the temperature in the external system. Consequently the external setpoint may thus not be reached.



Sense of a limit setting:

- Protects the bath fluid from overheating.
- Prevents an undesired alarm shut-down by the excess temperature protection - >ALARM CODE 14<.  
 Set >IntMax:< to a value at least 5 °C below the >SafeTemp:< value.
- Protects the pump motor from high viscosity of the bath fluid at low temperatures.
- For refrigerated Cryo-Compact Circulator s: Freezing protection when using water as bath fluid.

**Upper band:** and **Lower band:** Band limit  
 The band limit is active for external control.

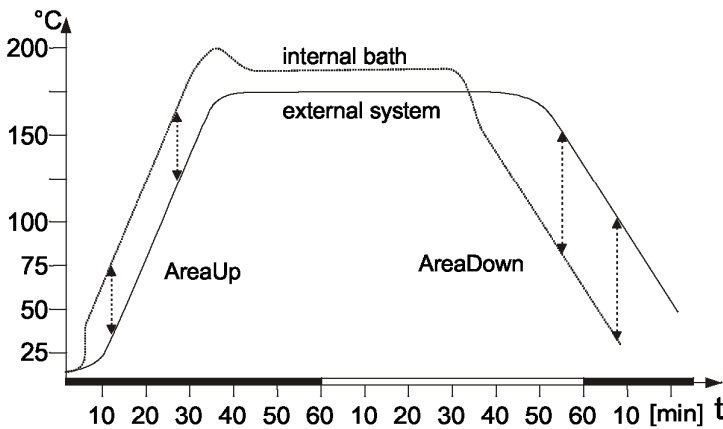

For the heat-up and cool-down phases different settings to conform to the requirements of the particular application are possible.

Setting range: 0 ... 200 °C

Using Upper band and Lower band, the difference between the temperatures in the internal bath and the external system can be limited to any maximum value for the heat-up or the cool-down phase.

During the heat-up phase the difference value always adds to the actual external temperature. During the cool-down phase, the difference value is subtracted.


Sense of a band limit:

- Ensures gentle heating for the objects requiring temperature application.
- Protects glass reactors, for example, from thermal shock.

ⓘ The **IntMax:** and **IntMin:** values are superordinated to this band limit.

## 10.7. Serial interface

For communication between Cryo-Compact Circulator and a PC or a superordinated process system the interface parameters of both units must be identical. The adjustment is usually carried out one time only in the menu >Interface serial<.

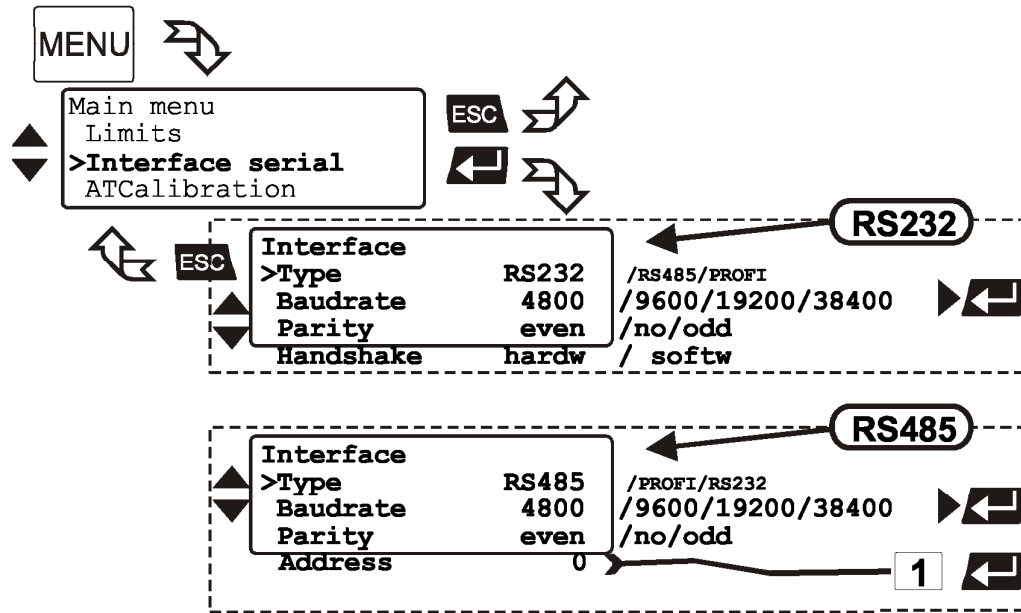

<b>Type</b>	RS232 / RS485
<b>Baud rate</b>	4800/9600/19200/38400
<b>Parity</b>	none/even/odd
<b>Handshake</b>	software handshake/hardware handshake


**Address** 0 to 127

Factory settings:  
 4800 Bauds  
 even  
 hardware handshake



## 10.8. ATC Absolute Temperature Calibration, 3-point calibration

ATC serves to compensate a temperature difference that might occur between Cryo-Compact Circulator and a defined measuring point in the bath tank because of physical properties.

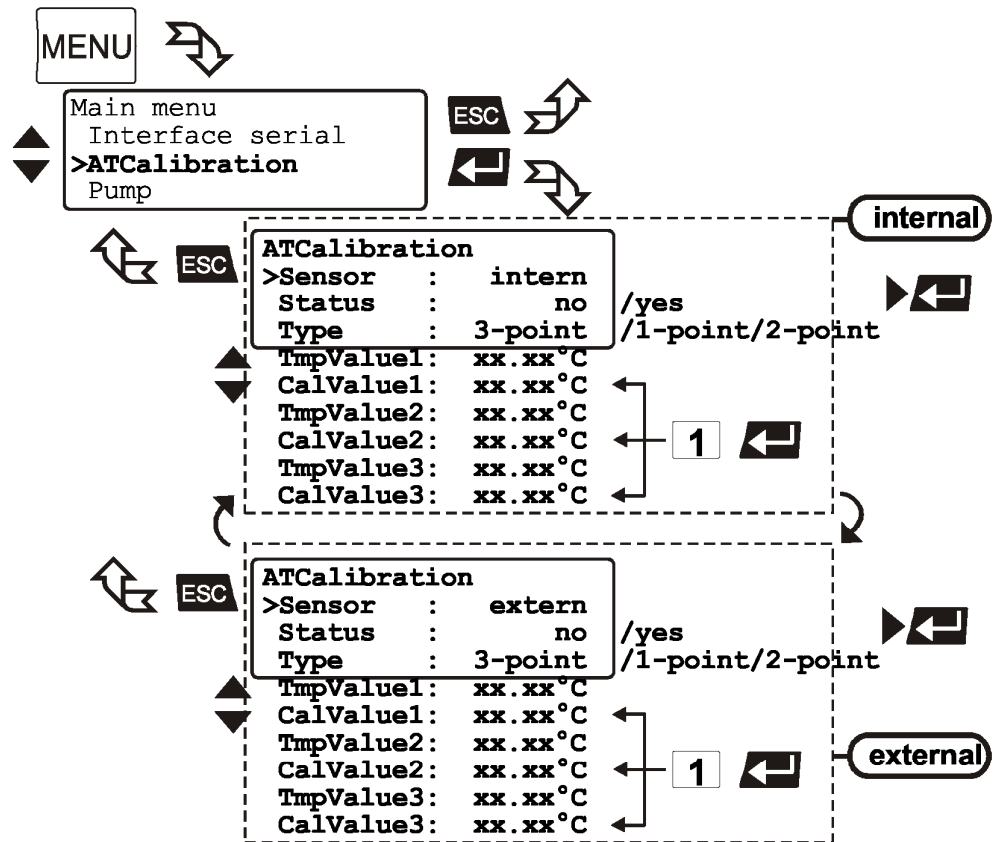

### Sensor: >intern<, >extern<

Calibration can be carried out for the internal temperature sensor and for the external temperature sensor connected to the socket „ext. Pt100“.

The Cryo-Compact Circulator is able to store both parameter sets. However, only the one set under this menu item is indicated.

### Status:

**>no<** The controller of the Cryo-Compact Circulator uses the original curve of the temperature sensor.  
Important: During the calibration process **>no<** needs to be set.

**>yes<** The controller of the Cryo-Compact Circulator uses the new calibration curve.

**Type:**

A >1-point<, >2-point < or >3-point < calibration can be carried out.

First geometrically define the location for calibration (measuring point  $T_M$ ), then determine the temperature values of the calibration points.

The calibration type also determines the number of pairs of values indicated on the LCD DIALOG-DISPLAY.

**TmpValue:** 1 or 2 or 3


Defined temperature value of the calibration point. This value is automatically stored with >CalValue< and can be indicated for control purposes.

**CalValue:** 1 or 2 or 3


The „Calibration value“ is determined with a temperature measuring device and stored under menu item >CalValue<.

**Example:**

**1-point calibration**

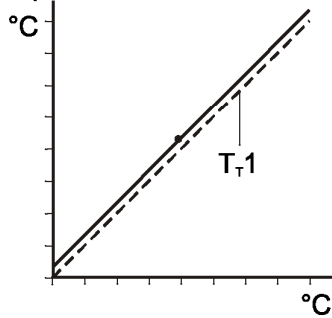


**3-point calibration**

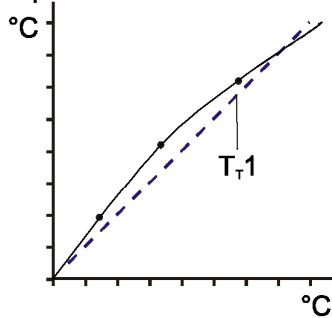

$T_T 1$  = Original curve

**Principle:**

For ATC calibration, in steady state the bath temperature at the location of the temperature sensor ( $T_M$ ) is determined at the respective adjusted working temperature. This value is then set on the Cryo-Compact Circulator in the menu >ATCalibration< under menu item >Calwert X<. This can be a 1-point, 2-point or 3-point calibration.

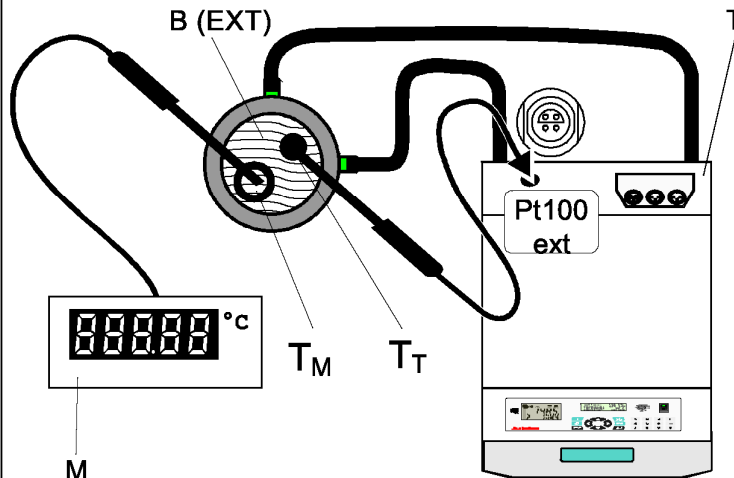

M = Temperature measuring instrument with temperature sensor

B = Bath tank (internal or external)

T = Cryo-Compact Circulator

$T_M$  = Temperature on measuring point

$T_T$  = Temperature on Cryo-Compact Circulator

1. Use the keys to select the menu item.
2. With the keys select the parameter (parameter blinks) or use the numeral keypad to enter the numbers (number segment blinks).
3. Press enter to store the new parameter or value.

## Examples:

$T_T = 80.00\text{ }^{\circ}\text{C}$   
 $T_M = 79.73\text{ }^{\circ}\text{C}$   
 $T_T = 120.00\text{ }^{\circ}\text{C}$   
 $T_M = 119.51\text{ }^{\circ}\text{C}$   
 $T_T = 160.00\text{ }^{\circ}\text{C}$   
 $T_M = 159.34\text{ }^{\circ}\text{C}$

```
Control
>ControlType: intern
Selftuning:  once
Dynamic :  aperio
```




```
Setpoints
>Setpoint1:  80.00°C
Setpoint2:  37.00°C
Setpoint3:  70.00°C
```

```
ATCalibration
>Sensor :  intern
Status  :  no
Type    :  3.point
TmpValue1: 80.00°C
CalValue1: 79.73°C
TmpValue2: xx.xx°C
CalValue2: xx.xx°C
TmpValue3: xx.xx°C
CalValue3: xx.xx°C
```




## Example:

3-point calibration for internal control.

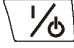In the temperature range of 80 °C to 160 °C the calibration curve of the temperature sensor ( $T_T$ ) should be assimilated to the actual temperatures on the measuring point ( $T_M$ ).

- Press the  button. Under the menu >Control< set >ControlType< to >intern< and press enter . (**ControlType**: Can only be adjusted in the **-OFF-** condition)
- Press start/stop .

**Calibration procedure:**

- Press  and set the first temperature value under >Setpoint1< for example (example 1<sup>st</sup> value = 80 °C).
- Wait until this temperature is maintained constant in the bath for about 5 minutes.
- Press the  button and open the menu >ATCalibration<.
- Set the following:  
Menu item >Sensor< to >intern< ,  
Menu item >Status< to >no< ,  
Menu item >Type< to >3-point<.
- Read the value of  $T_M$  on the temperature measuring device and enter the value under menu item >CalValue 1< (79.73 °C) using the numeral keypad.  
Press enter  and the Cryo-Compact Circulator also stores the value of  $T_T$  as value for >TmpValue 1< (80.00 °C) .  
The first of the 3 points is now calibrated.
- Repeat the calibration procedure for 120 °C and 160 °C.


## 10.9. Setting the pump pressure

Examples:



The pressure of the circulating pump is adjustable in four grades. After setting, the VFD-Info-Display indicates the corresponding value.


Adjustable pump capacity stage 1 ... 4

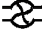




Illuminated display:  for pump pressure

- ① Display for the adjusted pump pressure stage in the –OFF- mode.
- ① Display for the effective pump pressure stage (rotation speed) after start.

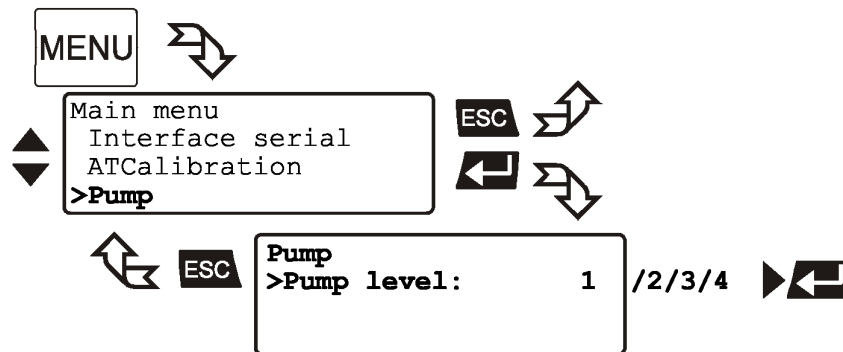For protecting the pump motor, the rotation speed changes with the load applied, for example depending on the viscosity of the bath fluid at different working temperatures.

Flow rate:	22 ... 26 l/min			
Pump capacity stage	1	2	3	4
Pump pressure:	[bar] 0,4	0,5	0,6	0,7
Suction pump	[bar] 0,2	0,26	0,33	0,4
Total capacity in a loop circuit	[bar] 0,6	0,76	0,93	1,1



1. Use the keys   to select the menu item.
2. With the keys   select the parameter (parameter blinks).
3. Press enter  to store the new parameter or value.

## 11. Troubleshooting guide / Error messages

**XXXXXX**  
ALARM  
CODE 01



**XXXXXX**  
WARNING  
CODE 40

**Alarm** with a complete shutdown of the unit

Whenever the microprocessor electronics registers a failure, a complete shutdown of the compressor, heater and circulating pump is performed.

The alarm light „“ illuminates and a continuous signal tone sounds.

The VFD-COMFORT-DISPLAY indicates the cause for the alarm in form of a code.

**Warning** without a complete shutdown of the unit

The VFD-COMFORT-DISPLAY indicates the cause for the warning in form of a code and an acoustic signal sounds in regular intervals.

These messages appear every 10 seconds.



Press enter to quit the audible signal.

Press **F** to indicate the help text in the LCD DIALOG-DISPLAY when alarm messages appear.

### ALARM

CODE 01 **F**

Insufficient liquid level or float is defective.

- The Cryo-Compact Circulator is operated without bath fluid, or the liquid level is insufficient. Replenish the bath tank with the bath fluid.
- Tube breakage has occurred (insufficient filling level due to excessive bath fluid pumped out). Replace the tubing and replenish the bath tank with the bath fluid.
- The float is defect (e. g., because damaged in transit). Repair by authorized JULABO service personnel.
- 

### WARNING

CODE 03 **F**

Actual temperature above high temperature value.

### ALARM

CODE 03 **F**

Actual temperature above high temperature value.

- Excess temperature warning or Excess temperature alarm

**Warn-Type:** Set to >Warning< or >Alarm<

### WARNING

CODE 04 **F**

Actual temperature below low temperature value.

### ALARM

CODE 04 **F**

Actual temperature below low temperature value.

- Low temperature warning or Low temperature alarm.

**Warn-Type:** Set to >Warning< or >Alarm<

### ALARM

CODE 05 **F**

Internal working sensor short-circuited or interrupted.

- Cable of the working temperature sensor interrupted or short-circuited.

**ALARM**

**CODE 06** **F** ↗

Difference between working- and safety-temp. Bigger than permissible value.

- Defect of the working or excess temperature sensor. Working temperature and excess temperature sensors report a temperature difference of more than 25 °C.

**ALARM**

**CODE 07** **F** ↗

The internal I2C bus-system shows a failure.

- Other errors (I<sup>2</sup>C-BUS errors)

**ALARM**

**CODE 12** **F** ↗

Error in A/D-converter of the internal measuring system.

- Error in A/D converter

**ALARM**

**CODE 14** **F** ↗

Working temperature exceeds the adjusted safety temperature.

- Excess temperature sensors defect.
- The excess temperature value lies below the working temperature setpoint. Set the excess temperature to a higher value.

**ALARM**

**CODE 15** **F** ↗

External control selected, but ext. sensor not connected or defective.

- External control selected, but external Pt100 sensor not connected or defect.

**WARNING**

**CODE 20** **F** ↗

Cooling is affected or ambient temperature to high  
Clean condenser

- Cooling of the condenser is affected. Clean air-cooled condenser. Check the flow rate and cooling water temperature on water-cooled condenser.

**WARNING**

**CODE 21** **F** ↗

Compressor does not work. Compressor will be automatically reconnected.

- Compressor does not work. After a short cooling interval, the compressor motor will be automatically reconnected and the message "CODE 21" no longer appears.

**ALARM**

**CODE 33** **F** ↗

Internal safety-temperature sensor short-circuited or interrupted.

- Cable of the excess temperature sensor interrupted or short-circuited

**ALARM**

**CODE 38** **F** ↗

Ext. Sensor input without signal but setpoint programming set to ext. Pt100.


- Ext. Pt100 sensor input without signal, but setpoint programming set to external Pt100.

**WARNING**

**CODE 40** **F** ↗

Niveau level warning  
Critical fluid level  
Replenish the bath tank with bath fluid

- The early warning system for low level signals a critical fluid level. Replenish the bath tank with the bath fluid.

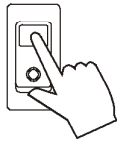


**Cancel the alarm state.**

Press the mains power switch off. After eliminating the malfunction, press the mains power on again to cancel the alarm state. If the unit cannot be returned to operation, contact an authorized service station.

**Disturbances that are not indicated.**

The electronic pump motor is overload-protected by an electronic current limiter. If viscosity of the bath fluid is or becomes too high, the motor stops running.

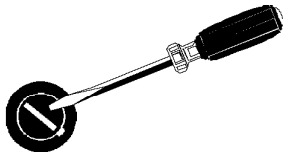

**Mains fuses:**

The mains fuses on the rear of the unit may easily be exchanged as shown on the left.  
Fine fuses (16a)– T 16 A, 250 V~ , D5 x 20 mm



**Warning:**

Before exchanging the fuses, turn off the mains power switch and disconnect the power plug from the mains socket!  
Only use fine fuses with a nominal value as specified.

Example:

Manufacturer	Supplier	Type	Order No.
Schurter	Schurter	G-fuse insert SPT T16A 5x20mm	No. 0001.2516

## 12. Electrical connections




**Notice:**

Use shielded cables only.

The shield of the connecting cable is electrically connected to the plug housing.

The unit ensures safe operation if connecting cables with a maximum length of 3 m are used. The use of longer cables does not affect proper performance of the unit, however external interferences may have a negative impact on safe operation.

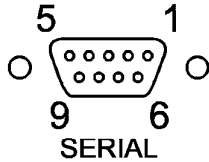

**RS232/RS485 serial interface**

This port can be used to connect a computer with an RS232 or RS485 cable for remote control of the Cryo-Compact Circulator .

**Pin assignments RS232:**

Pin 2	RxD	Receive Data
Pin 3	TxD	Transmit Data
Pin 4	DTR	Data terminal ready
Pin 5	0 V	Signal GND
Pin 7	RTS	Request to send
Pin 8	CTS	Clear to send

Pin 1; 6; 9      Reserved - do not use!

**Pin assignments RS485:**

Pin 3	B	
Pin 5	0 V	Signal GND
Pin 6	+5 V (50 mA max. load current)	
Pin 8	A	

Pins 1; 2; 7; 9 are reserved - do not use!




**Notice: RS485 serial interface**

The max. load current between Pin 5 and Pin 6 is 50 mA.


**Accessories:**

Order No.	Description
8 980 073	RS232 interface cable 9-pol./9-pol. , 2,5 m
8 900 110	USB interface adapter cable



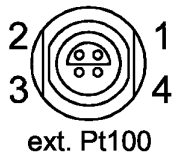


Socket for external Pt100 sensor

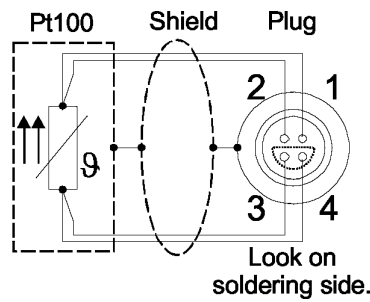

Pin assignment:

Pin	Signal
1	I+
2	U+
3	U-
4	I-

The shield of the connecting cable is electrically connected to the plug housing and the sensor tube.

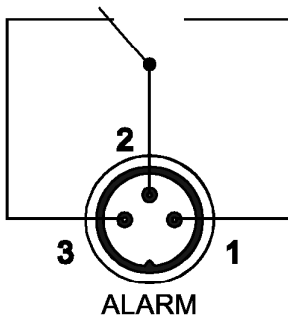


**Alarm output**

(for external alarm signal)

This potential-free change-over contact is activated in case of an alarm when pins 2 and 3 are connected.

Switching capacity	max.	30 W / 40 VA
Switching voltage	max.	125 V~/–
Switching current	max.	1 A

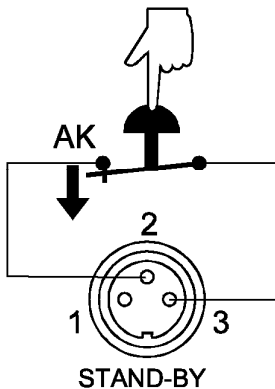


**STAND-BY input**

(for external emergency switch-off)

Pin assignment:	Pin	Signal
	1	not connected
	2	5 V / DC
	3	0 V

Functional description see page 52

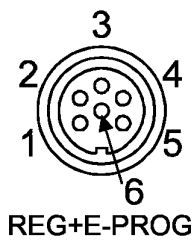


**Programmer input / temperature recorder output**

Pin	Signal
1 Voltage output	Channel 1 0 ... 10 V
2 Voltage output	Channel 2 0 ... 10 V
3 GND for outputs	0 V
4 Programmer input	EPROG 0 to 10 V / 0 to 20 mA
5 Current output	Channel 3 0 to 20 mA / 4 to 20 mA
6 GND for Programmer	0 V

Functional description see page 49

## 13. Remote control

### 13.1. Setup for remote control

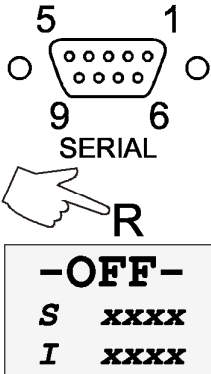

1. Check the interface parameters for both interfaces (on Cryo-Compact Circulator and PC) and make sure they match.  
In the menu >Interface< set the menu item >Type< to >RS232< or >RS485<.( Serial interface see page 56)
2. In the menu >Configuration< set the menu item >Setpoint< to >RS232< or >RS485<.  
(>Configuration< see page 30)
3. Connect both units with an interface cable..



Like all parameters which can be entered through the keypad, interface parameters are stored in memory even after the Cryo-Compact Circulator is turned off.

### 13.2. Communication with a PC or a superordinated data system

If the Cryo-Compact Circulator is put into remote control mode via the configuration level, the VFD COMFORT-DISPLAY will read „R -OFF-„ = REMOTE STOP. The Cryo-Compact Circulator is now operated via the computer.  
In general, the computer (master) sends commands to the Cryo-Compact Circulator (slave). The Cryo-Compact Circulator sends data (including error messages) only when the computer sends a query.



In remote control mode: After a power interruption the order to start and all values which have to be adjusted must be resent from the personal computer via the interface.  
AUTOSTART is not possible.

A transfer sequence consists of:

- address (RS485 interface only)
- command
- space (↔; Hex: 20)
- parameter (the character separating decimals in a group is the period)
- end of file (↵; Hex: 0D)
- The response (data string) after an in command is always followed by a line feed (LF, Hex: 0A).

#### Important times for a command transmission:



To ensure a safe data transfer, the time gap between two commands should be at least 250 ms.

The Cryo-Compact Circulator automatically responds to an in command with a data string followed by a LF (Line Feed). The next command should only be sent after 10 ms.

The commands are divided into **in** or **out** commands.

**in** commands: asking for parameters to be displayed

**out** commands: setting parameters



The **out** commands are valid only in remote control mode.

Examples:

When the RS485 interface is used, the three-digit instrument address stands in front of each command.

(example: address Ad32 = **A032** )

Command to set the working temperature >Setpoint1< to 55.5 °C

**out\_sp\_00** ↔ **55.5**↵

**A032\_out\_sp\_00** ↔ **55.5**↵

Command to ask for the working temperature >Setpoint1<

**in\_sp\_00**↵

**A032\_in\_sp\_00**↵

Response from the Cryo-Compact Circulator :

**55.5**↵ LF

**A032\_55.5**↵ LF



**Notice:** **out-commands**

Temperature values can be displayed on the Cryo-Compact Circulator in the unit °C or °F.

According to this setting the instructions via the interface also have to be made in the set unit.

### 13.3. List of commands

**out commands:** Setting parameters or temperature values.

Command	Parameter	Response of Cryo-Compact Circulator
out_mode_01	0	Use working temperature >Setpoint1<
out_mode_01	1	Use working temperature >Setpoint2<
out_mode_01	2	Use working temperature >Setpoint3<
out_mode_02	0	Selftuning „off“. Temperature control by using the stored parameters.
out_mode_02	1	Selftuning „once“ Single selftuning of controlled system after the next start.
out_mode_02	2	Selftuning „always“ Continual selftuning of controlled system whenever a new setpoint is to be reached.

## Remote control

Command	Parameter	Response of Cryo-Compact Circulator
out_mode_03	0	Set external programmer input to voltage. Voltage 0 V ... 10 V
out_mode_03	1	Set external programmer input to current. Current 0 mA ... 20 mA
out_mode_04	0	Temperature control of internal bath.
out_mode_04	1	External control with Pt100 sensor.
out_mode_05	0	Stop the unit = R -OFF-.
out_mode_05	1	Start the unit.
out_mode_08	0	Set the control dynamics - aperiodic
out_mode_08	1	Set the control dynamics - standard
out_sp_00	xxx.xx	Set working temperature. „Setpoint 1“
out_sp_01	xxx.xx	Set working temperature. „Setpoint 2“
out_sp_02	xxx.xx	Set working temperature. „Setpoint 3“
out_sp_03	xxx.xx	Set high temperature warning limit „OverTemp“
out_sp_04	xxx.xx	Set low temperature warning limit „SubTemp“
out_sp_06	xxx.xx	Set manipulated variable for the heater via serial interface -100 ... +100 [%]
out_sp_07	x	Set the pump pressure stage. (1 ... 4)
out_par_04	x.x	CoSpeed for external control 0 ... 5.0
out_par_06	xxx	Xp control parameter of the internal controller. 0.1 ... 99.9
out_par_07	xxx	Tn control parameter of the internal controller. 3 ... 9999
out_par_08	xxx	Tv control parameter of the internal controller. 0 ... 999
out_par_09	xxx	Xp control parameter of the cascade controller. 0.1 ... 99.9
out_par_10	xxx	Proportional portion of the cascade controller. 1 ... 99.9
out_par_11	xxx	Tn control parameter of the cascade controller. 3 ... 9999
out_par_12	xxx	Tv control parameter of the cascade controller. 0 ... 999
out_par_13	xxx	Maximum internal temperature of the cascade controller.
out_par_14	xxx	Minimum internal temperature of the cascade controller.
out_par_15	xxx	Band limit (upper) 0 ... 200 °C
out_par_16	xxx	Band limit (lower) 0 ... 200 °C
out_hil_00	-xxx	Set the desired maximum cooling power (0 % to 100 %). <b>Note:</b> Enter the value with a preceding negative sign! This adjustment only makes sense if CF41is used.
out_hil_01	xxx	Set the desired maximum heating power (10 % to 100 %).

**in commands:** Asking for parameters or temperature values to be displayed.

<b>Command</b>	<b>Parameter</b>	<b>Response of Cryo-Compact Circulator</b>
version	None	Number of software version (V X.xx)
status	none	Status message, error message (see page 70)
in_pv_00	none	Actual bath temperature.
in_pv_01	none	Heating power being used (%).
in_pv_02	none	Temperature value registered by the external Pt100 sensor.
in_pv_03	none	Temperature value registered by the safety sensor.
in_pv_04	none	Setpoint temperature („SafeTemp“) of the excess temperature protection
in_sp_00	none	Working temperature „Setpoint 1“
in_sp_01	none	Working temperature „Setpoint 2“
in_sp_02	none	Working temperature „Setpoint 3“
in_sp_03	none	High temperature warning limit „OverTemp“
in_sp_04	none	Low temperature warning limit „SubTemp“
in_sp_05	none	Setpoint temperature of the external programmer (REG+E-PROG) .
in_sp_06	none	Temperature indicatio in °C or °F
in_sp_07	none	1. Adjusted pump stage in the -OFF- condition. 2. Pump stage corresponding to the effective rotation speed after start.
in_sp_08	none	Value of a flowrate sensor connected to the E-Prog input
in_par_00	none	Temperature difference between working sensor and safety sensor
in_par_01	none	Te - Time constant of the external bath.
in_par_02	none	Si - Internal slope
in_par_03	none	Ti - Time constant of the internal bath.
in_par_04	none	CoSpeed - Band limit (max. difference between the temperatures in the internal bath and external system).
in_par_05	none	Factor pk/ph0: Ratio of max. cooling capacity versus max. heating capacity
in_par_06	none	Xp control parameter of the internal controller.
in_par_07	none	Tn control parameter of the internal controller.
in_par_08	none	Tv control parameter of the internal controller.
in_par_09	none	Xp control parameter of the cascade controller.
in_par_10	none	Proportional portion of the cascade controller.


## Remote control

Command	Parameter	Response of Cryo-Compact Circulator
in_par_11	none	Tn control parameter of the cascade controller.
in_par_12	none	Tv control parameter of the cascade controller.
in_par_13	none	Adjusted maximum internal temperature of the cascade controller.
in_par_14	none	Adjusted minimum internal temperature of the cascade controller.
in_par_15	none	Band limit (upper)
in_par_16	none	Band limit (lower)
in_mode_01	none	Selected setpoint: 0 = Setpoint 1 1 = Setpoint 2 2 = Setpoint 3
in_mode_02	none	Selftuning type: 0 = Selftuning „off“ 1 = Selftuning „once“ 2 = Selftuning „always“
in_mode_03	none	Type of the external programmer input: 0 = Voltage 0 V to 10 V 1 = Current 0 mA to 20 mA
in_mode_04	none	Internal/external temperature control: 0 = Temperature control with internal sensor. 1 = Temperature control with external Pt100 sensor.
in_mode_05	none	Cryo-Compact Circulator in Stop/Start condition: 0 = Stop 1 = Start
in_mode_08	none	Adjusted control dynamics 0 = aperiodic 1 = standard
in_hil_00	none	Max. cooling power (%).
in_hil_01	none	Max. heating power (%).

### 13.4. Status messages

Status messages	Description
<b>00 MANUAL STOP</b>	Cryo-Compact Circulator in „OFF“ state.
<b>01 MANUAL START</b>	Cryo-Compact Circulator in keypad control mode.
<b>02 REMOTE STOP</b>	Cryo-Compact Circulator in „r OFF“ state.
<b>03 REMOTE START</b>	Cryo-Compact Circulator in remote control mode.

### 13.5. Error messages

Error messages	Description
<b>-01 LOW LEVEL ALARM</b>	Low liquid level alarm
<b>-03 EXCESS TEMPERATURE WARNING</b>	High temperature warning
<b>-04 LOW TEMPERATURE WARNING</b>	Low temperature warning.
<b>-05 WORKING SENSOR ALARM</b>	Working temperature sensor short-circuited or interrupted.
<b>-06 SENSOR DIFFERENCE ALARM</b>	Sensor difference alarm. Working temperature and safety sensors report a temperature difference of more than 25 °C.
<b>-07 I<sup>2</sup>C-BUS ERROR</b>	Internal error when reading or writing the I <sup>2</sup> C bus.
<b>-08 INVALID COMMAND</b>	Invalid command.
<b>-09 COMMAND NOT ALLOWED IN CURRENT OPERATING MODE</b>	Invalid command in current operating mode.
<b>-10 VALUE TOO SMALL</b>	Entered value too small.
<b>-11 VALUE TOO LARGE</b>	Entered value too large.
<b>-12 TEMPERATURE MEASUREMENT ALARM</b>	Error in A/D converter.
<b>-13 WARNING : VALUE EXCEEDS TEMPERATURE LIMITS</b>	Value lies outside the adjusted range for the high and low temperature warning limits. But value is stored.
<b>-14 EXCESS TEMPERATURE PROTECTOR ALARM</b>	Excess temperature protector alarm 
<b>-15 EXTERNAL SENSOR ALARM</b>	External control selected, but external Pt100 sensor not connected.
<b>-20 WARNING: CLEAN CONDENSOR OR CHECK COOLING WATER CIRCUIT OF REFRIGERATOR</b>	Cooling of the condenser is affected. Clean air-cooled condenser. Check the flow rate and cooling water temperature on water-cooled condenser.
<b>-21 WARNING: COMPRESSOR STAGE 1 DOES NOT WORK</b>	Compressor stage 1 does not work.
<b>-26 WARNING: STAND-BY PLUG IS MISSING</b>	External standby contact is open. (see page 52 and 65)
<b>-33 SAFETY SENSOR ALARM</b>	Excess temperature sensor short-circuited or interrupted.
<b>-38 EXTERNAL SENSOR SETPOINT PROGRAMMING ALARM</b>	Ext. Pt100 sensor input without signal and setpoint programming set to external Pt100.
<b>-40 NIVEAU LEVEL WARNUNG</b>	Low liquid level warning in the internal reservoir.

## 14. JULABO Service – Online remote diagnosis

JULABO Cryo-Compact Circulator s of the HighTech series are equipped with a so-called black box. This box is implemented in the controller and records all significant data for the last 30 minutes.

In case of a failure, this data can be read out from the unit by using special software. The respective program is available for **free** download from [www.julabo.de \ EasyBlackBox](http://www.julabo.de \ EasyBlackBox).

- Installation is easy and carried out step by step. Please observe the instructions.
- Data read-out is possible in the conditions „OFF“, „R OFF“ or „ALARM“.
- Connect the Cryo-Compact Circulator to the computer using an interface cable.
- Start the EasyBlackBox program. The program asks for the used port (COM1, ..... ) and the baud rate of the unit. You do not have this information on hand? Simply try it out! The program keeps on sending this request until the actually used port and correct baud rate are entered.



- Data is read out and shown on the monitor divided in the sections >Einstellungen/Settings<, >Alarmspeicher/Alarms stored<, >Blackbox<   
 ← see example
- After pressing >Speichern/Save< a text file is compiled. The program proposes a filename - >C:\model description and barcode no.<. Modifications are possible.
- E-mail this file to [service@julabo.de](mailto:service@julabo.de), JULABO's service department. JULABO is thus able to provide rapid support.




## 15. Installation of electronic module



### Caution:

Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit. Configuration, installation, maintenance and repairs on the Cryo-Compact Circulator may only be carried out by qualified personnel.

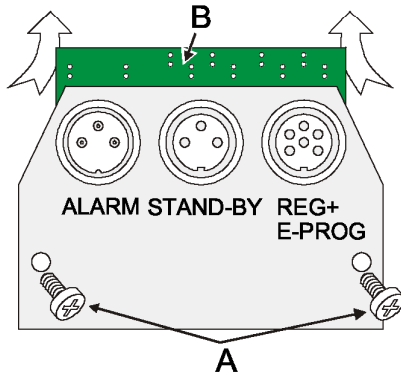


- Turn off the unit with the mains switch and disconnect the power plug.
- Open the screws (A) and remove the dummy plate.
- Carefully insert the upper edge (B) into the rear panel from below, align the electronic module at top and bottom and applying light pressure connect the 15-pole connector.
- Attach the electronic module using the screws (A).
- The Cryo-Compact Circulator is ready for use. It automatically recognizes the installed electronic module.

## 16. Maintaining the cooling performance



### Caution:

Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit. The opening of the Cryo-Compact Circulator must be performed by qualified personnel only.

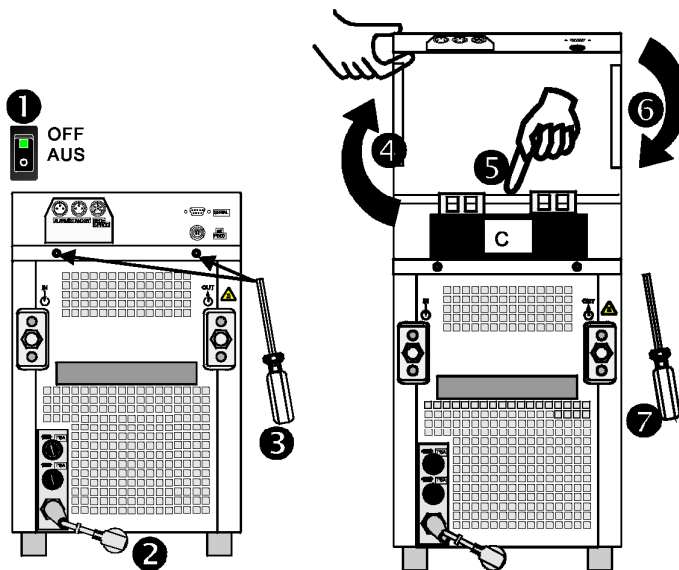

To maintain the full cooling performance, clean the condenser ( C ) from time to time.

1. Switch off device by pressing the main power switch and
2. disconnect mains cable from power source.
3. Remove 2 screws
4. Lift cover upwards.
5. Remove dirt at condenser by suction cleaning.
6. Close cover and
7. Fix by means of screws.
8. Unit is ready for operation.

## 16.1. Draining




**Notice:**

Exercise caution when emptying hot or cold bath fluids!

Check the temperature of the bath fluid prior to draining (by switching the unit on for a short moment, for example).

Store and dispose the used bath fluid according to the laws for environmental protection.

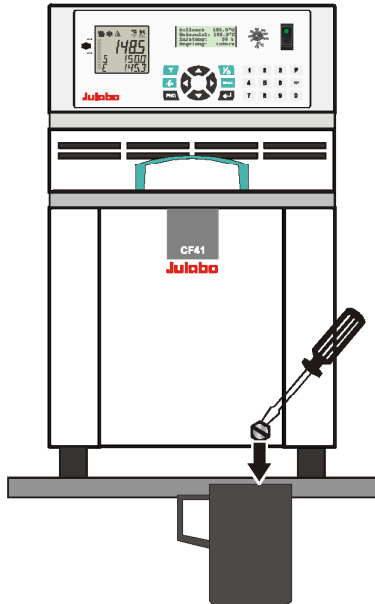

### Draining

- Turn off the unit and disconnect the mains cable from the power source.
- Place the Cryo-Compact Circulator near the rim of the table. Use a suitable vessel as recipient for the bath liquid
- Unscrew the drain tap and empty the unit completely.
- Tighten the drain tap.

## 17. Cleaning / repairing the unit

**Caution:**

Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.

Prevent humidity from entering into the Cryo-Compact Circulator .

Electrical connections and any other work must be performed by qualified personnel only.

**Cleaning:**

Clean the outside of the unit using a wet cloth and low surface tension water.

The Cryo-Compact Circulator is designed for continuous operation under normal conditions. Periodic maintenance is not required.

The tank should be filled only with a bath fluid recommended by JULABO. To avoid contamination, it is essential to change the bath fluid from time to time.

**Repairs:**

Before asking for a service technician or returning a JULABO instrument for repair, please contact an authorized JULABO service station.

When returning the unit:

- Clean the unit in order to avoid any harm to the service personnel
- Attach a short fault description.  
If you intend to return your JULABO unit to us, you will find a Service Return Form on our website [www.julabo.de](http://www.julabo.de) . Please use this as a delivery note and include it to the unit or send it in advance either by Fax or E-Mail.
- When returning a unit, take care of careful and adequate packing.
- JULABO is not responsible for damages that might occur from insufficient packing.



JULABO reserves the right to carry out technical modifications with repairs for providing improved performance of a unit.

## 18. Warranty conditions

JULABO Labortechnik GmbH warrants its products against defects in material or in workmanship, when used under appropriate conditions and in accordance with appropriate operating instructions

**for a period of ONE YEAR.**

**Extension of the warranty period – free of charge**



With the '1PLUS warranty' the user receives a free of charge extension to the warranty of up to 24 months, limited to a maximum of 10 000 working hours.

To apply for this extended warranty the user must register the unit on the JULABO web site [www.julabo.de](http://www.julabo.de), indicating the serial no. The extended warranty will apply from the date of JULABO Labortechnik GmbH's original invoice.

JULABO Labortechnik GmbH reserves the right to decide the validity of any warranty claim. In case of faults arising either due to faulty materials or workmanship, parts will be repaired or replaced free of charge, or a new replacement unit will be supplied.

Any other compensation claims are excluded from this guarantee.