## Super Clean<sup>™</sup> Gas Filters





- High Capacity
- High-purity output ensures
   99.9999% Pure gas
- "Quick connect" fittings for easy, leak-tight cartridge changes
- Glass inside to prevent diffusion, plastic outside for safety
- Easy to read indicators

#### **ABOUT SUPER CLEAN™ GAS FILTERS**

Cartridge systems make changing gas filters quick and easy. A base plate allows cartridges to be exchanged without introducing oxygen. Spring-loaded check valves seal when a filter is removed and open only when a new filter has been locked in place. There is no longer a need for loosening and tightening fittings every time a trap is changed and your system will not become contaminated during the process

#### WHY GAS PURITY IS IMPORTANT

Carrier gas should contain less than 1 ppm of oxygen, moisture, or other trace contaminants, to prevent column degradation, increase column lifetime, and decrease stationary phase bleed. The expense of using high-purity gases in combination with carrier gas line purifiers will be offset by longer column lifetime and less GC maintenance.

Contaminants cause ghost peaks to appear during temperature programming and degrade the validity of analytical data. Make-up gas also should be contaminant-free, or baseline fluctuations and excessive detector noise can occur; detector gases should be free of water and hydrocarbons, or excessive baseline noise can result. Gas purifiers remove these contaminants from gas sources, thereby improving system performance.

#### GAS FILTER SYSTEM COMPONENTS



#### **Base plate**

Connecting unit with in- and outlet connectors for the gas line and two spring-loaded check valves that automatically start the flow of gas once is filter is installed.



#### **Universal Ring Nut**

The Universal Ring Nut is used for mounting a filter cartridge or flush-cap to a base plate. It can also serve as a mounting point for the Electronic Maintenance Indicator device.



#### Filter Cartridge

The filter cartridges are made of glass to prevent diffusion, and protected by a plastic housing for safety. The PTFE seals at the base of the Filter will only be punctured during installation on the base plate.



#### **Electronic Indicator**

Optional Electronic Maintenance Indicator device warns when scheduled replacement or maintenance is due.



#### Wall mount bracket set

Optional wall mounting brackets to mount a Base plate on a wall.




#### GC / MS (ITD/MSD)




Product	Part No.	Qty
Triple Filter	SGT-F0301	1
1 Position Base plate	SGT-B0010-XX	1

#### GC / TCD

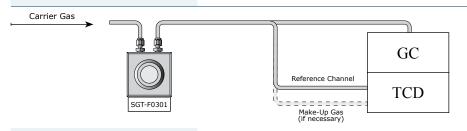

Product	Part No.	Qty
Triple Filter	SGT-F0301	1
1 Position Base plate	SGT-B0010-XX	1

#### GC / ECD

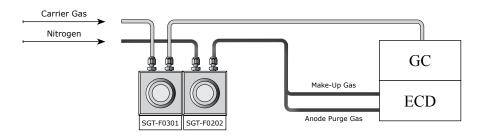


Product	Part No.	Qty
Triple Filter	SGT-F0301	1
Oxygen/Moisture Filter	SGT-F0202	1
2 Position Base plate	SGT-B0020-XX	1

#### **ICP-OES / ICP-MS**



Product	Part No.	Qty
High Flow Oxygen/ Moisture Filter Bundle	SGT-F0724	1
Particle Filter	B0060	1
High Flow Base plate	SGT-B0021-XX	1


#### **ICP-OES**

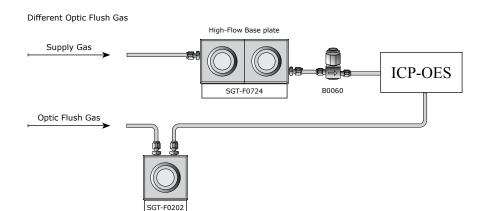

Product	Part No.	Qty
High Flow Oxygen/ Moisture Filter Bundle	SGT-F0724	1
Oxygen/Moisture Filter	SGT-F0202	1
Partile Filter	B0060	1
High Flow Base plate	SGT-B0021-XX	1
1 Position Base plate	SGT-B0010-XX	1


#### GC / FID

GC/FID Solution 1 (Carrier Gas = Make-Up Gas)

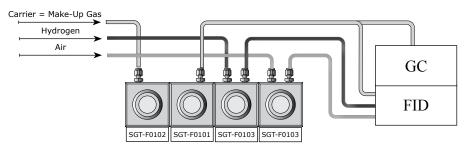


Product	Part No.	Qty
Oxygen Filter	SGT-F0102	1
Moisture Filter	SGT-F0101	1
Hydrocarbon Filter	SGT-F0103	2
4 Position Base plate	SGT-B0040-XX	1

GC/FID Solution 2 (Carrier Gas = Make-Up Gas)

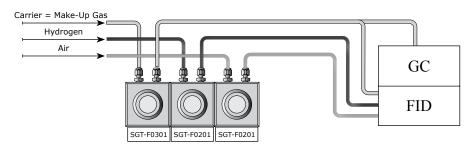

Product	Part No.	Qty
Triple Filter	SGT-F0301	1
Hydrocarbon/Moisture Filter	SGT-F0201	2
3 Position Base plate	SGT-B0030-XX	1


GC/FID Solution 3 (Carrier Gas differs from Make-Up Gas)

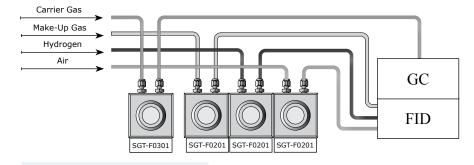

Product	Part No.	Qty
Triple Filter	SGT-F0301	1
1 Position Base plate	SGT-B0010-XX	1


Product	Part No.	Qty
Hydrocarbon/Moisture Filter	SGT-F0201	3
3 Position Base plate	SGT-B0030-XX	1

#### GC (FPD/PFPD)

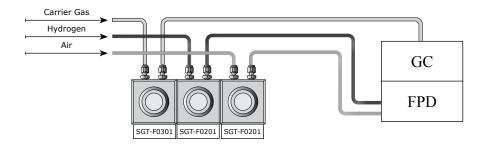


Product	Part No.	Qty
Triple Filter	SGT-F0301	1
Hydrocarbon/Moisture Filter	SGT-F0201	2
3 Position Base plate	SGT-B0030-XX	1

#### GC (TSD/TID/NPD/PND)

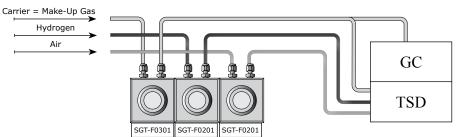


Product	Part No.	Qty
Triple Filter	SGT-F0301	1
Hydrocarbon/Moisture Filter	SGT-F0201	2
3 Position Base plate	SGT-B0030-XX	1

#### LC/MS

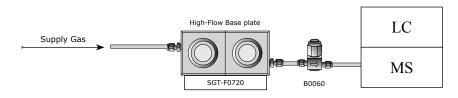

Product	Part No.	Qty
High Flow Hydrocarbon Filter Bundle	SGT-F0720	1
Particle Filter	B0060	1
High Flow Base plate	SGT-B0021-XX	1

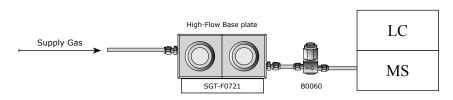

Part No.	Qty
SGT-F0721	1
B0060	1
SGT-B0021-XX	1
	SGT-F0721 B0060




Product	Part No.	Qty
High Flow Hydrocarbon/ Moisture Filter Bundle	SGT-F0723	1
Particle Filter	B0060	1
High Flow Base plate	SGT-B0021-XX	1

# Carrier Gas TOC Analyzer

SGT-F0205

Product	Part No.	Qty
CO <sub>2</sub> / Moisture Filter	SGT-F0205	1
1 Position Base plate	SGT-B0010-XX	1

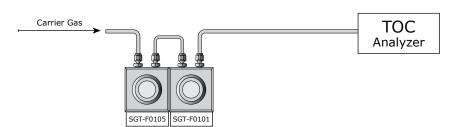


Product	Part No.	Qty
CO <sub>2</sub> Filter	SGT-F0105	1
Moisture Filter	SGT-F0101	1
2 Position Base plate	SGT-B0020-XX	1

#### **Zero-Air Generator**



Product	Part No.	Qty
CO <sub>2</sub> / Moisture Filter	SGT-F0205	1
1 Position Base plate	SGT-B0010-XX	1

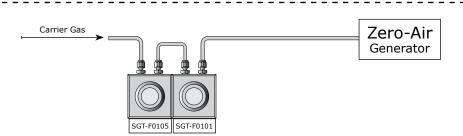

Product	Part No.	Qty
CO <sub>2</sub> Filter	SGT-F0105	1
Moisture Filter	SGT-F0101	1
2 Position Base plate	SGT-B0020-XX	1

Catalog No. *	SGT-B1010-B8 SGT-B1011-B8 (Helium) SGT-B1012-B8 (Hydrogen)
Usable for	Benefit
GC/MS	Higher data accuracy and less maintenance
GC/ECD	Greater sensitivity
GC/TCD	Greater sensitivity and less maintenance



#### TRIPLE FILTER KIT

The triple combination filter kit is ideal for purifying GC/MS carrier gases. It contains oxygen, moisture and hydrocarbon scrubbers in one easy to change economical cartridge.

Catalog No. *	SGT-B1015-B8
Usable for	Benefit
GC/TCD	Greater sensitivity and less maintenance



#### COMBI (OXYGEN/MOISTURE) KIT

The oxygen/moisture combination filter kit is ideal for purifying GC/ECD carrier gases. It contains oxygen and moisture scrubbers in one easy to change economical cartridge.

Catalog No. *	SGT-B1014-B8
Usable for	Benefit
TOC Analyzer	Greater sensitivity
Zero-Air generator	Cleaner gas



## COMBI (CO2/MOISTURE) FILTER KIT

The carbon dioxide/moisture combination filter kit is ideal for TOC analysers and Zero-Air generators. It removes both carbon dioxide and moisture from gas streams.

Catalog No. *	SGT-B1030-B8
Usable for	Benefit
GC/FID	Greater sensitivity



#### **3 POSITION KIT**

The FID Filter kit is the perfect all-in-one solution for purifying flame ionization detector (FID) fuel gases together with the carrier gas.

This kit removes hydrocarbons, moisture and oxygen from the carrier gas and removes both moisture and hydrocarbons from the Hydrogen and Air fuel gases.

Catalog No. *	SGT-B1040-B8
Usable for	Benefit
GC/FID	Greater sensitivity



#### **4 POSITION KIT**

The high capacity FID Filter kit is the perfect all-in-one solution for purifying flame ionization detector (FID) fuel gases together with the carrier gas.

This kit removes both moisture and oxygen from the carrier gas and removes hydrocarbons from the Hydrogen and Air fuel gases.

<sup>\*</sup> Standard fittings are 1/8" Brass. All Kits are available in 1/4" to 1/8" Brass and Stainless steel fittings.

To order, replace Part No. suffix into '-B4' for 1/4" Brass, '-S8' for 1/8" Stainless steel and '-S4' for 1/4" Stainless steel fittings.

Catalog No. \*\*

**Usable for** 

LC/MS

COMBI (MOISTURE/HYDRO-**CARBON) FILTERS KIT** 

The Fuel Gas Filter kit is perfect for

detector (FID) fuel gases, removing both moisture and hydrocarbons.

Using the Fuel Gas Filter for FID Hydrogen and air will produce a stable baseline, improving overall reproducibility and sensitivity.

purifying flame ionization

Catalog No. *	SGT-B1020-B8
Usable for	Benefit
FID	Greater sensitivity

SGT-B1021-B4

Greater sensitivity

**Benefit** 



### **HYDROCARBON KIT** Up to 20 Liter of hydrocarbon-free



Catalog No. **	SGT-B1022-B4	
Usable for	Benefit	



**MOISTURE KIT** Up to 20 Liter of moisture-free nitrogen per minute with this LC/MS High Flow Kit.

**HIGH FLOW** 





**HIGH FLOW** COMBI (HYDROCARBON/ **MOISTURE) KIT** Removes Moisture and

Hydrocarbons from high flow gas streams.

Catalog No. **	SGT-B1024-B4  Benefit	
Usable for		
ICP-MS	Greater sensitivity	
ICP-OES	Greater sensitivity	



**HIGH FLOW** COMBI (OXYGEN/MOISTURE)

Removes Oxygen and Moisture from high flow gas streams.

Standard fittings are 1/8" Brass. All Kits are available in 1/4" to 1/8" Brass and Stainless steel fittings. To order, replace Part No. suffix into '-B4' for 1/4" Brass, '-S8' for 1/8" Stainless steel and '-S4' for 1/4" Stainless steel fittings.

<sup>\*\*</sup> Standard fittings are 1/4" Brass. All Kits are available in 1/4" Brass and Stainless steel fittings. To order, replace Part No. suffix into '-S4' for 1/4" Stainless steel fittings.

Catalog No.	SGT-F0101
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	7 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR, Air
Dimensions	24 cm x Ø 4.4 cm
Weight	0.26 Kg
Estimated Lifetime	> 2 years



#### **MOISTURE FILTER**

Moisture in carrier gas lines will prematurely degrade oxygen and hydrocarbon traps and increase detector noise.

As a precaution, we highly recommend installing a moisture trap before the hydrocarbon and oxygen traps on all carrier gas lines.

_			
Ca	pa	CI	t۷

H <sub>2</sub> O	72σ	
п2О	7.2 g	

Catalog No.	SGT-F0102
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	7 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR
Dimensions	24 cm x Ø 4.4 cm
Weight	0.26 Kg
Estimated Lifetime	> 2 years




#### **OXYGEN FILTER**

Oxygen is a column killer. It is present even in UHP gases, as minutes leaks at fittings allow oxygen to influx against the concentration gradient. Because oxygen can enter a gas line at any fitting or during gas bottle exchange, the oxygen trap should be the last connection before the gas line enters the chromatograph.

Capacity	
----------	--

	-		
$O_2$		150 mL	

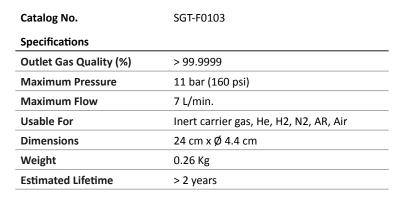




#### **HYDROCARBON FILTER**

Use a hydrocarbon trap if your gas has a potential source of hydrocarbon contaminants or if you suspect you are observing carrier gas ghost peaks.

Install the hydrocarbon trap after the moisture trap, to prevent moisture from degrading the hydrocarbon-trapping ability of the activated carbon in the hydrocarbon trap.

#### Capacity

HC	12 g (as <i>n</i> -butane)
----	----------------------------

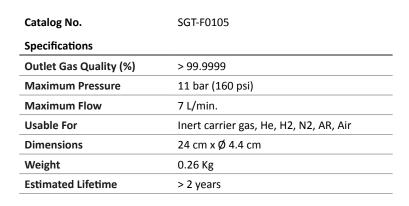



#### **CARBON DIOXIDE FILTER**

Removes carbon dioxide and sulfur from gas streams. To be used in combination with a Moisture Filter.

•	•		
co <sub>2</sub>		12 g	

Catalog No.	SGT-F0201
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	7 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR, Air
Dimensions	24 cm x Ø 4.4 cm
Weight	0.26 Kg
Estimated Lifetime	> 2 years



#### COMBI (HYDROCARBON/ MOISTURE) FILTER

The Fuel Gas Filter is perfect for purifying flame ionization detector (FID) fuel gases, removing both moisture and hydrocarbons. Using the Fuel Gas Filter for FID Hydrogen and air will produce a stable baseline, improving overall reproducibility and sensitivity.

Capacity	
H <sub>2</sub> O	3.5 g
НС	6 g (as <i>n</i> -butane)

Catalog No.	SGT-F0202
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	7 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR
Dimensions	24 cm x Ø 4.4 cm
Weight	0.26 Kg
Estimated Lifetime	> 2 years



## COMBI (OXYGEN/MOISTURE) FILTER

This Combi trap is ideal for purifying carrier gas. It contains oxygen and moisture scrubbers in one easy to change economical cartridge.

Capacity	
H <sub>2</sub> O	3.5 g
02	75 mL

Catalog No.	SGT-F0205
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	7 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR, Air
Dimensions	24 cm x Ø 4.4 cm
Weight	0.26 Kg
Estimated Lifetime	> 2 years



#### COMBI (CARBON DIOXIDE/ MOISTURE) FILTER

Removes carbon dioxide, sulfur and moisture from gas streams.

Capacity	
H <sub>2</sub> O	3.5 g
co <sub>2</sub>	6 g

Catalog No.	SGT-F0301 SGT-F0302 (Helium Specfic) SGT-F0304 (Hydrogen Specific)
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	7 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR
Dimensions	24 cm x Ø 4.4 cm
Weight	0.26 Kg
Estimated Lifetime	> 2 years



#### TRIPLE (OXYGEN/MOISTURE/ HYDROCARBON) FILTER

The Triple trap is ideal for purifying carrier gas. It contains oxygen, moisture and hydrocarbon scrubbers in one easy to change economical cartridge.

Capacity	
H <sub>2</sub> O	1.8 g
02	75 mL
НС	4 g (as <i>n</i> -butane)

Catalog No.	SGT-F0720
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	20 L/min.
Usable For	Nitrogen
Dimensions	24 cm x Ø 4.4 cm
Weight	0.5 Kg
Estimated Lifetime	3 to 6 months



## HIGH FLOW HYDROCARBON FILTER BUNDLE

Up to 20 Liter of hydrocarbon-free nitrogen per minute. To be used in combination with a high flow base plate.

#### Capacity

нс	24 g (as <i>n</i> -butane)
пс	24 g (as ii-butaile)

Catalog No.	SGT-F0721
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	20 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR, Air
Dimensions	24 cm x Ø 4.4 cm
Weight	0.5 Kg
Estimated Lifetime	3 to 6 months



## HIGH FLOW MOISTURE FILTER BUNDLE

Removes moisture from high flow gas streams. To be used in combination with a high flow base plate.

#### Capacity

Catalog No.	SGT-F0723
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	20 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR, Air
Dimensions	24 cm x Ø 4.4 cm
Weight	0.5 Kg
Estimated Lifetime	3 to 6 months



#### HIGH FLOW COMBI (HYDRO-CARBON/MOISTURE) FILTER BUNDLE

Removes Moisture and Hydrocarbons from high flow gas streams. To be used in combination with a high flow base plate.

#### Capacity

HC	12 g (as <i>n</i> -butane)
H <sub>2</sub> O	7.2 g

Catalog No.	SGT-F0724
Specifications	
Outlet Gas Quality (%)	> 99.9999
Maximum Pressure	11 bar (160 psi)
Maximum Flow	20 L/min.
Usable For	Inert carrier gas, He, H2, N2, AR
Dimensions	24 cm x Ø 4.4 cm
Weight	0.5 Kg
Estimated Lifetime	3 to 6 months



#### HIGH FLOW COMBI (OXYGEN/ MOISTURE) FILTER BUNDLE

Removes Oxygen and Moisture from high flow gas streams. To be used in combination with a high flow base plate.

#### Capacity

H <sub>2</sub> O	7.2 g
O <sub>2</sub>	150 mL

Catalog No.	Fitting Type
SGT-B0010-B8	1/8" Brass
SGT-B0010-B4	1/4" Brass
SGT-B0010-S8	1/8" SS
SGT-B0010-S4	1/4" SS




#### **1 POSITION BASE PLATE**

Single position base plate



Catalog No.	Fitting Type
SGT-B0020-B8	1/8" Brass
SGT-B0020-B4	1/4" Brass
SGT-B0020-S8	1/8" SS
SGT-B0020-S4	1/4" SS

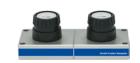

#### **2 POSITION BASE PLATE**

Double position base plate




Catalog No.	Fitting Type
SGT-B0030-B8	1/8" Brass
SGT-B0030-B4	1/4" Brass
SGT-B0030-S8	1/8" SS
SGT-B0030-S4	1/4" SS



#### **3 POSITION BASE PLATE**

Three position base plate

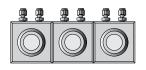

Catalog No.	Fitting Type
SGT-B0040-B8	1/8" Brass
SGT-B0040-B4	1/4" Brass
SGT-B0040-S8	1/8" SS
SGT-B0040-S4	1/4" SS



#### **4 POSITION BASE PLATE**

Four position base plate




Catalog No.	Fitting Type
SGT-B0021-B4	1/4" Brass
SGT-B0021-S4	1/4" SS



#### **HIGH FLOW BASE PLATE**

High Flow double position base plate (in parallel)

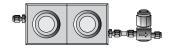

Catalog No.	Clip Type
SGT-E0040	Filter Cartridge Clip
SGT-E0010	Universal Ring Nut Clip
Usable for	
All base plates with Universal Ring Nut	





#### **ELECTRONIC INDICATOR**

Electronic Maintenance Indicator device warns when scheduled filter replacement or base plate maintenance is due.

Catalog No.	B0140
Usable for	
All base plates	



#### **UNIVERSAL RING NUT**

Universal Ring Nut to mount a filter or flush-cap on a base plate.

All base plates		
Catalog No.	B0110	
Usable for		
All base plates		



#### **O-RING REPLACEMENT SET**

A set of 10 + 10 replacement O-rings for replacing the O-rings on the in- and outlet valves on a base plate.





#### **WALL-MOUNT BRACKET SET**

Wall mounting brackets to mount a Base plate on a wall.

Catalog No.	B0130
Usable for	
All base plates	



#### **UNIVERSAL FLUSH-CAP SET**

Flush-cap that mounts on a base plate, and allows the gas to pass through the base plate without a filter attached.

Catalog No. Connection Typ		
B0120	1/4" Brass	
B0121	1/8" Brass	
B0122	1/4" SS	
B0123	1/8" SS	
Usable for		
1, 2, 3 and 4 position base plates		



## STANDARD BASE PLATE CONNECTOR SET

Replacement connectors for standard base plates.

Catalog No.	Connection Type
B0124	1/4" Brass
B0125	1/4" SS
Usable for	
High flow base plates	



#### HIGH FLOW BASE PLATE CONNECTOR SET

Replacement connector for high flow base plates.

Catalog No.	Description
B0060	0.5 Micron Particle Filter (1/4" Brass)
B0065	0.5 Micron Particle Filter Cup

Replacement Pack



#### **PARTICULATE FILTER**

Particulate filter for high flow base plates.

#### **Installation Kits**

Description	qty.
SGT 1 Triple Filter Kit - $^1/8''$ Brass Includes (1) 1 position base plate with $^1/8''$ Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Trap	ea.
SGT 1 Triple Helium Specific Filter Kit - $^{1}$ /8" Brass Includes (1) 1 position base plate with $^{1}$ /8" Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Helium Specific Trap	ea.
SGT 1 Triple Hydrogen Specific Filter Kit - $^{1}/_{8}$ " Brass Includes (1) 1 position base plate with $^{1}/_{8}$ " Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Hydrogen Specific Trap	ea.
SGT Co2 Filter Kit - $^1$ /8" Brass Includes (1) 1 position base plate with $^1$ /8" Brass inlet/outlet fittings and (1) Carbon dioxide Trap	ea.
SGT Combi (Co2/Moisture) Filter Kit - 1/8" Brass Includes (1) 1 position base plate with 1/8" Brass inlet/outlet fittings and (1) Carbon dioxide/Moisture Trap	ea.
SGT Combi (Oxygen/Moisture) Kit - $^1/_8$ " Brass Includes (1) 1 position base plate with $^1/_8$ " Brass inlet/outlet fittings and (1) Oxygen/Moisture Trap	ea.
SGT 2 Combi Filters Kit - $^{1}/_{8}$ " Brass Includes (1) 2 position base plate with $^{1}/_{8}$ " Brass inlet/outlet fittings and (2) Hydrocarbon/Moisture Traps	ea.
SGT 3 Filters Kit - $^1$ /8" Brass Includes (1) 3 position base plate with $^1$ /8" Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Trap and (2) Hydrocarbon/Moisture Traps	ea.
SGT 4 Filters Kit - $^1$ /8" Brass Includes (1) 4 position base plate with $^1$ /8" Brass inlet/outlet fittings and (1) Moisture Trap, (1) Oxygen Trap, and (2) Hydrocarbon Traps	ea.
	SGT 1 Triple Filter Kit - 1/8" Brass Includes (1) 1 position base plate with 1/8" Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Trap  SGT 1 Triple Helium Specific Filter Kit - 1/8" Brass Includes (1) 1 position base plate with 1/8" Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Helium Specific Trap  SGT 1 Triple Hydrogen Specific Filter Kit - 1/8" Brass Includes (1) 1 position base plate with 1/8" Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Hydrogen Specific Trap  SGT Co2 Filter Kit - 1/8" Brass Includes (1) 1 position base plate with 1/8" Brass inlet/outlet fittings and (1) Carbon dioxide Trap  SGT Combi (Co2/Moisture) Filter Kit - 1/8" Brass Includes (1) 1 position base plate with 1/8" Brass inlet/outlet fittings and (1) Carbon dioxide/Moisture Trap  SGT Combi (Oxygen/Moisture) Kit - 1/8" Brass Includes (1) 1 position base plate with 1/8" Brass inlet/outlet fittings and (1) Oxygen/Moisture Trap  SGT 2 Combi Filters Kit - 1/8" Brass Includes (1) 2 position base plate with 1/8" Brass inlet/outlet fittings and (2) Hydrocarbon/Moisture Traps  SGT 3 Filters Kit - 1/8" Brass Includes (1) 3 position base plate with 1/8" Brass inlet/outlet fittings and (1) Oxygen/Moisture/Hydrocarbon Trap and (2) Hydrocarbon/Moisture Traps  SGT 4 Filters Kit - 1/8" Brass Includes (1) 4 position base plate with 1/8" Brass inlet/outlet fittings and (1)

#### **High Flow Installation Kits**

Part No. **	Description	qty.
SGT-B1021-B4	SGT High Flow Hydrocarbon Filter Kit - $^1/_4$ " Brass Includes (1) 2 position high flow base plate with $^1/_4$ " Brass inlet/outlet fittings, (2) High Flow Hydrocarbon Traps, and a $^1/_4$ " Brass Particle Filter	ea.
SGT-B1022-B4	SGT High Flow Moisture Filter Kit - $^1/_4$ " Brass Includes (1) 2 position high flow base plate with $^1/_4$ " Brass inlet/outlet fittings, (2) High Flow Moisture Traps, and a $^1/_4$ " Brass Particle Filter	ea.
SGT-B1023-B4	SGT High Flow Combi (Hydrocarbon/Moisture) Filter Kit - $^1$ / $_4$ " Brass Includes (1) 2 position high flow base plate with $^1$ / $_4$ " Brass inlet/outlet fittings, (2) High Flow Combi (Hydrocarbon/Moisture) Traps, and a $^1$ / $_4$ " Brass Particle Filter	ea.
SGT-B1024-B4	SGT High Flow Combi (Oxygen/Moisture) Filter Kit - $^1/_4$ " Brass Includes (1) 2 position high flow base plate with $^1/_4$ " Brass inlet/outlet fittings, (2) High Flow Combi (Oxygen/Moisture) Traps, and a $^1/_4$ " Brass Particle Filter	ea.

<sup>\*</sup> Standard fittings are 1/8" Brass. All Kits are available in 1/4" to 1/8" Brass and Stainless steel fittings.

To order, replace Part No. suffix into '-B4' for 1/4" Brass, '-S8' for 1/8" Stainless steel and '-S4' for 1/4" Stainless steel fittings.

<sup>\*\*</sup> Standard fittings are 1/4" Brass. All Kits are available in 1/4" Brass and Stainless steel fittings. To order, replace Part No. suffix into '-S4' for 1/4" Stainless steel fittings.

#### Replacement Filters

Part No.	Description	qty.
SGT-F0101	SGT Moisture Filter	ea.
SGT-F0102	SGT Oxygen Filter	ea.
SGT-F0103	SGT Hydrocarbon Filter	ea.
SGT-F0105	SGT Carbon dioxide Filter	ea
SGT-F0201	SGT Combi (Hydrocarbon/Moisture) Filter	ea.
SGT-F0202	SGT Combi (Oxygen/Moisture) Filter	ea
SGT-F0205	SGT Combi (Carbon dioxide/Moisture Filter)	ea
SGT-F0301	SGT Triple (Oxygen/Moisture/Hydrocarbon) Filter	ea.
SGT-F0302	SGT Triple (Oxygen/Moisture/Hydrocarbon) Helium-specific Filter	ea.
SGT-F0304	SGT Triple (Oxygen/Moisture/Hydrocarbon) Hydrogen-specific Filter	ea.

#### **Replacement Filter Bundles**

Part No.	Description	qty.
SGT-F0730	SGT Filter Bundle of 3 Includes (1) Oxygen/Moisture/Hydrocarbon Trap and (2) Hydrocarbon/ Moisture Traps	ea.
SGT-F0740	SGT Filter Bundle of 4 Includes (1) Moisture Trap, (1) Oxygen Trap, and (2) Hydrocarbon Traps	ea.

#### High Flow Replacement Filter Bundles

Part No.	Description	qty.
SGT-F0720	SGT High Flow Hydrocarbon Filter Bundle Includes (2) High Flow Hydrocarbon Filters	ea.
SGT-F0721	SGT High Flow Moisture Filter Bundle Includes (2) High Flow Moisture Filters	ea.
SGT-F0723	SGT High Flow Combi (Hydrocarbon/Moisture) Filter Bundle Includes (2) High Flow Combi (Hydrocarbon/Moisture) Filters	ea.
SGT-F0724	SGT High Flow Combi (Oxygen/Moisture) Filter Bundle Includes (2) High Flow Combi (Oxygen/Moisture) Filters	ea.

#### Base plates

Part No. *	Description	qty.
SGT-B0010-B8	SGT Base plate 1 Position - 1/8" Brass	ea.
SGT-B0020-B8	SGT Base plate 2 Position - 1/8" Brass	ea.
SGT-B0030-B8	SGT Base plate 3 Position - 1/8" Brass	ea.
SGT-B0040-B8	SGT Base plate 4 Position - 1/8" Brass	ea.

#### High Flow Base plates

Part No. **	Description	qty.
SGT-B0021-B4	SGT High Flow Base plate 2 Position - $^1/_4$ " Brass Includes $^1/_4$ " Brass Particle Filter	ea.

<sup>\*</sup> Standard fittings are 1/8" Brass. All base plate are available in 1/4" to 1/8" Brass and Stainless steel fittings.

To order, replace Part No. suffix into '-B4' for 1/4" Brass, '-S8' for 1/8" Stainless steel and '-S4' for 1/4" Stainless steel fittings.

<sup>\*\*</sup> Standard fittings are 1/4" Brass. All high flow base plates are available in 1/4" Brass and Stainless steel fittings. To order, replace Part No. suffix into '-S4' for 1/4" Stainless steel fittings.

#### **Electronic Maintenance Indicator**

Part No.	Description	qty.
SGT-E0010	Electronic Maintenance Indicator Kit for Universal Ring Nut For use with Universal Ring Nut. Includes Electronic Indicator Device, Battery, USB Cable and Mounting clip.	ea.
SGT-E0040	Electronic Maintenance Indicator Kit for Super Clean Gas Filters For use with Super Clean Gas Filters. Includes Electronic Indicator Device, Battery, USB Cable and Mounting clip.	ea.
B0140	Universal Ring Nut	ea.

#### Replacement O-Rings for all Cartridge base plates

Part No.	Description	qty.
B0110	Base plate O-ring Replacement Set	20/pk.

#### Wall mounting Bracket for all Cartridge base plates

Part No.	Description	qty.
B0050	Wall-mounting Bracket Set	ea.

#### Base plate Flush Cap Replacement Set

Part No.	Description	qty.
B0130	Universal Flush Cap Replacement Set	2-pk.
	Compatible with normal and high flow base plates	

#### Base plate fittings

Part No.	Description	qty.
B0120	Connector Set - 1/4" Brass	6/pk
B0121	Connector Set - 1/8" Brass	6/pk
B0122	Connector Set - 1/4" Stainless steel	6/pk.
B0123	Connector Set - 1/8" Stainless steel	6/pk.

#### High Flow Base plate fittings

Part No.	Description	qty.
B0124	High Flow Connector Set - 1/4" Brass	6/pk.
B0125	High Flow Connector Set - 1/4" Stainless steel	6/pk.

#### Particle Filter for High flow Base plates

Part No.	Description	qty.
B0060	0.5 Micron Particle Filter $(1/4"$ Brass)	ea.
B0065	0.5 Micron Particle Filter Cup Replacement Pack	12/pk.





Management System ISO 9001:2008

www.tuv.com ID 9108620930

Distributed by:



Tallaght Business Park Whitestown, Dublin 24, Ireland D24 RFK3

Tel: (01) 4523432

Web: www.labunlimited.com

Quatro House, Frimley Road,

Camberley, United Kingdom **GU16 7ER** 

Tel: 08452 30 40 30

Fax: (01) 4523967 Fax: 08452 30 50 30 E-mail: info@labunlimited.com E-mail: info@labunlimited.co.uk

Web: www.labunlimited.co.uk