Data sheet

Saliva Sample Collection & stabilization Kit

Cat. No: SC001 Cat. No: SC002 Cat. No: SC003


Description

The Saliva sample collection & stabilization kit provide a safe and rapid all-in-one procedure for the collection, stabilization and transportation of saliva sample at ambient temperatures.

Advantages

- Easily self-collect. •
- Ship saliva samples at room temperature safely.
- The stabilized samples can be stored for over 2 years.
- Allows for subsequent isolation of DNA of excellent quality and integrity.

PROCEDURE:

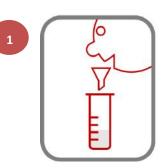

Saliva sample are collected spitting inside the bv collection funnel which has been assembled with the collection tube.



After collecting 2 ml of saliva the contents of saliva preservation solution are added and mixed with the collected saliva.




ITEM	001	002	003
SalivaCollection Tubes	20	100	250
Saliva Preservation Solution (2.5 mL)	20	100	250
Funnels	20	100	250

Kit Storage:



Store at room temperature.

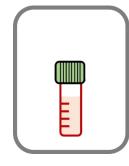

The saliva collection tube is sent to the laboratory for DNA isolation and analysis using **HigherPurity™ Buccal** Saliva Genomic DNA Extraction Kit (ref.AN0037).

Distributed by:



Tallaght Business Park Whitestown, Dublin 24, Ireland D24 RFK3

Tel: (01) 4523432 Fax: (01) 4523967

2

Quatro House, Frimley Road, Camberley, United Kingdom **GU16 7ER**

Tel: 08452 30 40 30 Fax: 08452 30 50 30 E-mail: info@labunlimited.com E-mail: info@labunlimited.co.uk Web: www.labunlimited.com Web: www.labunlimited.co.uk





Notes

PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively for research purposes and in vitro use only. The product was not tested for use in diagnostics or for drug development. Please refer to www.canvaxbiotech.com for Material Safety Data Sheet of the product.

Distributed by:



Tallaght Business Park Whitestown, Dublin 24, Ireland D24 RFK3

Tel: (01) 4523432 Fax: (01) 4523967 Web: www.labunlimited.com

Quatro House, Frimley Road, Camberley, United Kingdom **GU16 7ER**

Tel: 08452 30 40 30 Fax: 08452 30 50 30 E-mail: info@labunlimited.com E-mail: info@labunlimited.co.uk Web: www.labunlimited.co.uk

🛆 canvax