Distributed by:

**Lab Unlimited**  
CARL STUART GROUP

Tallaght Business Park  
Whitestown, Dublin 24,  
Ireland  
D24 RFK3

Tel: (01) 4523432  
Fax: (01) 4523967  
E-mail: [info@labunlimited.com](mailto:info@labunlimited.com)  
Web: [www.labunlimited.com](http://www.labunlimited.com)

Quatro House, Frimley Road,  
Camberley,  
United Kingdom  
GU16 7ER

Tel: 08452 30 40 30  
Fax: 08452 30 50 30

E-mail: [info@labunlimited.co.uk](mailto:info@labunlimited.co.uk)  
Web: [www.labunlimited.co.uk](http://www.labunlimited.co.uk)

## ULT / Low Temperature Freezers

GB Instructions for use  
DE Bedienungsanleitung  
FR Mode d'emploi

## Warning.

As the appliance contains a flammable refrigerant, it is essential to ensure that the refrigerant pipes are not damaged.

**The quantity and type of the refrigerant used in your appliance is indicated on the rating plate.**

Standard EN378 specifies that the room in which you install your appliance must have a volume of 1m<sup>3</sup> per 8 g of hydrocarbon refrigerant used in the appliances. This is to avoid the formation of flammable gas/air mixtures in the room where the appliance is located in the event of a leak in the refrigerant circuit.

### **WARNING:**

Keep ventilation openings in the appliance's cabinet or in the built-in structure clear of obstruction

### **WARNING:**

Do not use other **mechanical devices** or other means to accelerate the defrosting process than those recommended by the manufacturer

### **WARNING:**

Do not damage the refrigerant system

### **WARNING:**

Do not use **electrical appliances** inside the refrigerated storage compartment, unless they are of a type recommended by the manufacturer

### **WARNING:**

Do not expose the appliance to rain, and secure not splashing water when cleaning the floor

### **WARNING:**

This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack experience and knowledge, unless they have been given supervision or instructions concerning use of the appliance by a person responsible for their safety. Children should be supervised to ensure that they do not play with the appliance

**WARNING:**  
If a not compatible USB pen drive is used it can cause a reset of the controller

See page 18

**WARNING:**

Do not store explosives, such as aerosol cans with flammable propellants in the unit.

**WARNING:**

Danger risk of fire or explosion if flammable refrigerant are used. To be repaired only by trained personnel.

**WARNING:**

Sharp edges on cabinet, compressor compartment, evaporator, ventilation cover and on internal equipment can occur. Please be aware to avoid injury.

**WARNING:**

The condenser on the back of the appliance will in some cases have a hot surface. Please be aware to avoid injury.

- Always keep the keys in a separate place and out of reach of children
- Do not step on the lower panel to reach medicine in the top of the appliance.
- Before servicing or cleaning the appliance, unplug the appliance from the mains or disconnect the electrical power supply
- If the supply cord is damaged, it must be replaced by the manufacturer, its service agent, or similarly qualified persons in order to avoid a hazard
- **Relevant for Australia:** Supply cord fitted with a plug complies with AS/NZS 3112.
- Please note that changes to the appliance construction will cancel all warranty and product liability
- This device is intended to be used exclusively for medical products.
- If the instructions is lost please contact your supplier of the appliance to have a new instruction for use.

# Contents

- If service needed to this device, please be aware of only using service personnel with education in handling medical devices.

Warning .....2  
Unpacking the appliance ..... 4  
Get to know your freezer .....5  
Before use.....6  
Electrical connection..... 6  
Installation and start-up ..... 7  
Battery backup system at start-up ..... 10  
Interior fitting..... 11  
Port Hole..... 12  
Operation and function ..... 13  
Chart Recorder. Operating instruction ..22  
Defrosting, cleaning and maintenance .25  
Fault finding .....27  
Warranty, sparepart and service..... 28  
Disposal.....29

## Unpacking the appliance

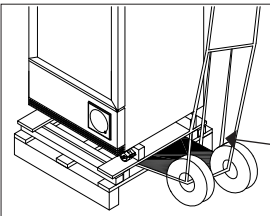
fig. 1

### Do not lay down the appliance!

Cut the band around the appliance and remove it.  
Remove the packing material of plastic and carton.  
The appliance is tipped from *from the side (handle side)*, by means of a sack truck.

Then the lower wooden frame can be removed.


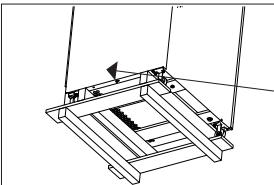
fig. 2

The lower wooden frame right under the appliance is bolted into the appliance.

Remove the 4 bolts using a wrench.

The appliance is tipped *from the side (handle side)*, by means of a sack truck. Then the wooden frame can be removed.

# Get to know your ULT/Low Temperature freezer
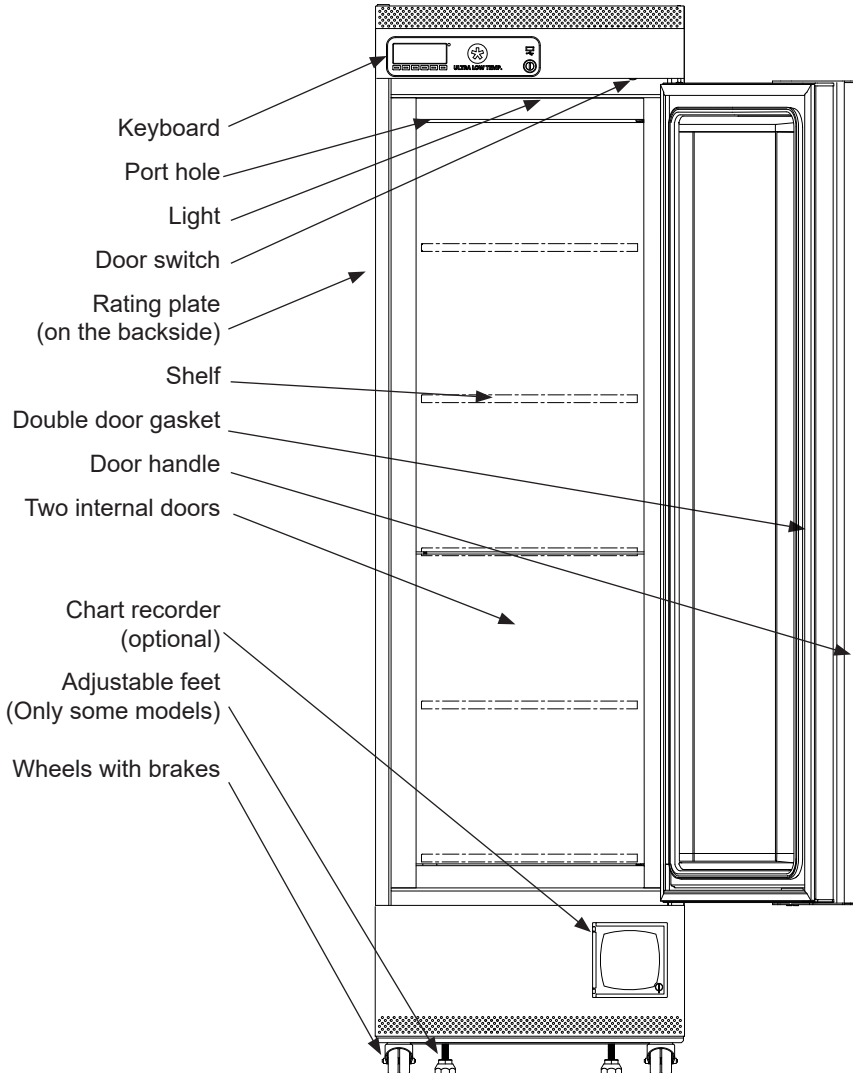
## Before use

Before operating your new appliance, please read the following instructions carefully as they contain important information on safety, installation, operation and maintenance. Keep the instructions for future reference.

On receipt, check to ensure that the appliance has not been damaged during transport. Transport damage should be reported to the local distributor before the appliance is put to use.

The foil on the shelves must be removed before cleaning and before using the unit. Clean the inside of the cabinet using warm water with a mild detergent. Rinse with clean water and dry thoroughly (see cleaning instructions). Use a soft cloth.

If during transport the appliance has been laid down, or if it has been stored in cold surroundings (colder than +5°C), it must be allowed to stabilise in an upright position for at least an hour before being switched on. Note: If the appliance is to be laid down, the door must face upwards and the appliance must be enclosed in the original packaging.

## Electrical connection

Wiring and connections in power supply systems must be all applicable (local and national) electrical codes. Consult these codes lengths and sizes prior to cabinet installation.

This device complies with relevant EU directives including Low Voltage Directive 2014/35/EU and Electromagnetic Compatibility Directive 2014/30/EU

The socket should be freely accessible.

Connect the appliance only to 220/240 V / 50Hz alternating current via a correctly installed earthed socket.

The socket must be fused with a 10 A or 13A fuse.

If the appliance is to be operated in a non-European country, check on the rating plate whether the indicated voltage and current type correspond to the values of your mains supply.

**Data regarding voltage and absorbed power / current are given on the rating plate.**

The power cord may be replaced by a technician only.

The rating plate provides various technical information as well as type and serial number.

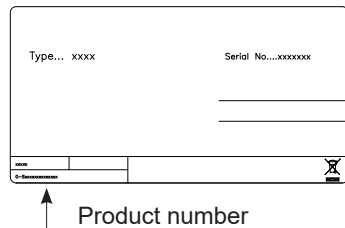
fig. 1

# Installation and start-up

## Placement

For safety and operational reasons, the appliance shall not be exposed to rain.

The appliance should be placed on a level surface in a dry, well ventilated room (max. 70% relative air humidity). Never place the appliance close to sources of heat and avoid placing it in direct sunlight.

The surface on which the appliance is to be placed must be level. Do not use a frame or similar.

The appliance can be installed as a free-standing unit against a wall or side-by-side with other appliances (see fig. 5 on page 7)

## Room temperature

The climate class is stated on the name plate. This specifies the optimum room temperature.

The appliance operates best at room temperatures:

Model	Optimum room temperature
ULT U250 -86°C	+16 °C to +20 °C
XLT U250 -60°C	+16 °C to +25 °C
LT U250 -45°C	+18 °C to +25 °C

## Ventilation

Keep ventilation openings, in the appliance enclosure or in the built-in structure, clear of obstruction. The figures below illustrate the clearance necessary to achieve the required air circulation. Fig 2 (Minimum 150 mm rear)

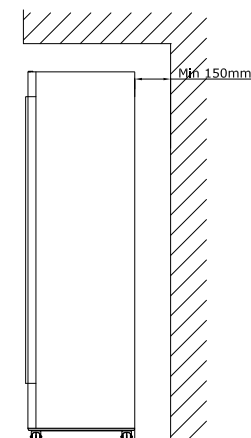
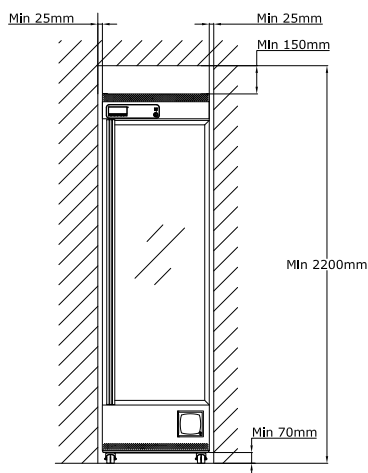
fig. 2

## Moving the unit

Unlock brakes on wheels.

Before moving the unit, ensure that the door is closed and locked. When moving the unit, always face the door and take a firm grip on either side of the door as shown. Fig 3

When the unit is installed, lock all the wheels (those that can be locked) and keep them locked until the unit shall be moved.

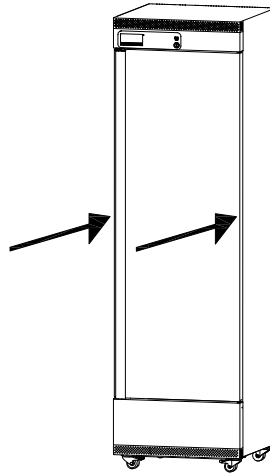
fig. 3

## Levelling the appliance

It is important that the appliance absolutely level.

If the appliance is to be placed on a soft surface, e.g. floorboards or a carpet, it is best to recheck whether the appliance is still level after a period of time as the underlying surface may yield under the weight of the appliance.

Some models have adjustment feet.  
(See fig 4)

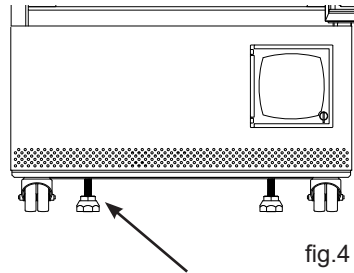
fig.4

## Side by side

In side-by-side arrangement, there must be at least 25 mm between and at the sides of the cabinets, so the doors can be opened freely (See fig. 5)

## Daily use

Please always wear gloves due to the very low temperatures inside the unit.

Always thoroughly remove defrosted ice/water on the front of the cabinet before closing the doors to prevent the gasket from freezing the door and cabinet together.

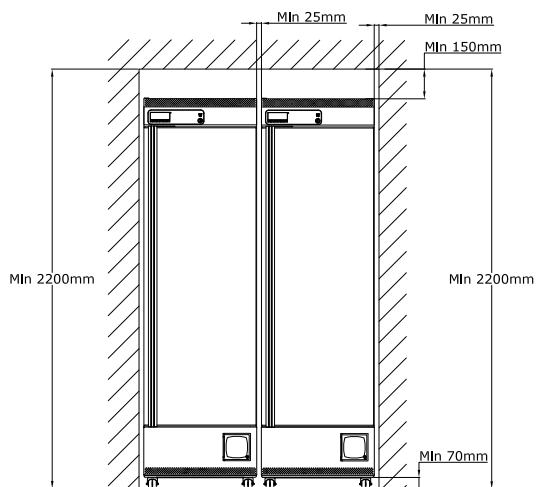
fig.5

## Installation of ventilation guard

The guards function is to secure an absolute minimum of ventilation, if the appliance is pushed too close up against a wall.

There are two different variants.

Fig 6.1 or 6.2

fig. 6.1

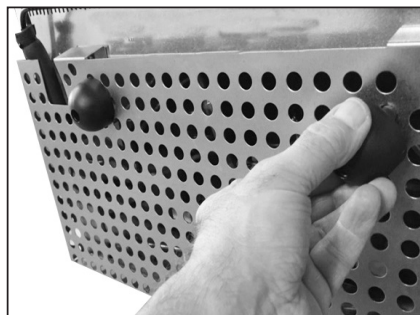
fig. 6.2

## Battery backup system at start-up

Some models are equipped with a battery back up system.

1. Open the door
2. Use a pen to switch on the battery backup system in the hole shown on fig. 7
3. When turned on the switch it will have a green light

### NOTE!

The battery back up system does not supply the cooling system with power to run.

When starting up the appliance for the first time it is necessary to switch on the battery back up system.

### NOTE!

The battery for back up should be changed every third year to secure 48 hours of back up. Please put this change in the maintenance schedule for every third year.

## Battery backup system

Some models are equipped with a battery back up system. The system supplies the controller and keyboard with power at power failure. This makes it possible to supervise the temperatures in the unit during the power failure.

The battery backup system makes it possible to supervise the temperatures for 50 hours. After a power failure and at the first start up the battery needs to be recharged. To regain the full capacity the battery will be reloading for 10 days

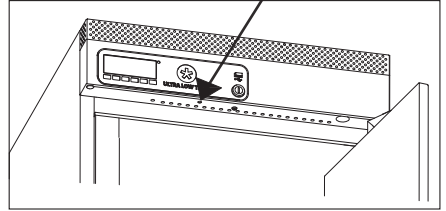
fig. 7

# Interior fitting

## Shelves

The shelves can be repositioned. To remove the shelves the door must be opened at least 90°.

1. Remove shelf brackets from shelf rails.
2. Replace shelf brackets in desired position (See fig. 8)




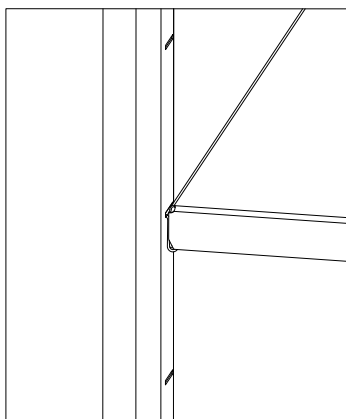
fig.8

The foil on the shelves must be removed before cleaning and before using the unit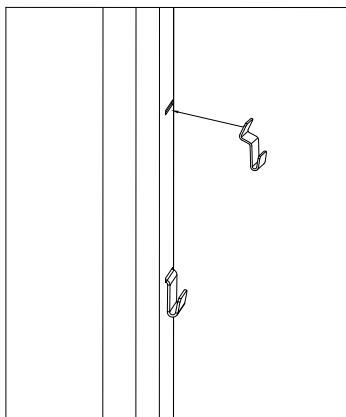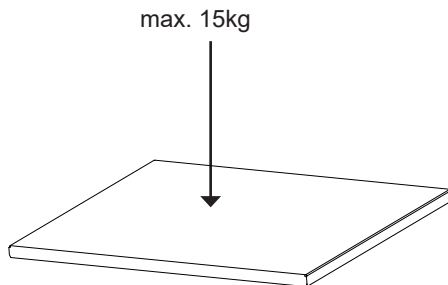


## Port Hole


Some models are equipped with an extra port hole. Ex. for use of an external sensor.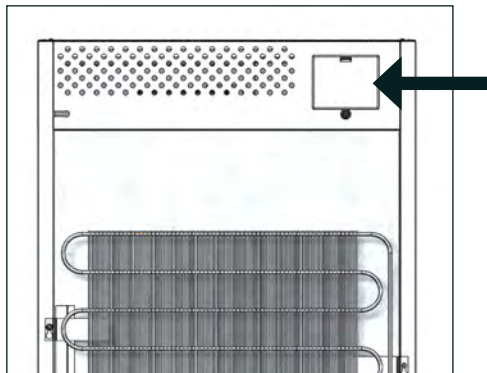

1. Open the service hatchway in the top on the back. Use a screwdriver.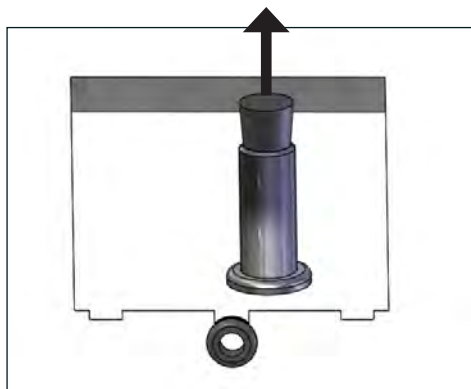

2. Remove the plug in the port hole tube.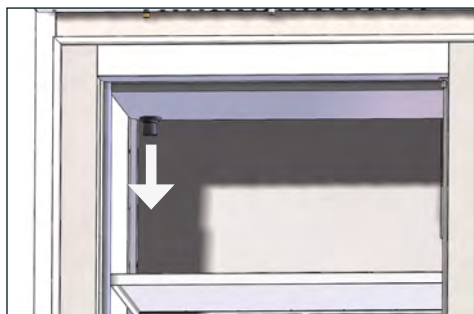

3. Remove the plug in the port hole inside the freezer.

4. Insert the sensor in the port hole.

5. Close the hole in the tube with putty – both inside and outside.

7. Close the hatchway

## Operation and function

### Start up:


When the appliance is connected to the power supply, the keyboard will automatically start up.

The start up view on the keyboard will show the different software installed on the controller of the appliance.


Press enter to return from the start up view.

### Operation – main view: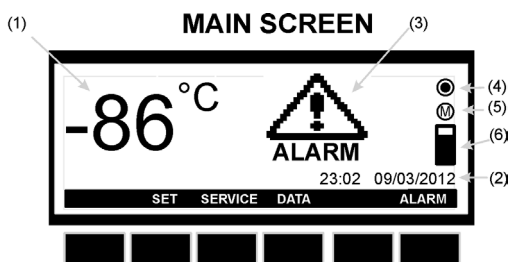

- (1) The temperature in the appliance (measured by the TR3 probe)
- (2) Time and date
- (3) ALARM icon. Flashing by alarm. On when there has been an alarm, but the alarm is no longer active.
- (4) Logging icon. Of if no logging. On when logging
- (5) Memory icon. On when the memory is 90% full. Flashing when the memory is full, and the controller is deleting the oldest logging data.
- (6) Memory bar. Shows the status of the memory.


Keys – main view: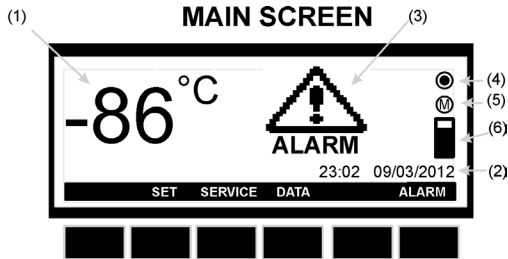

- SET:** Enter the Set Point menu of the temperature
- SERVICE:** Enter the Service menu
- DATA:** Enter the Data Logging menu
- ALARM:** Enter the Alarm menu


How to see and modify the temperature Set Point:

1. Push and immediately release the **SET** key: the display will show the Set Point value.
2. To modify the value push the **SET** key, the Set Point start flashing.
3. Use the **UP and DOWN** keys to modify the value.
4. To memorize the new Set Point value push the **SET** key again or wait 30sec.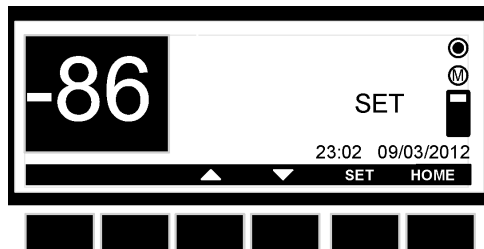

## Service menu:


From the main view push the **SERVICE** key and the SERVICE menu is entered.  
See below picture: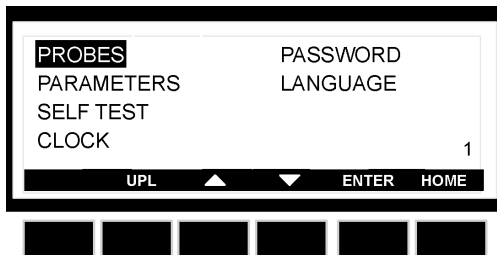

- PROBES:** Enter the probes, to see the measured temperatures. 1-4 probes is available depending on model.
- PARAMETERS:** Enter the setting of the parameters. Please note that changes made to the parameters should only be made of a technician. The code to enter the Pr2 parameters is standard set to 12.
- SELF TEST:** Enter the Self Test program of the controller.
- CLOCK:** Enter clock menu, where it is possible to change date and time.
- PASSWORD:** Enter the password menu, where it is possible to change the password.
- LANGUAGE:** Enter the language menu, where it is possible to change language.

## How to set time and date:


1. Enter the SERVICE menu
2. Select CLOCK sub-menu
3. Push the ENTER key.
4. Set the day by means of the UP and DOWN keys.
5. Push the SET key, to confirm and pass to the setting of time.
6. Use the same procedure as for the date.
7. Then confirm the selection by means of the SET key.

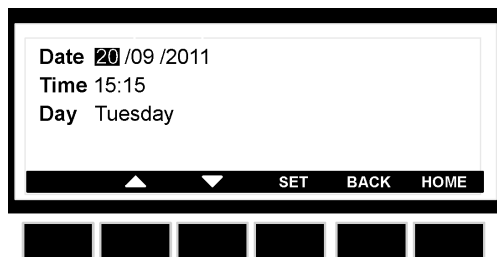
## How to use the Self Test:

1. Enter the **SERVICE** menu
2. Select **SELF TEST** sub-menu
3. Push the **ENTER** key,
4. If **PASSWORD** is required, insert the password,
5. OTHERWISE the SELF TEST menu is entered directly.
6. Push the **START** key
7. Then push the keys to activate the correspondent loads:
  - a. Compressor
  - b. Light
  - c. Fan
  - d. Alarm
  - e. Buzzer
8. The display shows the status of the digital input (ON or OFF)
9. Wait 30s or push the **BACK** key to come back to the previous screen.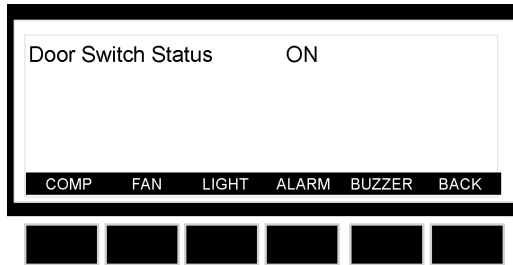

## How to change language:

1. Enter the **SERVICE** menu
2. Select **LANGUAGE** sub-menu
3. Push the **ENTER** key and the **LANGUAGE** menu is entered.
4. Push the **SET** key and then use the **UP** and **DOWN** keys to select the language and then the **SET** key to confirm it.

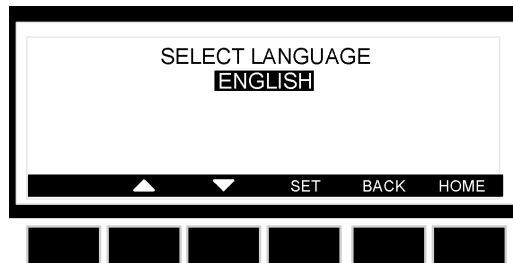## Data menu:

From the main view push the **DATA** key and the DATA menu is entered.  
See below picture:

**LOG:** Enter the data logged by the controller.


**EXPORT:** Export the data to an USB pen drive delivered with the appliance.

**GRAPH:** Enter the graph showing the temperature logged the last 24h (with a logging interval of 15min.)

## How to enter the log:

1. Enter the **DATA** menu
2. Select **LOG** sub-menu
3. Push the **ENTER** key and the LOG menu is entered.
4. By **UP and DOWN** keys chose the data interval to display
5. Push the **ENTER** key to display the selected data.

**NOTE: THE  KEY: IS USED TO STOP AND START LOGGING**

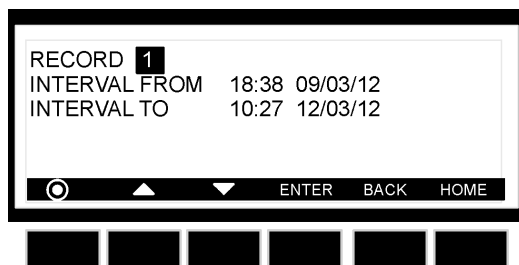Logged data will have this layout:

		TR3	EVP	TL1	TL2	ST
19/05/08	11.34	25.4	25.8	-	25.5	C
19/05/08	11.35	25.4	25.8	-	25.5	C
19/05/08	11.36	25.4	25.8	-	25.5	-
19/05/08	11.37	25.4	25.8	-	25.5	D

EXIT

Where TR3, EVP, TL1, TL2 = Value of probes.

With probe failure or absence:” - “ symbol is displayed. Please note that 1-4 probes are available depending on model.

ST: status of the controller/load

- - = operating, without any load activated;
- D = defrost running (if automatic defrost is available)
- C = compressor working

## How to export data:


1. Enter the **EXPORT** menu
2. Insert the USB pen drive supplied by **Nordic Freeze**.
3. Select **ALARM** or **DATA**, the controller starts sending data to the pen drive, when the export is finished the message: **EXPORT - Copy completed** is displayed.

The exported data will be exported as a CSV-file (Comma Separated Values). This file can be used in ex. Excel for making graphs.

**IMPORTANT:** during the download don't remove the USB pen drive: this action could damage the data files and USB pen drive itself.

**WARNING:** leave the USB pen inserted only for the time necessary to export data then remove.


**WARNING:** if a not compatible USB pen drive is used it can cause a reset of the controller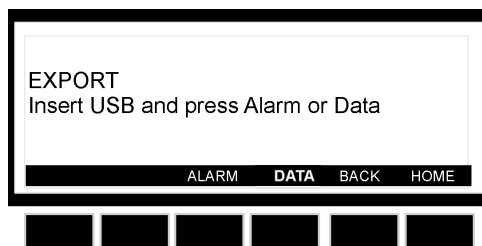

## How to enter the Graph:

1. Enter the DATA menu
2. Select GRAPH sub-menu
3. Push the ENTER key and the GRAPH menu is entered.
4. By UP and DOWN keys chose the probe that has to be displayed.
5. Push the HOME key to get back to the main view.

**NOTE: A graph is erased when the controller is switched off.**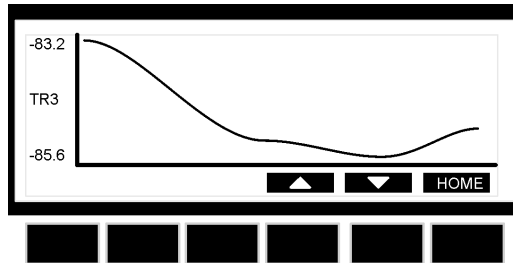

## Alarm menu:


If the alarm icon is **flashing** on the main display, an alarm is occurring.

If the alarm icon is **displayed but not flashing** on the main display, an alarm is occurred and recovered.

Once the alarm signal is detected the buzzer can be silenced by pressing any key.

### Active alarms:

1. Push the **ALARM** key to enter the alarm menu.
2. The alarm menu displays the active alarm with the following layout:
  - a. **First column** = alarm code
  - b. **Second column** = alarm description

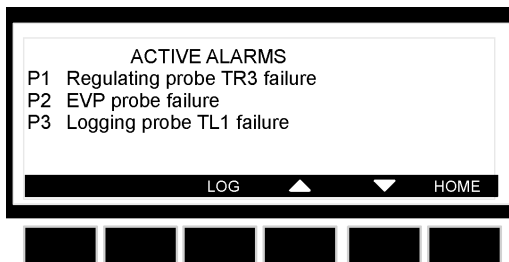

3. Push the **LOG** button to enter the **ACTIVE ALARM LOG**.

This menu contains all the information concerning the active alarms. In the first line, it is displayed how many alarms are happening.

4. It's possible to move through the alarms by the UP and DOWN keys.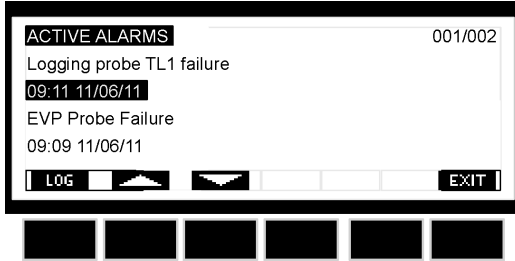

5. Push the **LOG** button to enter the **ALARM LOG**. This menu contains all the memorized alarms. For each alarm the starting time and date and the finish time and date are recorded.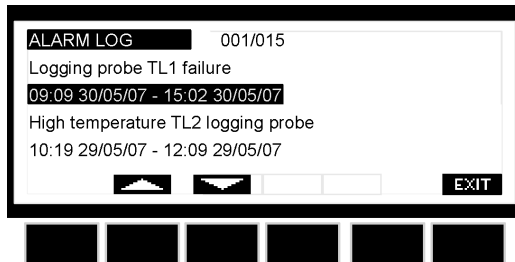

"P1"	Regulating probe TR3 failure	Alarm output ON; Compressor output according to parameters <b>Con</b> and <b>CoF</b>
"P2"	EVP Probe Failure	Alarm output ON; Other outputs unchanged
"P3"	Logging probe TL1 failure	Alarm output ON; Other outputs unchanged
"P4"	Logging probe TL2 failure	Alarm output ON; Other outputs unchanged
"HA1"	TR3 High Alarm	Alarm output ON; Other outputs unchanged
"LA1"	TR3 Low Alarm	Alarm output ON; Other outputs unchanged
"HA3"	High temperature alarm probe TL1	Alarm output ON; Other outputs unchanged
"LA3"	Low temperature alarm probe TL1	Alarm output ON; Other outputs unchanged
"HA4"	High temperature alarm probe TL2	Alarm output ON; Other outputs unchanged
"LA4"	Low temperature alarm probe TL2	Alarm output ON; Other outputs unchanged
"dA"	Door Open Alarm	Compressor and fans depend on " <b>rrd</b> "
"EA"	External Alarm	Output unchanged
"CA"	Serious Alarm	All outputs OFF
	"Real Time Clock Error".	Please set the date and time. See page 15

# Chart Recorder

## Operating instruction

### Maintenance:



- Regularly replace the chart and the fibre point pen.
- Replace the alkaline battery every year.

### Putting into operation

1. Open the door with the key.
2. Lift the steel pen (3) into upright position away from the chart (7).
3. Push the chart holder (6) sideways and lift it up, and remove the chart (7)
4. Remove the plastic strip to activate the battery.
5. Reinstall the chart (7), lower the chart holder (6), and push it sideways.
6. Lower the steel pen carefully (3).
7. Turn the chart manually in correct starting point and note the date on the chart.


**It is advisable to wait almost 6 hours to see the trail of the fibre point pen on the chart.**

8. Remove the protection of the pen-nib.
9. Close the door with the key.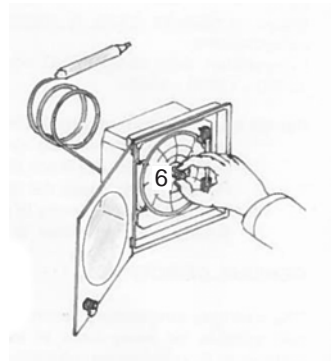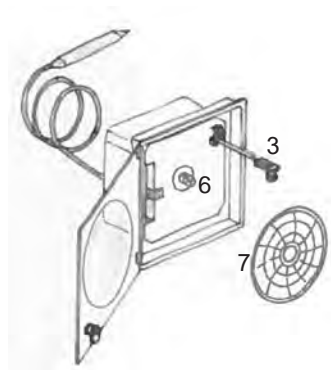


## Substituting the chart

1. Push the chart holder (6) sideways and lift it up.
2. Remove the chart (7).
3. Insert the new chart.
4. Lower the chart holder (6) and push it sideways.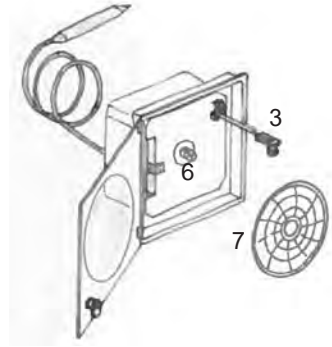

## Substituting the fiber tipped pen

1. Pull out the fibre pen (4) taking care not to damage the pen.
2. Gently place the new fibre pen on the pen arm.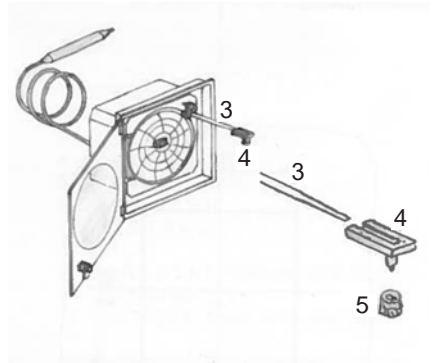

## Controll working - Pen arm calibration


The chart recorder is calibrated from the factory. However, it is important to check that the temperature trail is showing the right temperature.

We suggest to control the recorder's calibration every 12 months.

Messure the temperature near by the installed temperature sensor (probe) with a reliable thermometer.

If there is a difference of  $\pm 2^{\circ}\text{C}$  between the data provided by the reliable thermometer and those of the recorder, proceed as follows:

- Gently take the pen's arm B.
- With a small screwdriver, calibrate the regulation screw A, until the fibre pen C is writing the same temperature as the reliable thermometer shows.

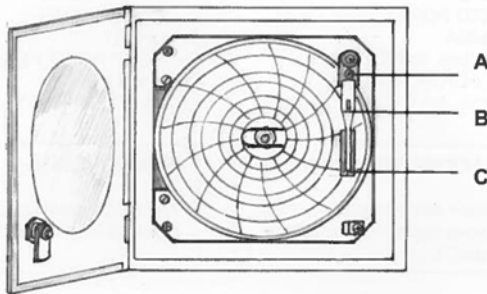# Defrosting, cleaning and maintenance

## Defrosting:

The items stored in the freezer and the humidity of the surrounding air gradually cause frost and ice to form inside the freezer. A thin layer of frost or ice does not affect freezer performance, but the thicker the layer, the higher the energy consumption.

Never use electrical apparatus or sharp implements as these may damage the inside of the appliance.

## Quick defrosting:

Small amounts of loose frost can be removed using a plastic scraper.

Defrosting has to be done regularly to prevent ice building.

The ice build depends on the use of the freezer. The more daily door openings – the more manual defrosting.

Door open/ close per 24h	Advised monthly manual defrosting
<5	1-2
<10	2-4
<15	4-6

## Thoroughly defrosting:

When the layer of frost and ice exceeds 4-5 mm the freezer should be completely defrosted. It is recommended that defrosting be performed when there are no items stored in the freezer.

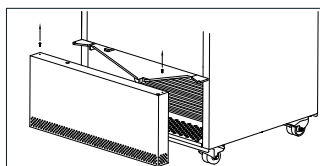
fig. 9

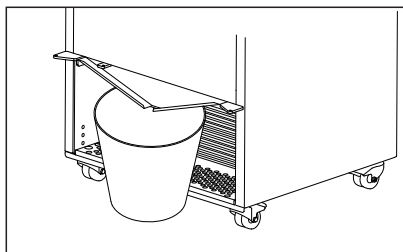
fig. 10

## Proceed as follows:

1. Disconnect the power supply.
2. Remove the lower panel from beneath the door by loosening the two screws. Use a 2mm Hexagon key. See fig 9
3. Position a bucket or a suitable container under the defrost water drain. /see fig 10)
4. Place a bowl of hot (but not boiling) water in the freezer, close the door and wait approx. 30 minutes before removing loose frost with the plastic scraper.
5. To prevent water from damaging the floor during defrosting, it is advisable to check the bucket/container from time to time and to cover the floor around the defrost water drain to protect against ice and splashes. When all the frost and ice have melted, clean the inside of the appliance using water with a mild, perfume-free detergent. Dry the freezer with a soft cloth.
6. Mounts the lower panels again and reconnect the power supply.
7. Return items to the freezer when the required temperature has been reached.

## Cleaning and maintenance

Before cleaning, disconnect the power supply and remove all loose fittings.

The appliance is best cleaned using warm water (max. 65°C) with a little mild, perfume-free detergent. Never use cleaning agents that scour. Use a soft cloth. Rinse with clean water and dry thoroughly. It is important to prevent water from entering the control panel.

The sealing strips around the door must be cleaned regularly to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strip, check that it continues to provide a tight seal.

It is recommended that the wire and tubular condensers on the rear of the appliance be cleaned.

The ventilation grille, the condenser, the compressor and the compressor compartment must also be kept free of dust and dirt. Clean it with a vacuum cleaner regularly. Remove the lower panel from beneath the door by loosening the two screws (see fig 9) Do the same on the back (see fig. 11)

**OBS:** If the appliance is not to be used for any length of time, switch it off, disconnect the power supply, empty it, clean the inside and leave the door open to allow air circulation and prevent smells.

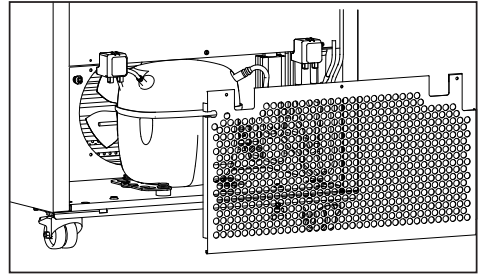
fig. 11

## Fault finding

Fault	Possible cause	Remedy
Screen of keyboard is not lit.	Power failure; the fuse is blown; the appliance is not plugged in correctly, the power is cut.	Check that power is connected. Reset the fuse
Temperature in freezer too high	<p>The ventilation grille is blocked.</p> <p>The door is not closed properly.</p> <p>The quantity of items placed in the freezer at a time is greater than the capacity of the appliance.</p> <p>Thick layer of frost and ice.</p> <p>The temperature setting is too high</p>	<p>Ensure unhindered air circulation</p> <p>Close the door and wait a few hours</p> <p>Wait a few hours</p> <p>Defrost the freezer.</p> <p>Lower the temperature setting.</p>
Temperature in freezer too low.	The temperature setting is too low	Raise the temperature setting
Vibration or bothersome noise	The appliance is not level.	Level the appliance (For model with adjustable feet: use a spirit level)
Compressor runs continuously.	<p>The temperature setting is too low.</p> <p>Too high room temperature.</p>	<p>Raise the temperature setting.</p> <p>Ensure adequate ventilation</p>
Display or alarm log shows "Real Time Clock Error".	The date and time is not set.	Please set the date and time See page 15

# Warranty, spare parts and service

## Warranty disclaimer

Faults and damage caused directly or indirectly by incorrect operation, misuse, insufficient maintenance, incorrect building, installation or mains connection. Fire, accident, lightning, voltage variation or other electrical interference, including defective fuses or faults in mains installations.

Repairs performed by others than approved service centres and any other faults and damage that the manufacturer can substantiate are caused by reasons other than manufacturing or material faults are not covered by the warranty.


Please note that changes to the construction of the appliance or changes to the component equipment of the appliance will invalidate warranty and product liability, and the appliance cannot be used lawfully. The approval stated on rating plate will also be invalidated.

Transport damage discovered by the buyer is primarily a matter to be settled between the buyer and the distributor, i.e. the distributor must ensure that such complaints are resolved to the buyer's satisfaction.

Before calling for technical assistance, please check whether you are able to rectify the fault yourself. If your request for assistance is unwarranted, e.g. if the appliance has failed as a result of a blown fuse or incorrect operation, you will be charged the costs incurred by your call for technical assistance.

## Spare parts

When ordering spare parts, please state the type, serial and product numbers of your appliance. This information is given on the rating plate. The rating plate contains various technical information, including product-type and serial numbers.

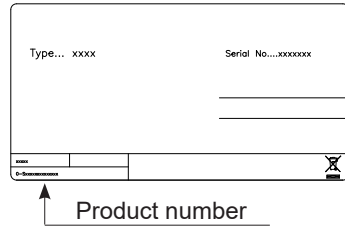# Disposal

## *Information for Users on Collection and Disposal of Old Equipment and used Batteries*

These symbols on the products, packaging, and/or accompanying documents mean that used electrical and electronic products and batteries should not be mixed with general household waste. For proper treatment, recovery and recycling of old products and used batteries, please take them to applicable collection points, in accordance with your national legislation and the Directives 2002/96/EC and 2006/66/EC.

By disposing of these products and batteries correctly, you will help to save valuable resources and prevent any potential negative effects on human health and the environment which could otherwise arise from inappropriate waste handling.

For more information about collection and recycling of old products and batteries, please contact your local municipality, your waste disposal service or the point of sale where you purchased the items.

Penalties may be applicable for incorrect disposal of this waste, in accordance with national legislation.

## **For business users in the European Union.**

If you wish to discard electrical and electronic equipment, please contact your dealer or supplier for further information.

## **Information on Disposal in other Countries outside the European Union**

These symbols are only valid in the European Union. If you wish to discard this product, please contact your local authorities or dealer and ask for the correct method of disposal.

### **Note for the battery symbol:**

This symbol might be used in combination with a chemical symbol. In this case it complies with the requirement set by the Directive for the chemical involved.

### **NOTE:**

It is very important to dismantle the door before disposal. This to prevent any accident by children playing with the cabinet and door.

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